

Advanced sample preparation by broad ion beam milling for EBSD analyses





Dr. Laurie Palasse



Senior Application Scientist EBSD
Bruker Nano, Berlin

Dr. Pawel Nowakowski

Applications Scientist
Fischione Instruments, USA

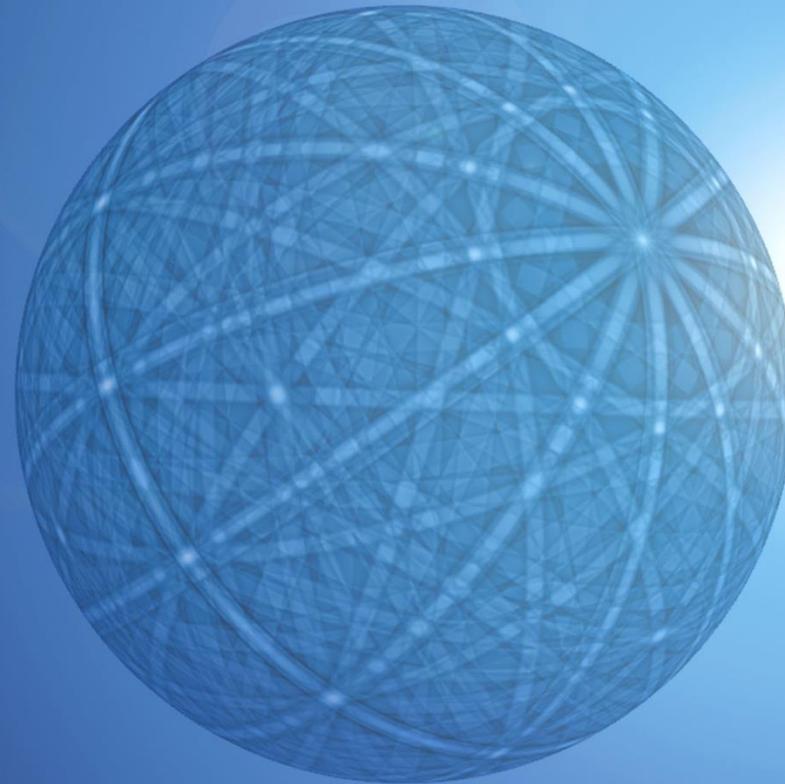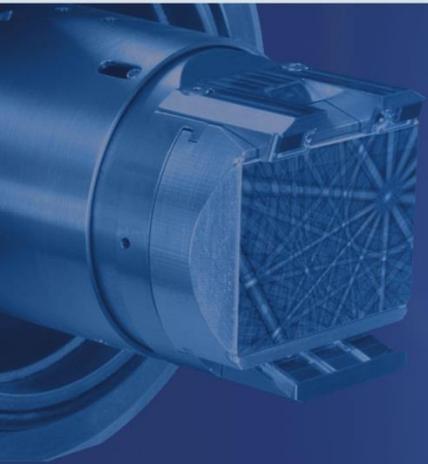

EBSD

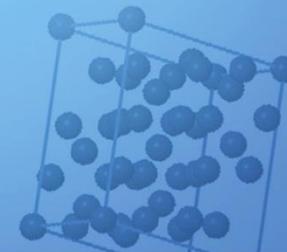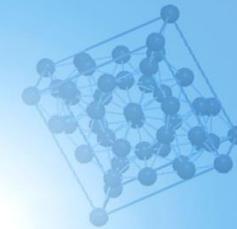

Outline



- EBSD technique overview
- ARGUS™ imaging system
- Broad ion beam milling: Model 1061 SEM Mill
- Application examples:
 - › Austenitic steel
 - › Zirconium alloy
 - › Ti CP
 - › Ti alloy
 - › Solder bumps

EBSD technique



Introduction

What is EBSD?



- **EBSD** - Electron **B**ack**S**catter **D**iffraction
- SEM based technique to measure **crystal orientations**
- Applicable to any **crystalline material** (in theory)
- Provides the absolute crystal orientation with sub-micron spatial resolution:
as fine as 300 nm grain size for bulk sample, 10 nm with Transmission Kikuchi Diffraction (TKD)
- Complementary technique between X-ray diffraction and TEM studies

➔ *The only technique that correlates local texture (orientation) with microstructure (grain metrics, grain boundary, ...)*

Introduction

EBSD: What is it used for?




To understand the material properties by:

- Analysing the grain structure, crystal orientation, phase ID, and distribution
- Residual stress, quality control, phase transformation, fracture analysis, strain analysis, tectonophysics studies...
- **Texture analysis:** Assessing the effects of the thermomechanical processes
- Complement the chemical information; **maximise your knowledge of the sample** and the processes involved

EBSD technique


EBSP formation

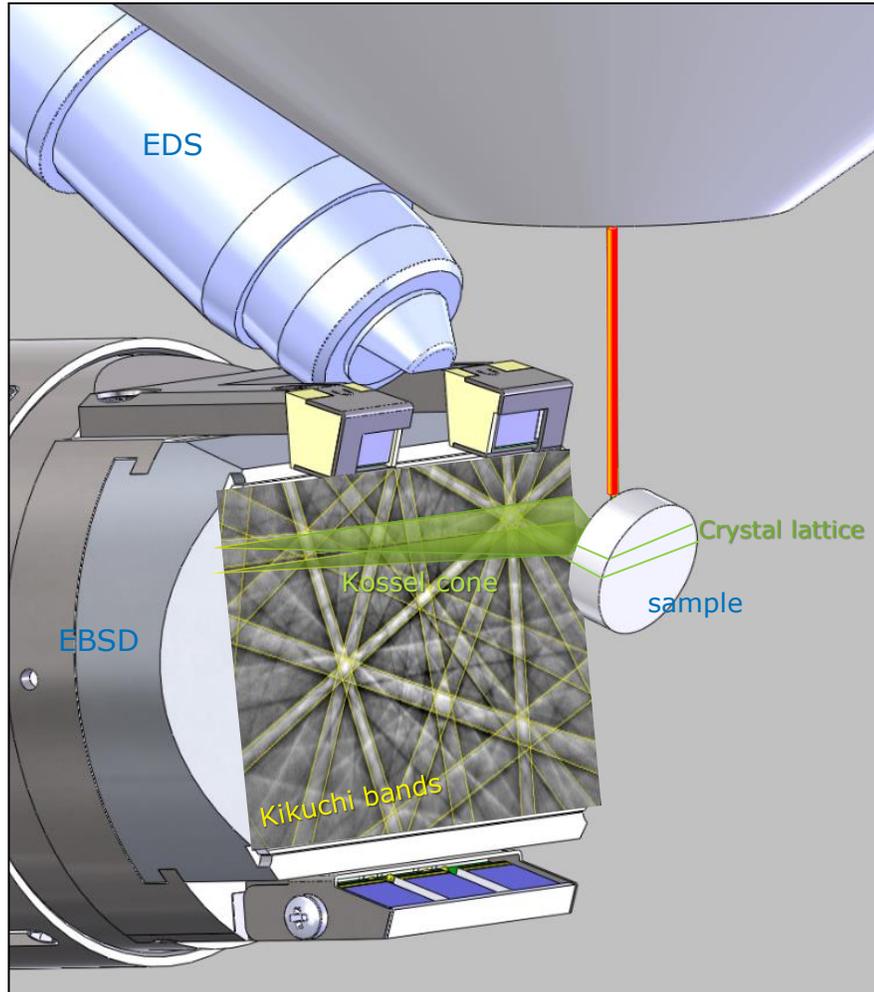

- e- beam strikes specimen
- Scattering produces e- travelling in all directions
- e⁻ that satisfy the Bragg condition ($n\lambda=2d.\sin\theta$) for a plane (hkl) are channeled \Rightarrow 2 hyperbolas/diffraction cones (**Kikuchi bands**) corresponding to the various diffracting planes
- Backscatter volume: Top few nanometers below the surface
- e⁻ strike the phosphor screen and produce light (gnomonic projection)
- Which is detected by a CCD or CMOS camera and digitised
- The resulting diffraction pattern "**EBSP**" is automatically analysed and indexed...

EBSD technique

EBSD measurement phase identification & distribution

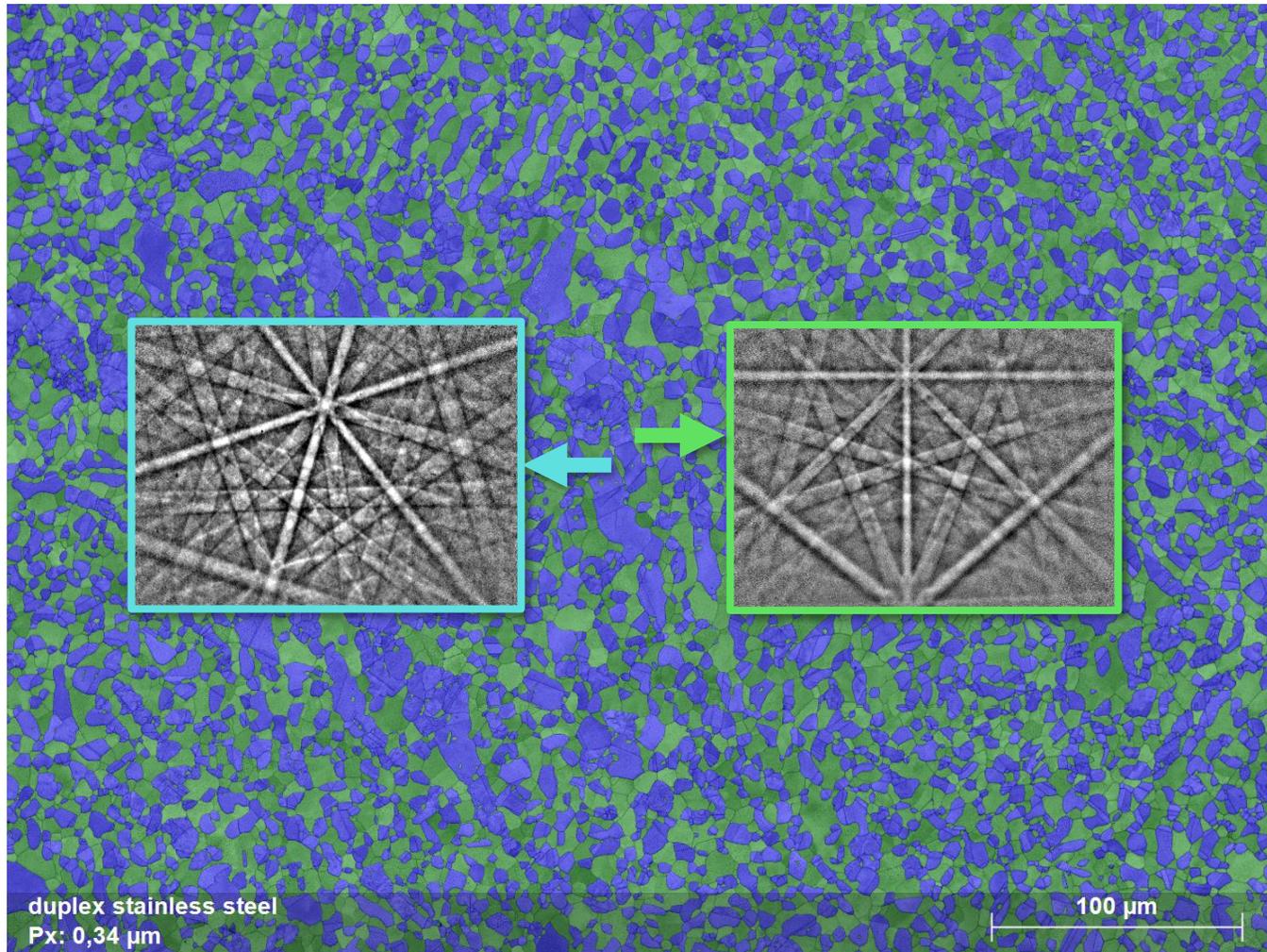


EBSD phase map

Hit rate 99,4%

	Fe fcc	49,1%
	Fe bcc	50,2%

EBSD technique

EBSD measurement phase identification & distribution

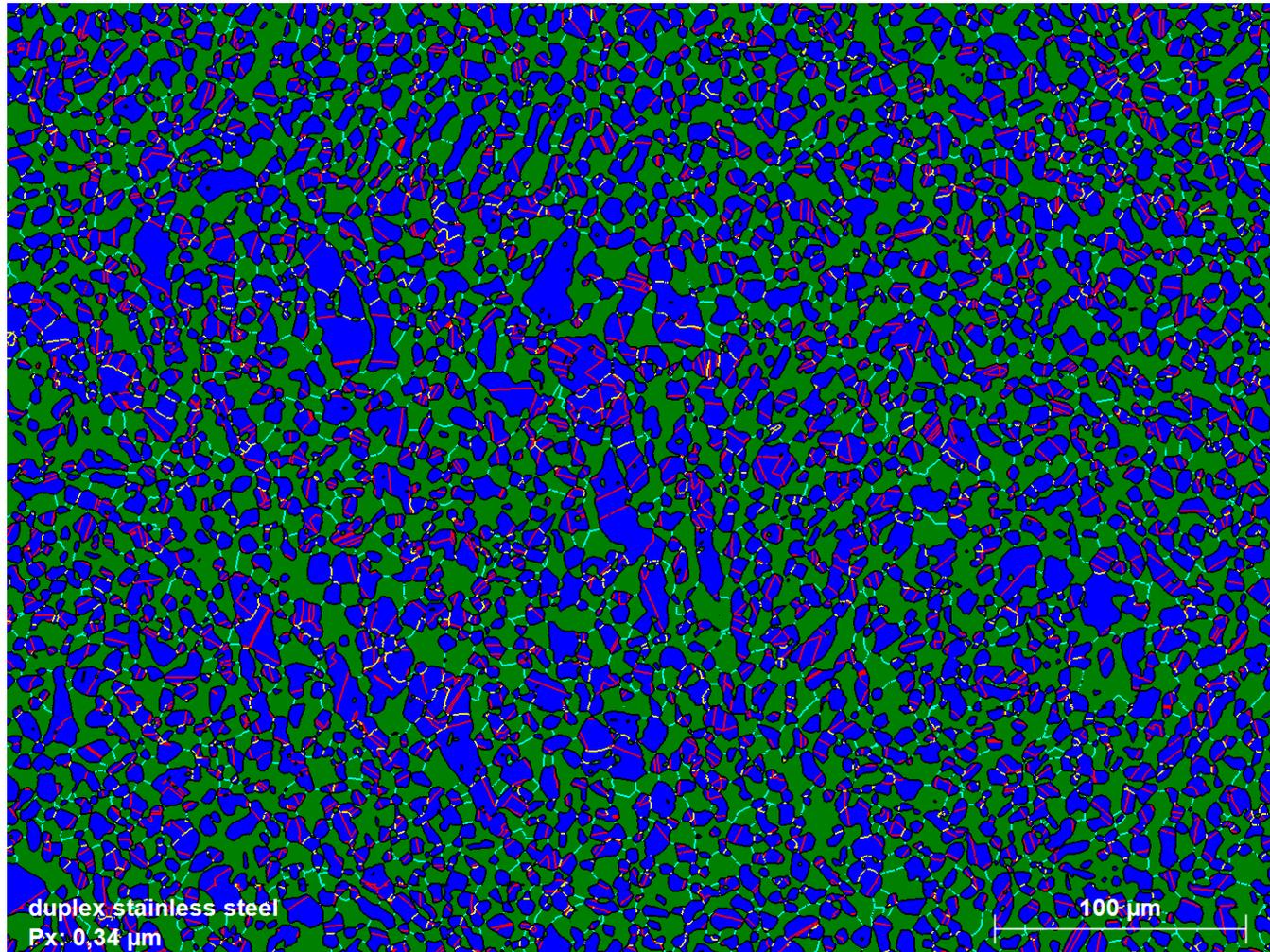


EBSD phase map

Hit rate 99,4%

- Fe fcc 49,1%
- Fe bcc 50,2%
- Grain boundaries (13,1 mm)
- Phase boundaries (63,4 mm)
- $\Sigma 3$ (17 mm)



EBSD technique

EBSD measurement – grain metrics

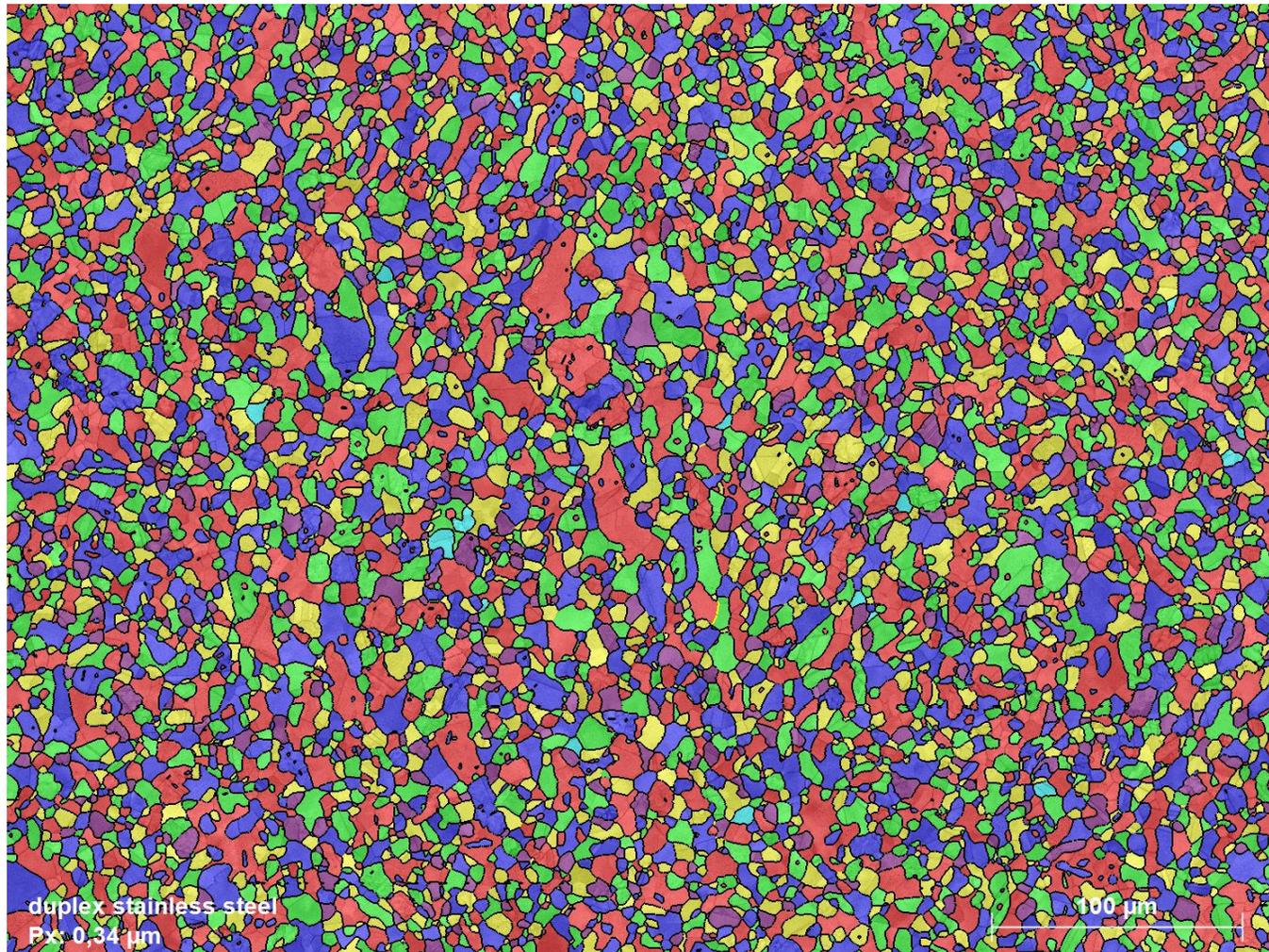

Grain distribution map

1,7M points in 30 min
5631 grains measured
Average grain size: 11 μm

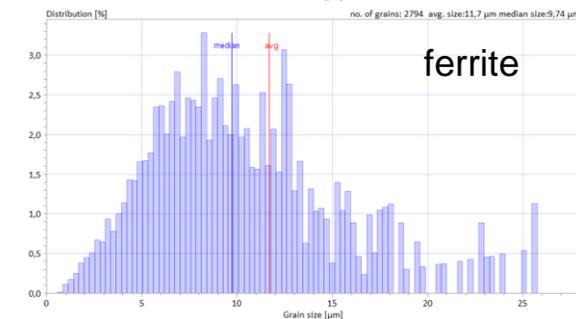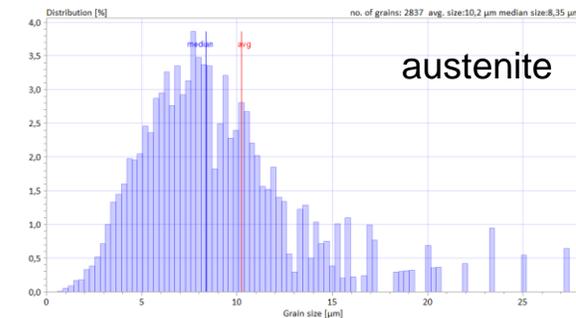


Correlation between local texture, microstructure, grain boundary and grain size analysis, ...

EBSD technique


EBSD measurement – other applications





Transmission Kikuchi Diffraction

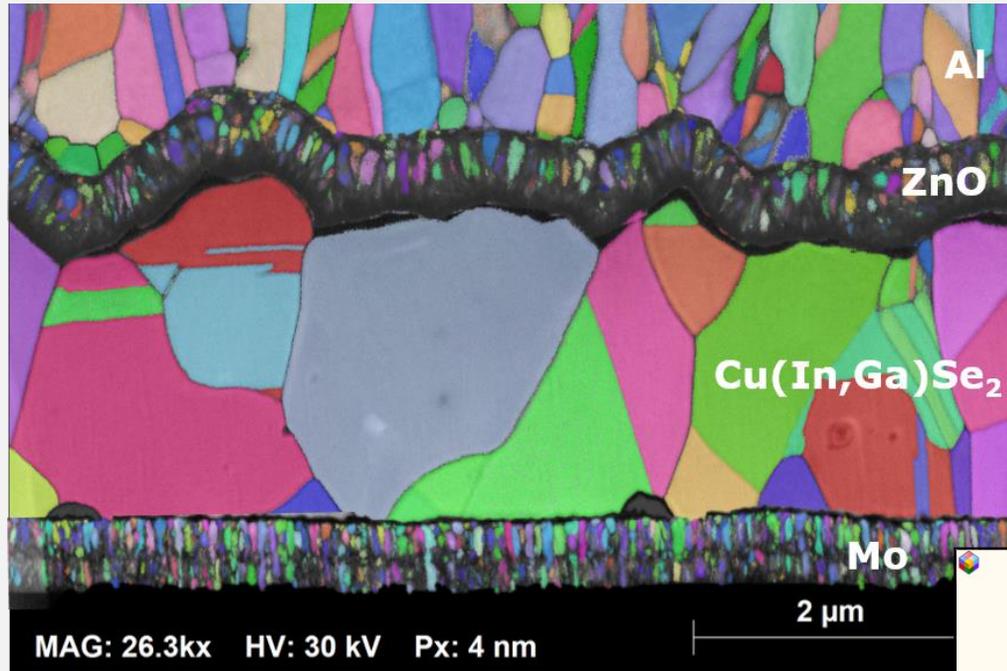

In situ EBSD/EDS & mechanical testing

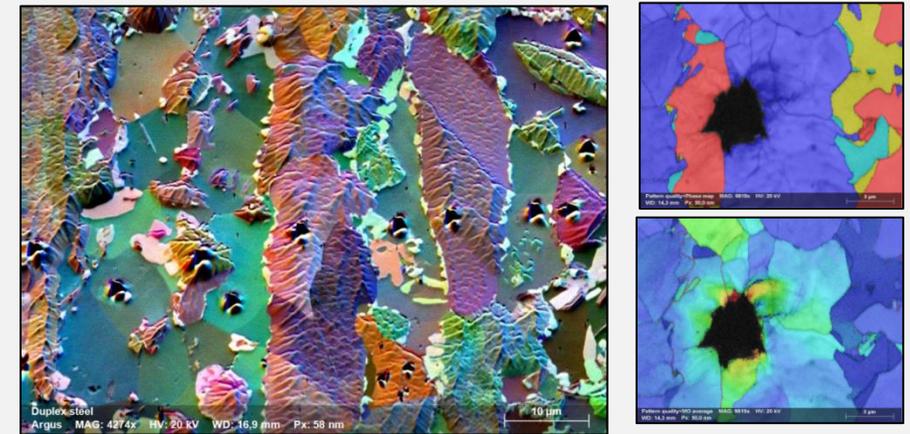



3D EBSD/EDS

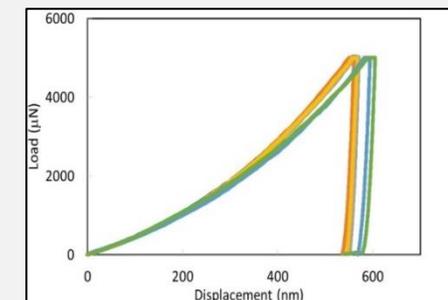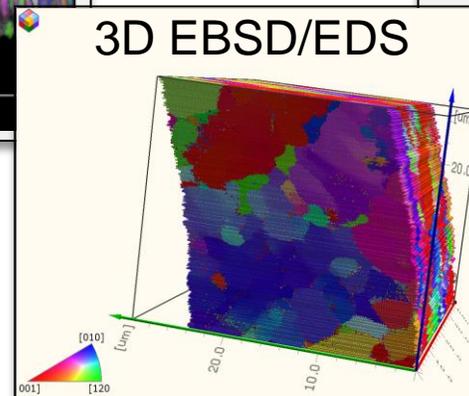


High-quality EBSD measurement requires high-quality sample preparation



Highly deformed copper sample prepared by ion milling

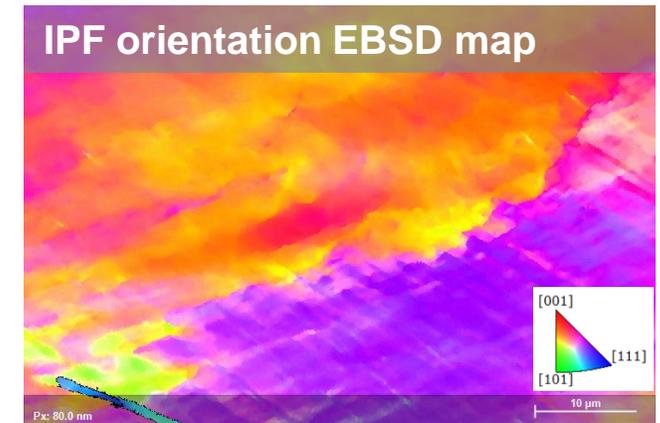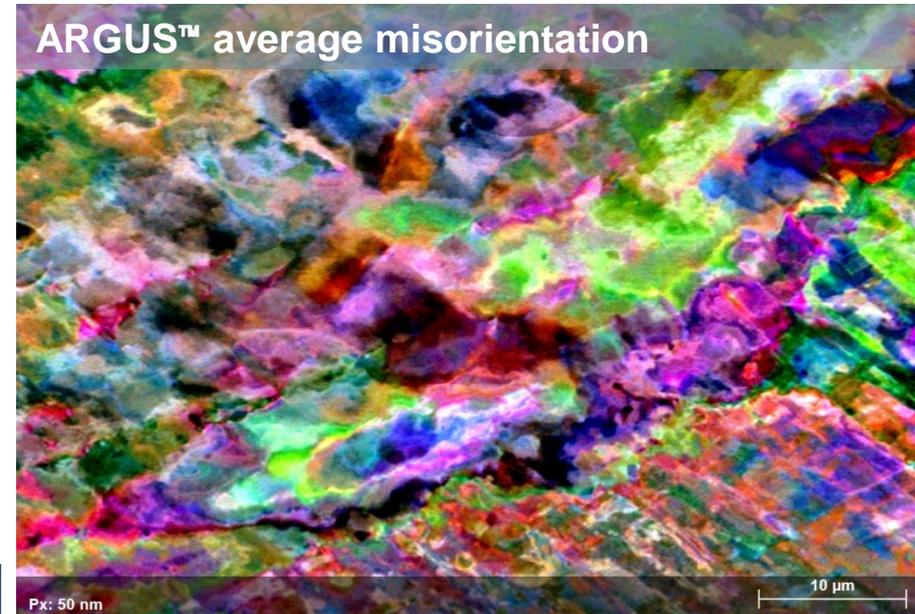

- Ion milling removes topography and mechanical preparation artifacts from the sample. An iterative workflow between the SEM Mill and the ARGUS™ imaging system results in optimal sample preparation.
- ARGUS™ orientation contrast imaging displays the quality of the sample preparation; orientation (IPF) and misorientation (KAM) maps reveal the microstructure of the material.

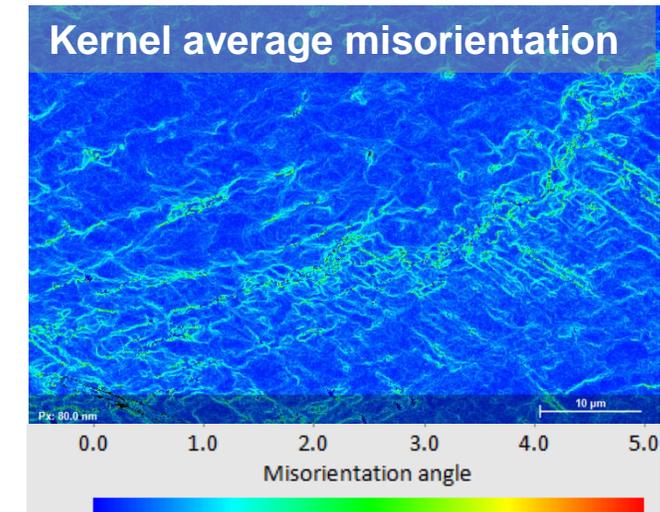



ARGUS™ imaging system

FSE imaging system


Color-coded orientation contrast



- SE images have very low contrast; not easy to find region of interest

FSE imaging system


Color-coded orientation contrast

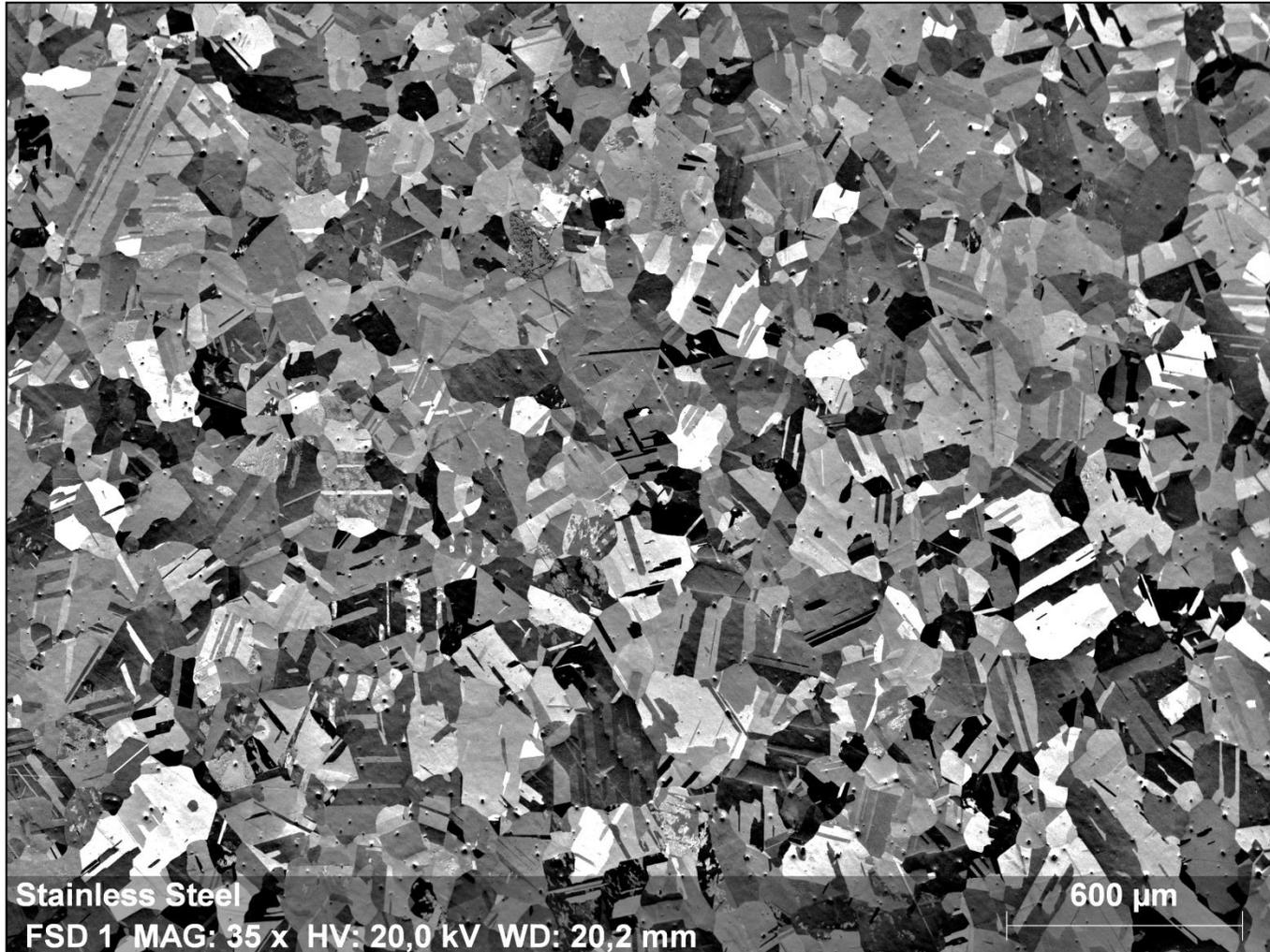

- SE images have very low contrast; not easy to find region of interest
- Common Fore Scattered Electron (FSE) detectors acquire images with better contrast

FSE imaging system


Color-coded orientation contrast

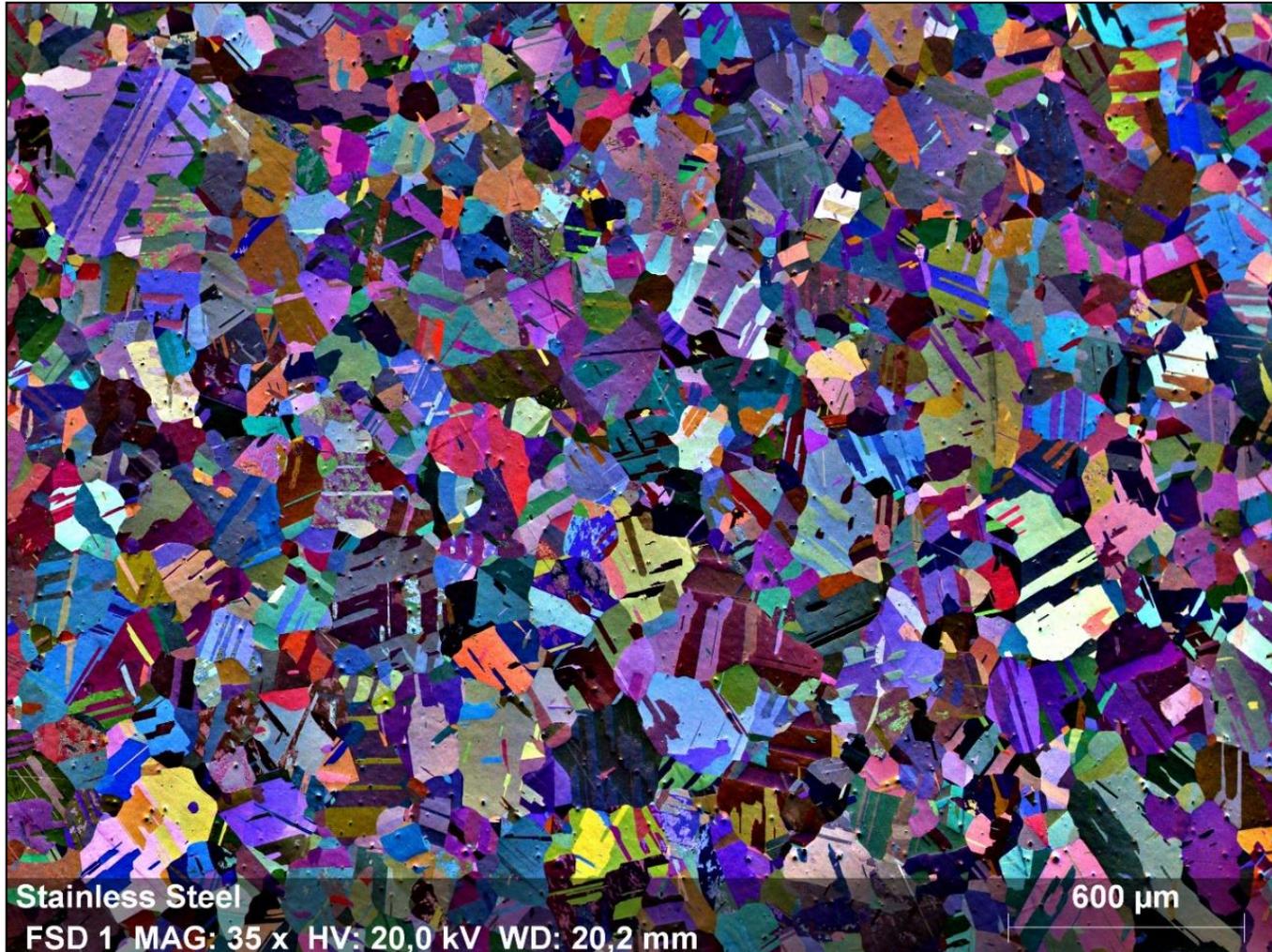

- SE images have very low contrast; not easy to find region of interest
- Common Fore Scattered Electron (FSE) detectors acquire images with better contrast
- However, ARGUS™ color-coded FSE images can reveal much more detail with **acquisition speed of 125 000 pps**

FSE imaging system with ARGUS™


How does it work?

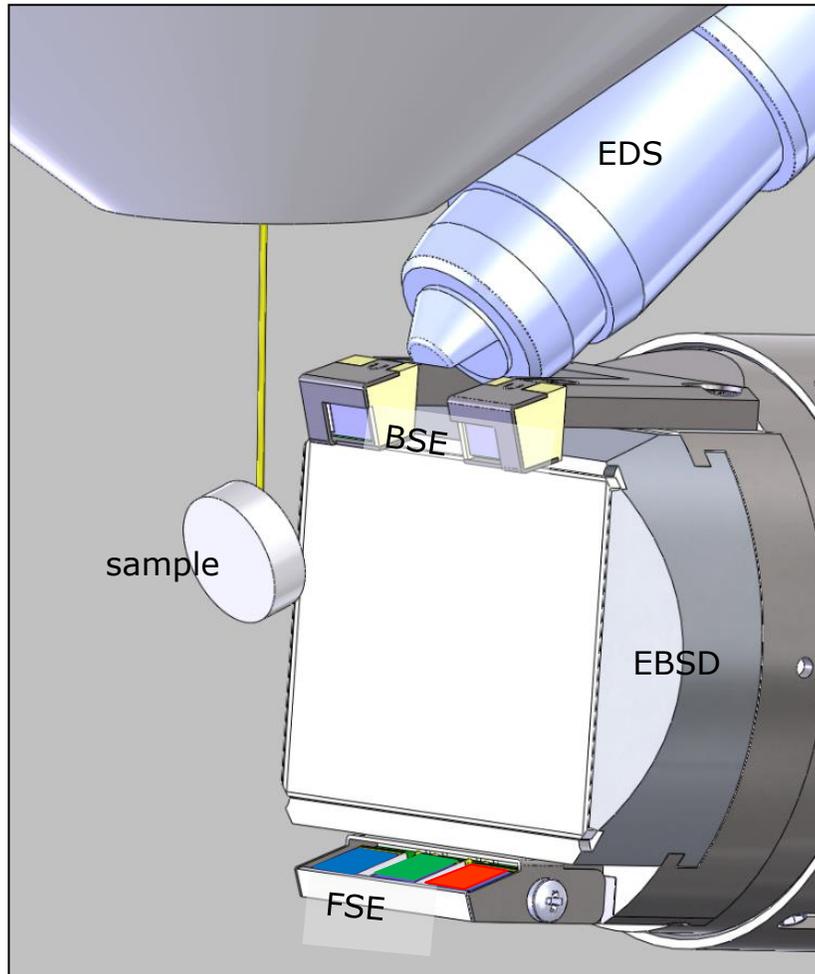

- Each diode captures a similar amount of noisy backscattered electrons and a different small part of the diffracted backscattered electrons, i.e., EBSD signal
- Signal (e^- counts) is transferred using a separate channel for each diode
- When scanning, for each pixel the system will obtain three numbers which will be transformed into three RGB levels
- When the three signals/RGB levels are mixed we obtain...


FSE imaging system with ARGUS™

How does it work?

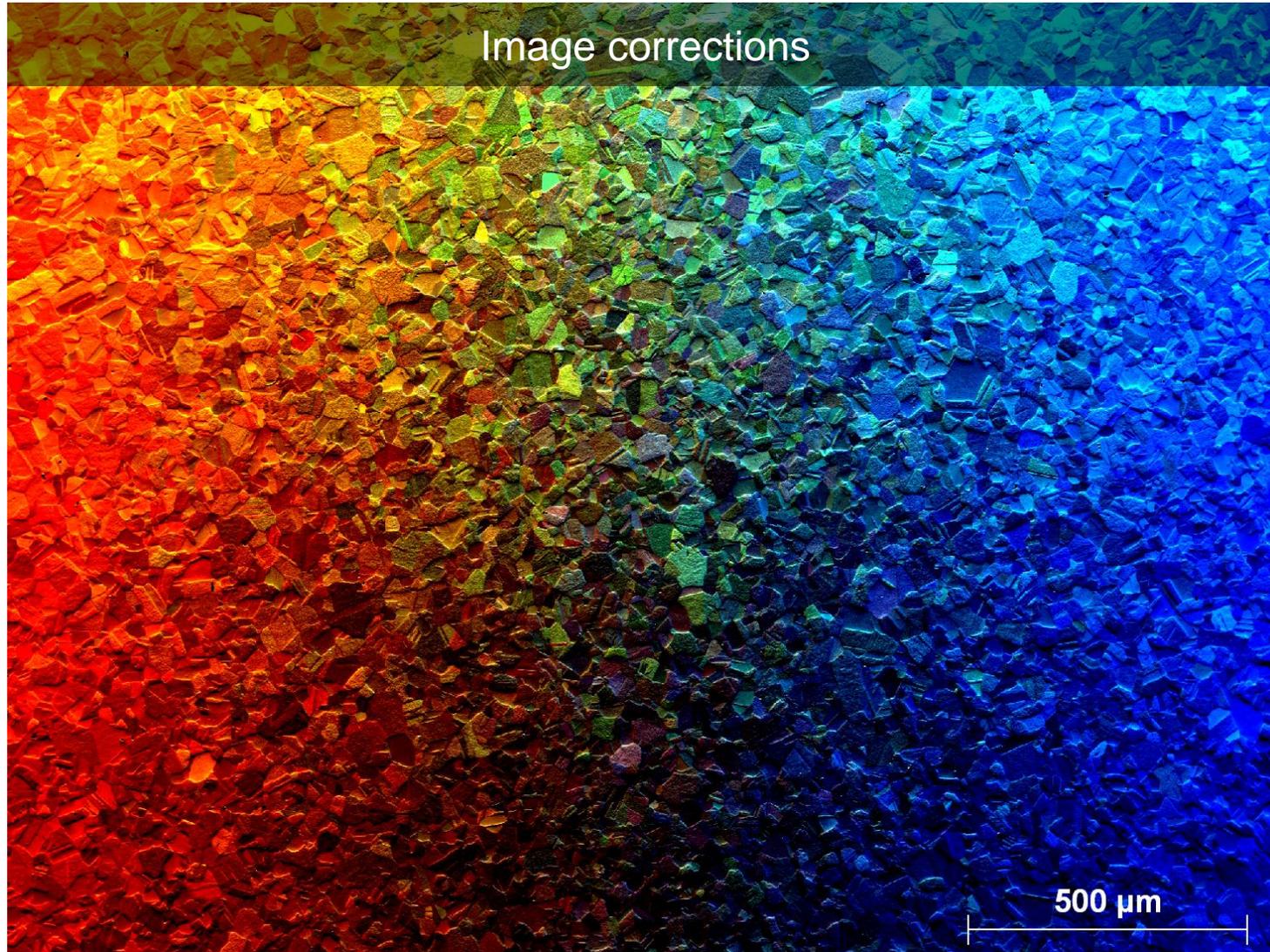

FSE imaging system with ARGUS™

How does it work?

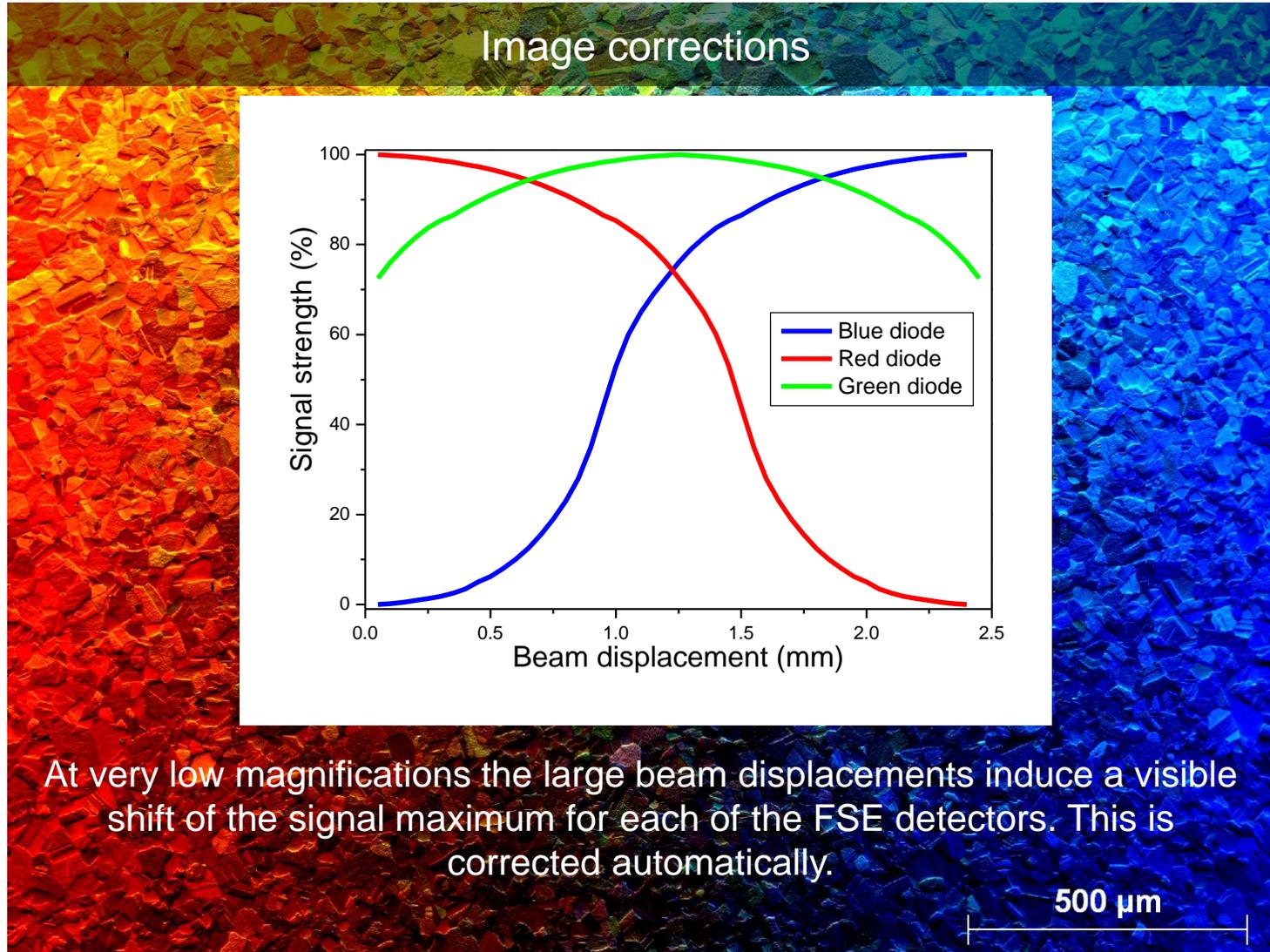


FSE imaging system with ARGUS™

How does it work?



FSE imaging system with ARGUS™

How does it work?

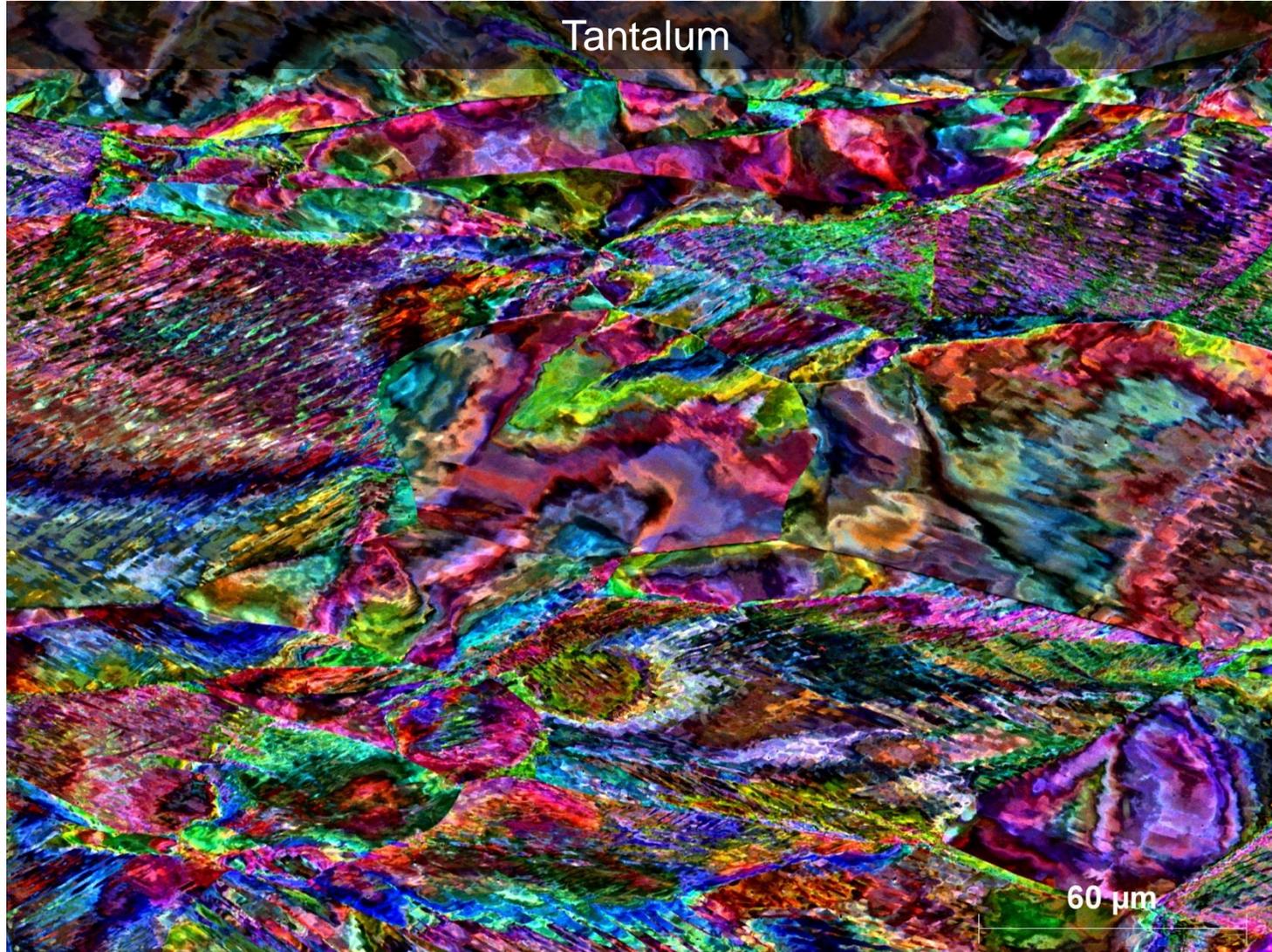



Broad Ion Beam Milling Model 1061 SEM Mill

Model 1061 SEM Mill

A versatile argon broad ion beam mill for SEM applications




- Produces ideal results for both **planar and cross-section** samples
- **Wide-ranging ion energies** allow either rapid milling or gentle polishing on a broad variety of sample materials
- **Artifact-free samples** are readily produced
- **Easy-to-use interface**
- **Fully automated**, including precise sample height detection, for high-throughput applications
- Liquid nitrogen **stage cooling**
- **Vacuum / inert gas transfer** to SEM capability

Model 1061 SEM Mill


Milling geometries



Planar milling

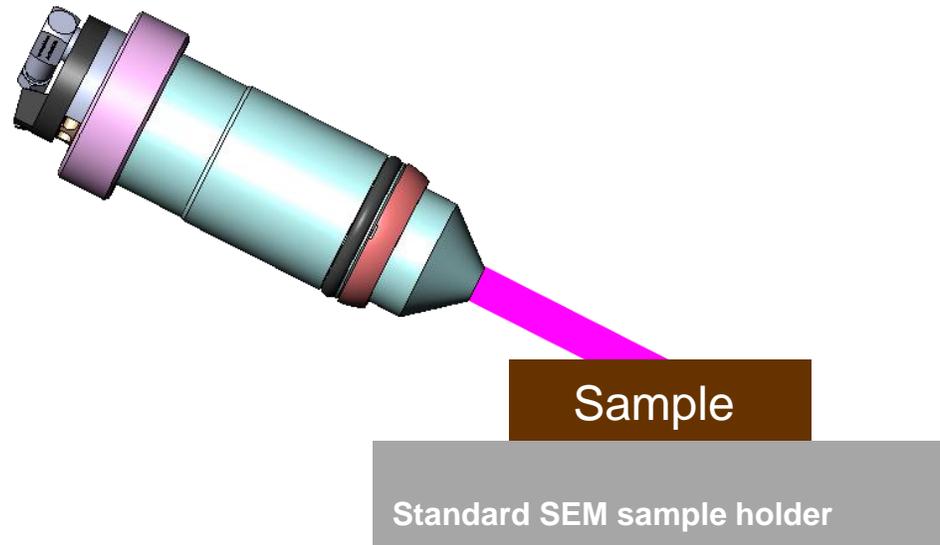

Cross-section milling

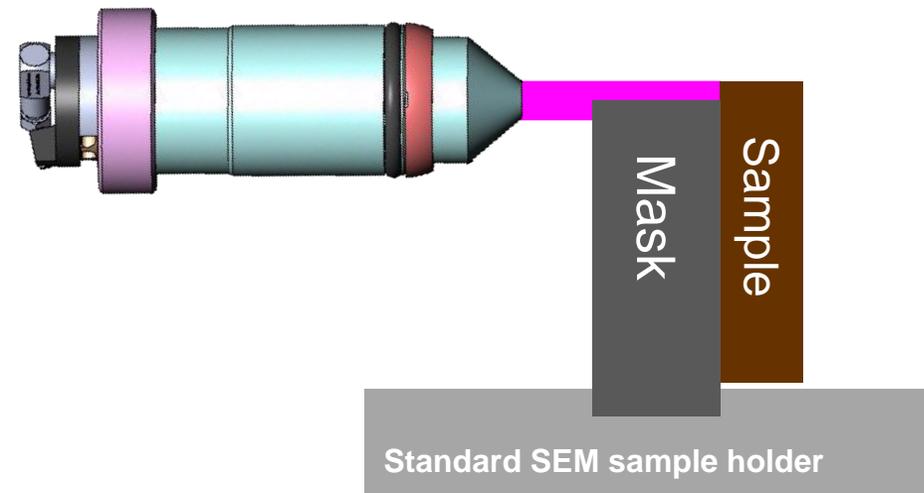


Ion energies range from 0.1 to 10 keV
Variable milling angle range: 0 to 10°
Stage movement: Continuous rotation or rocking

Model 1061 SEM Mill


Milling geometries



Planar milling

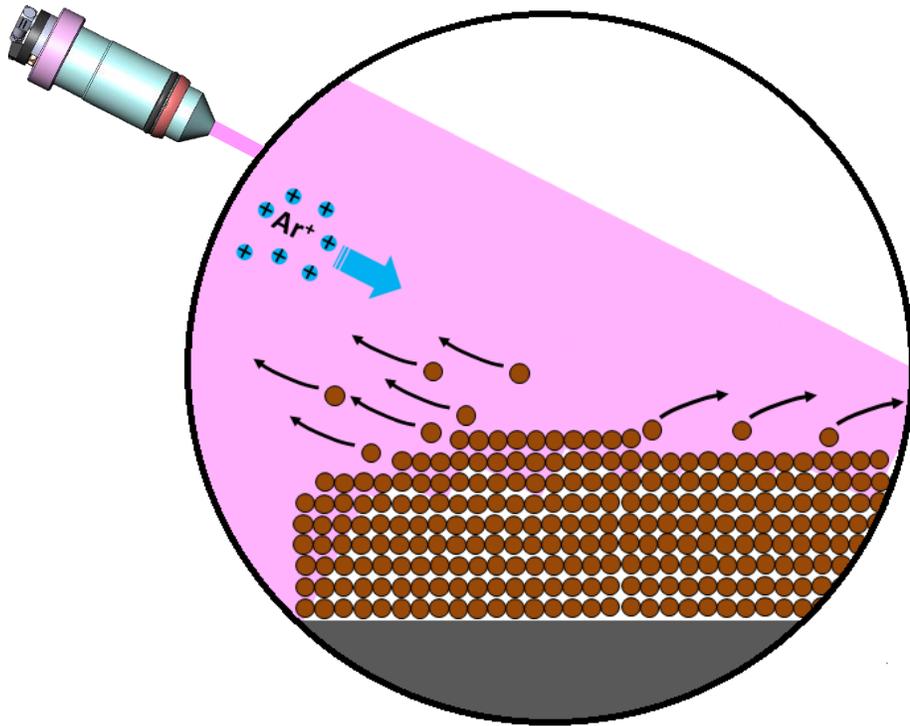


Cross-section milling

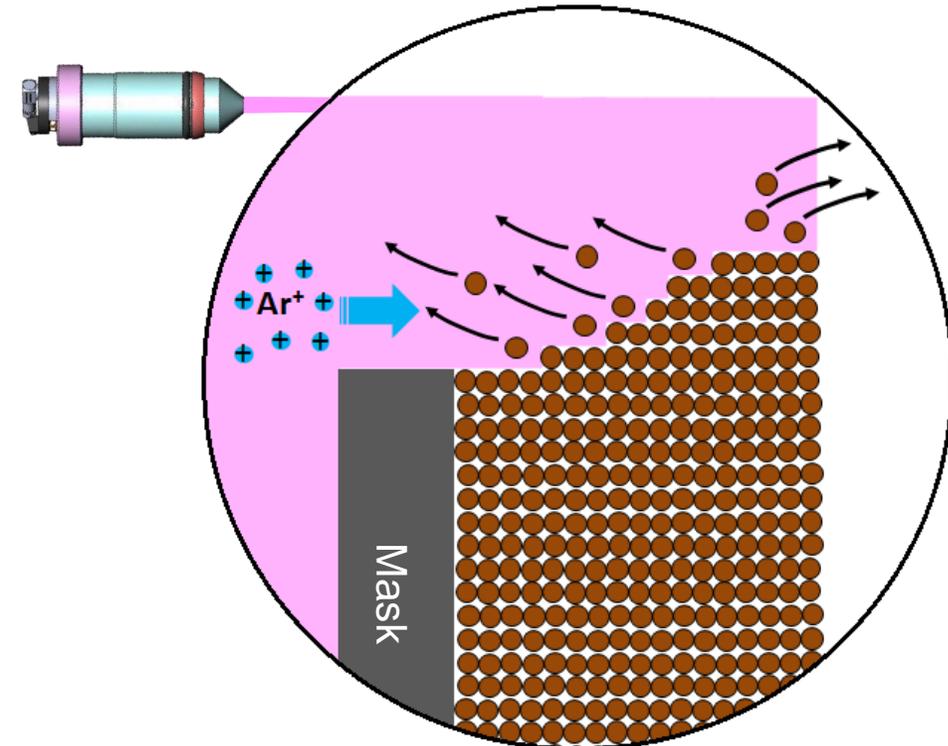


Model 1061 SEM Mill

Planar sample milling

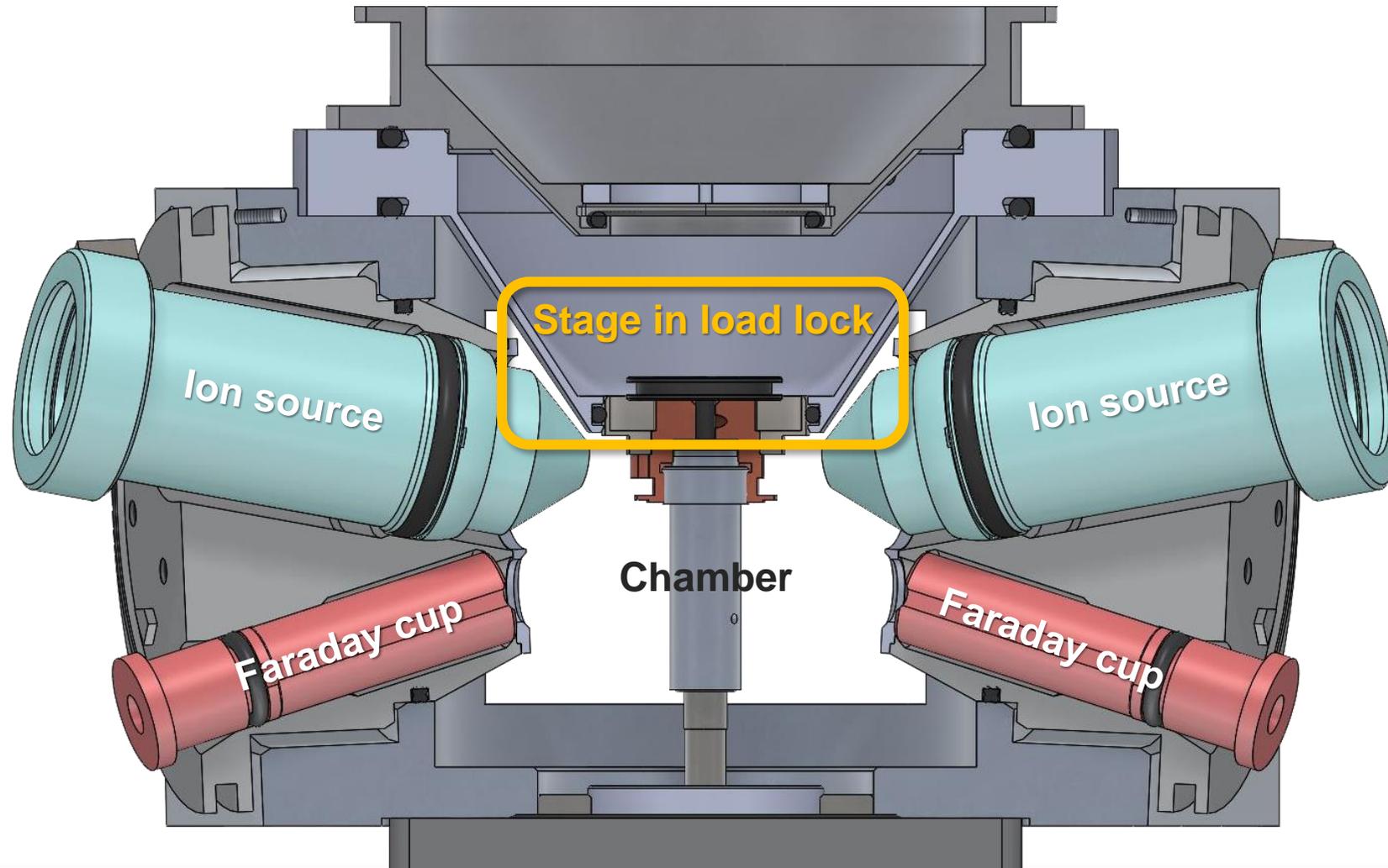


Model 1061 SEM Mill

Planar sample milling

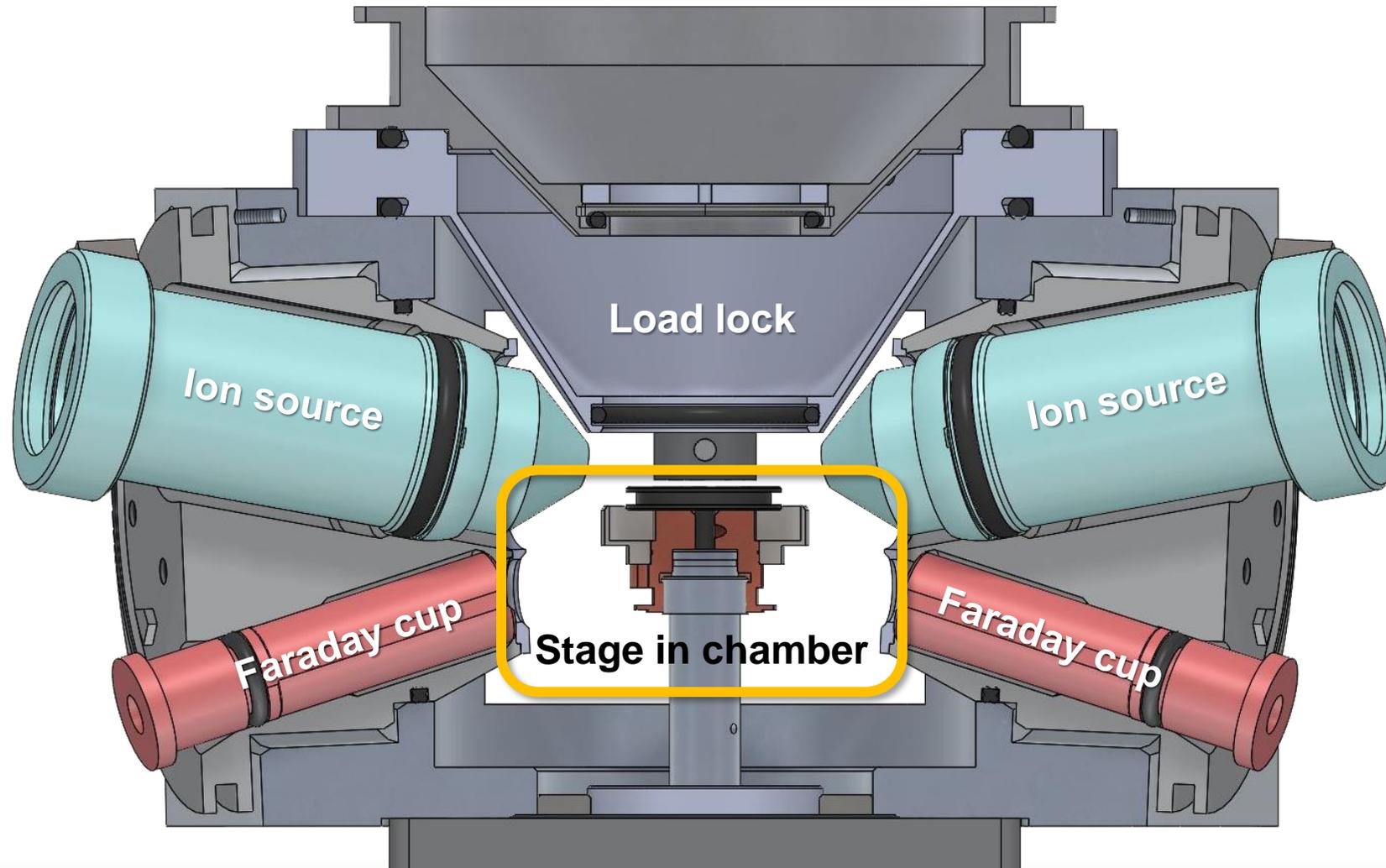


Model 1061 SEM Mill

Sample height detection

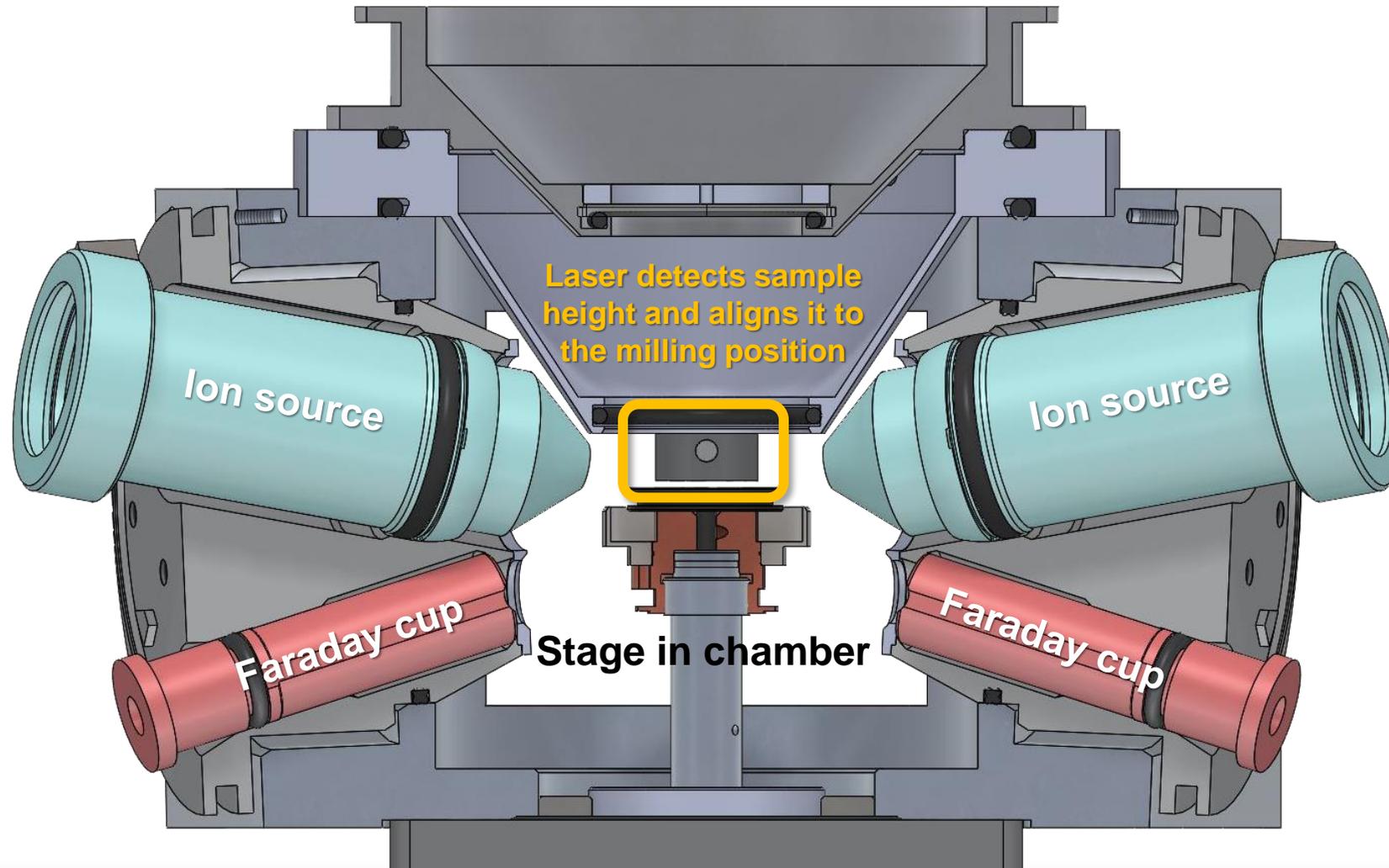

Model 1061 SEM Mill

Cross-section sample milling

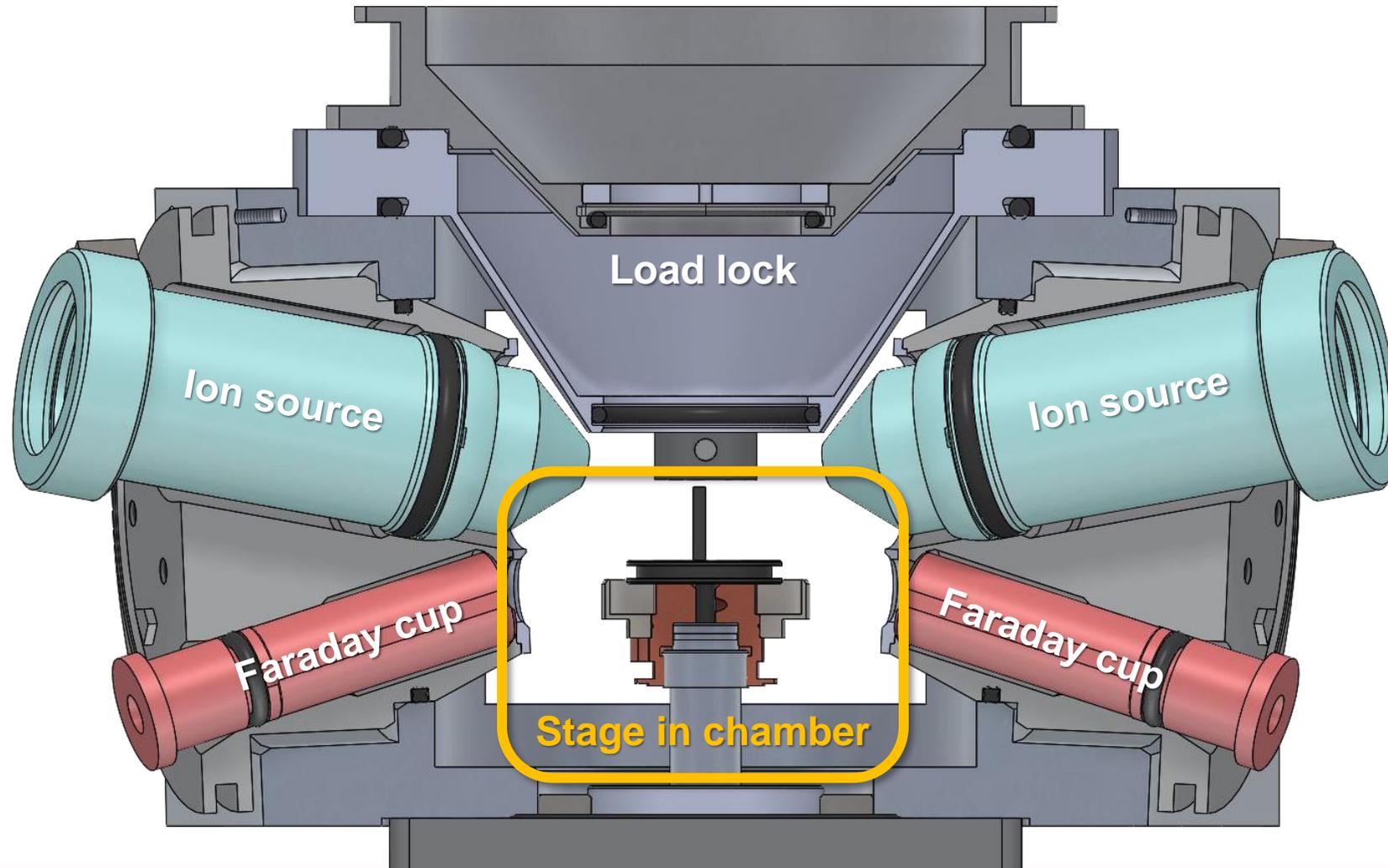

Model 1061 SEM Mill

TrueFocus ion sources



5 keV; 44 μ A; 0% focus



5 keV; 44 μ A; 50% focus




5 keV; 44 μ A; 100% focus




- Adjustable energy range from 100 eV to 10 keV
- Ion beam spot size ranges from 300 μ m to 5 mm
- Consistent beam current
- Motorized ion sources enable touch screen control of beam angle adjustments

Broad ion milling

Sample preparation

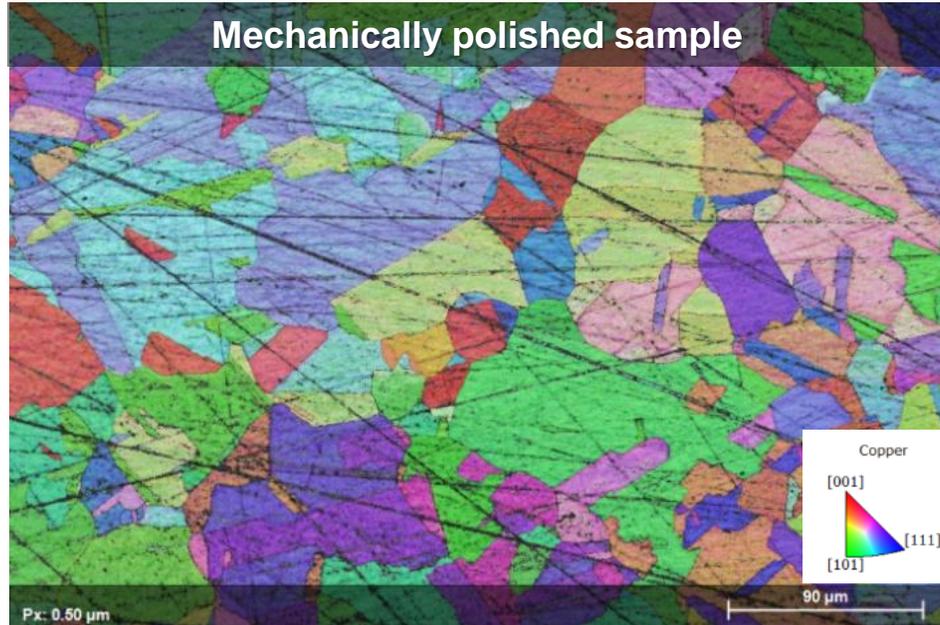

The specimen surface must be deformation-free to conduct EBSD measurements.

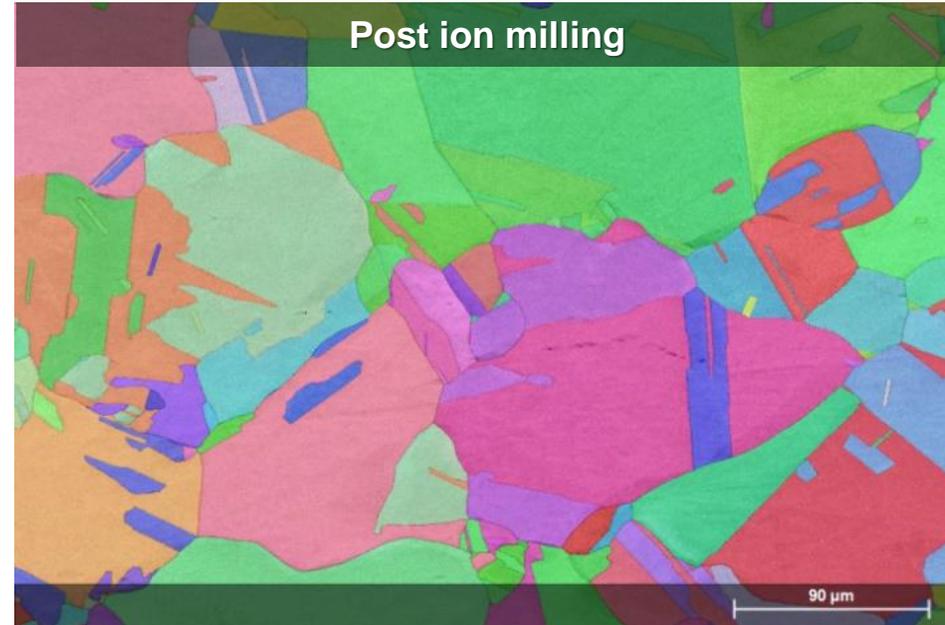

Scratches and damages caused by mechanical polishing are removed by the ion milling process.

As a result, the material's microstructure is revealed, and statistical data can be acquired.




Austenitic steels

Austenitic 300 series

Orientation contrast imaging with ARGUS™



Mechanically polished using colloidal silica


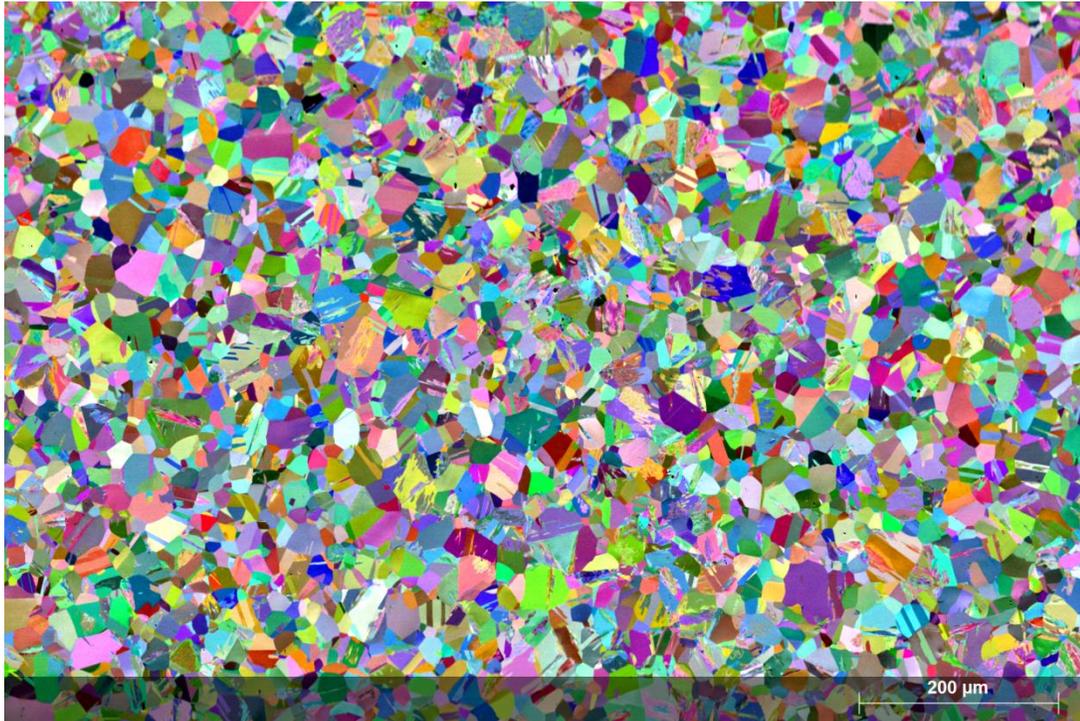


Austenitic 300 series

Orientation contrast imaging with ARGUS™






Mechanically polished using colloidal silica

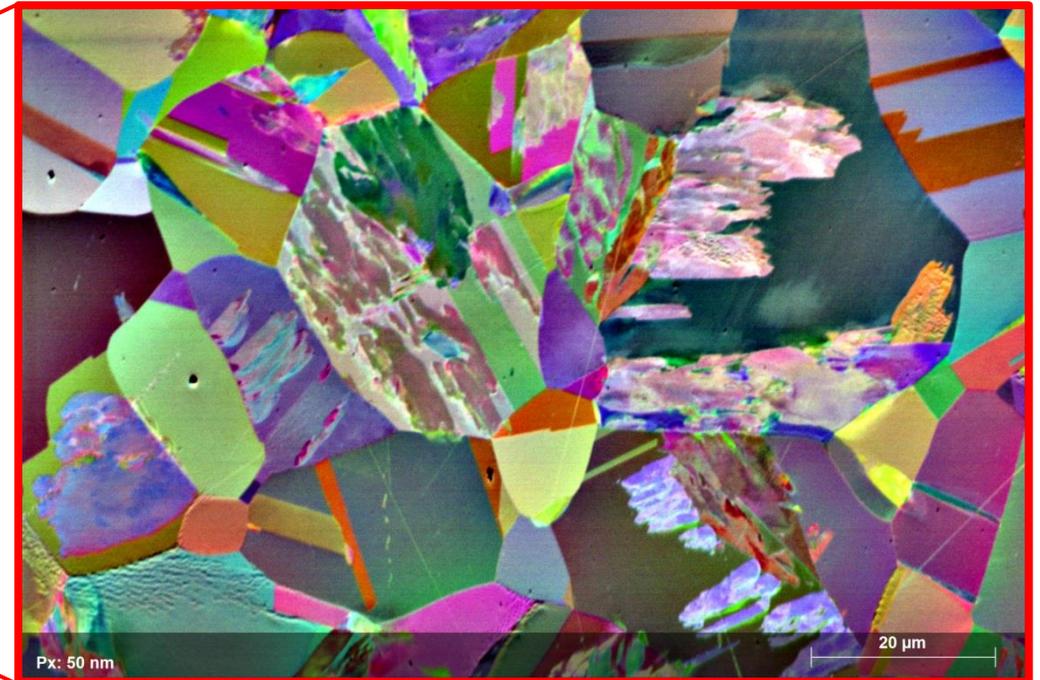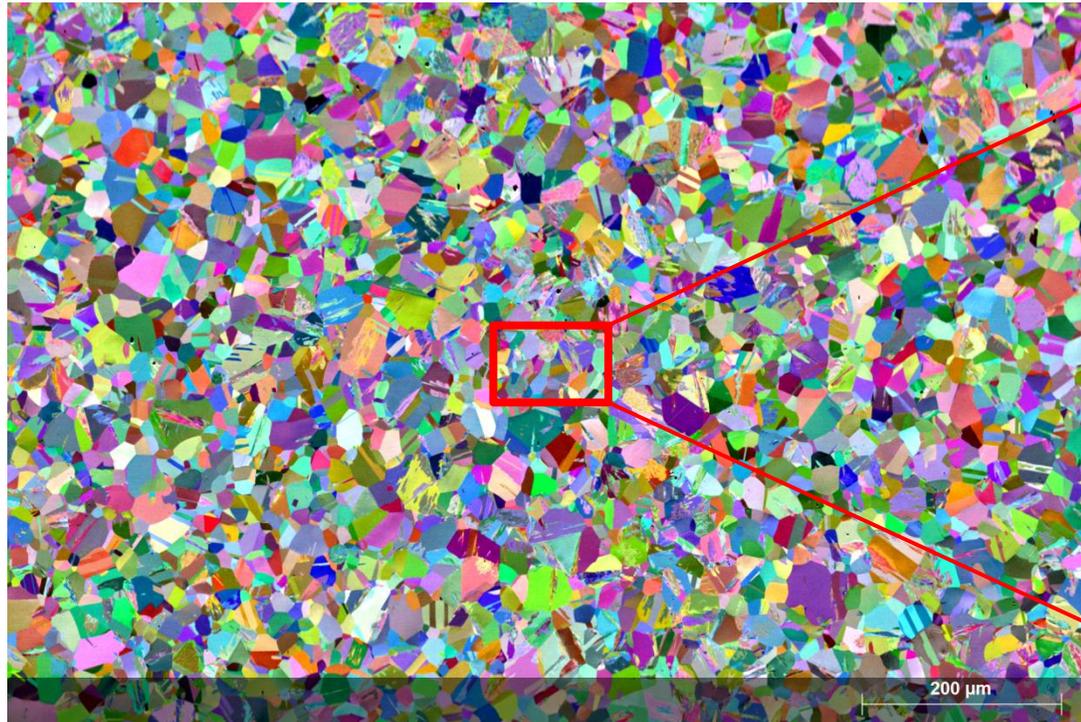

Austenitic 300 series

Orientation contrast imaging with ARGUS™

Mechanically polished using colloidal silica

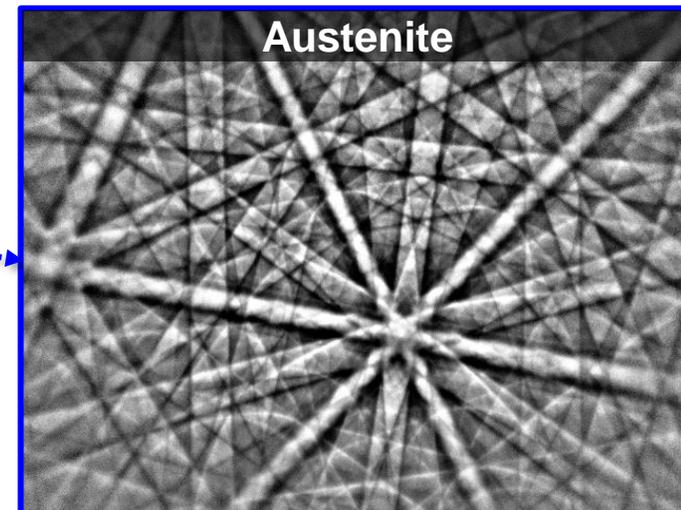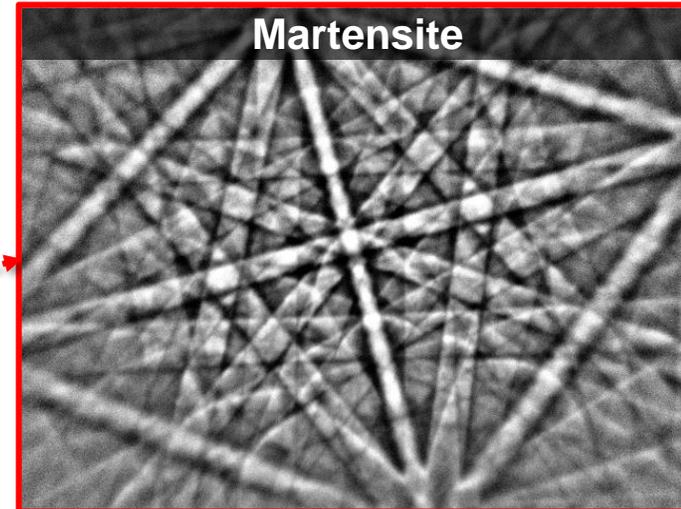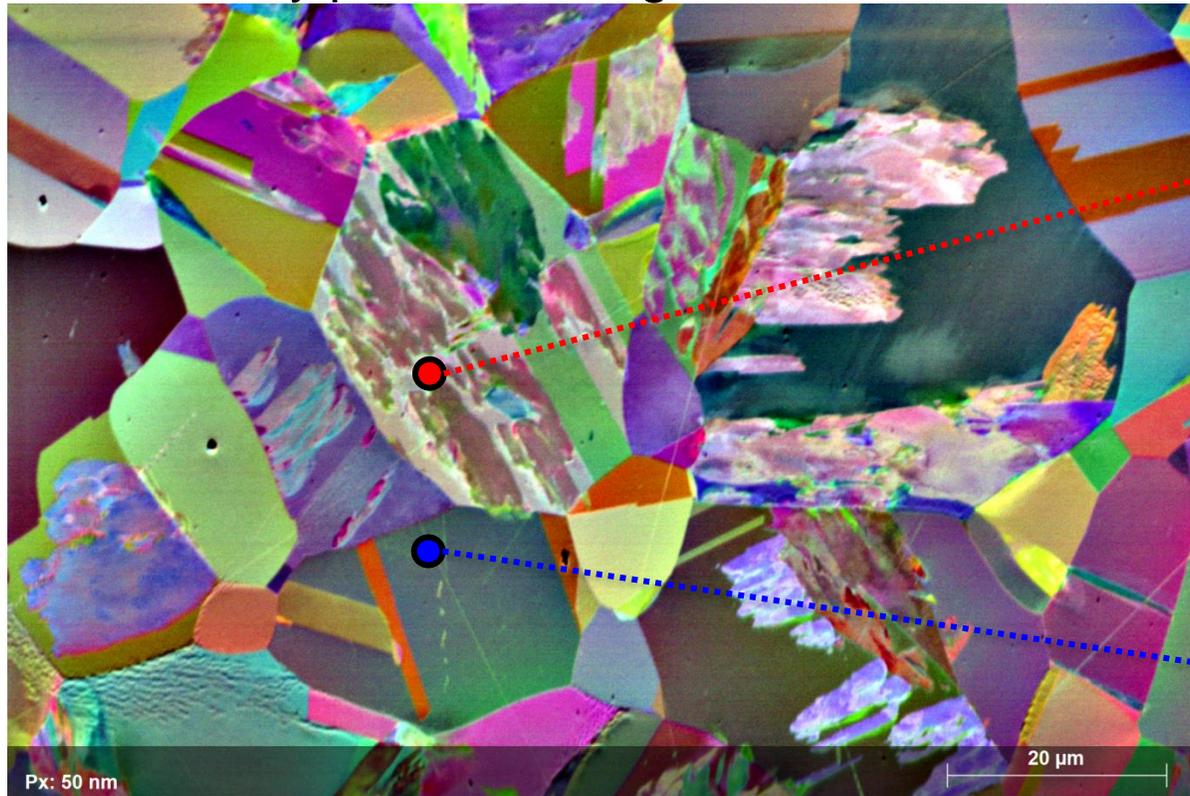


Dynamic strain-induced phase transformation of austenite to martensite during mechanical polishing

Austenitic 300 series

EBSD phase distribution map



Mechanically polished using colloidal silica

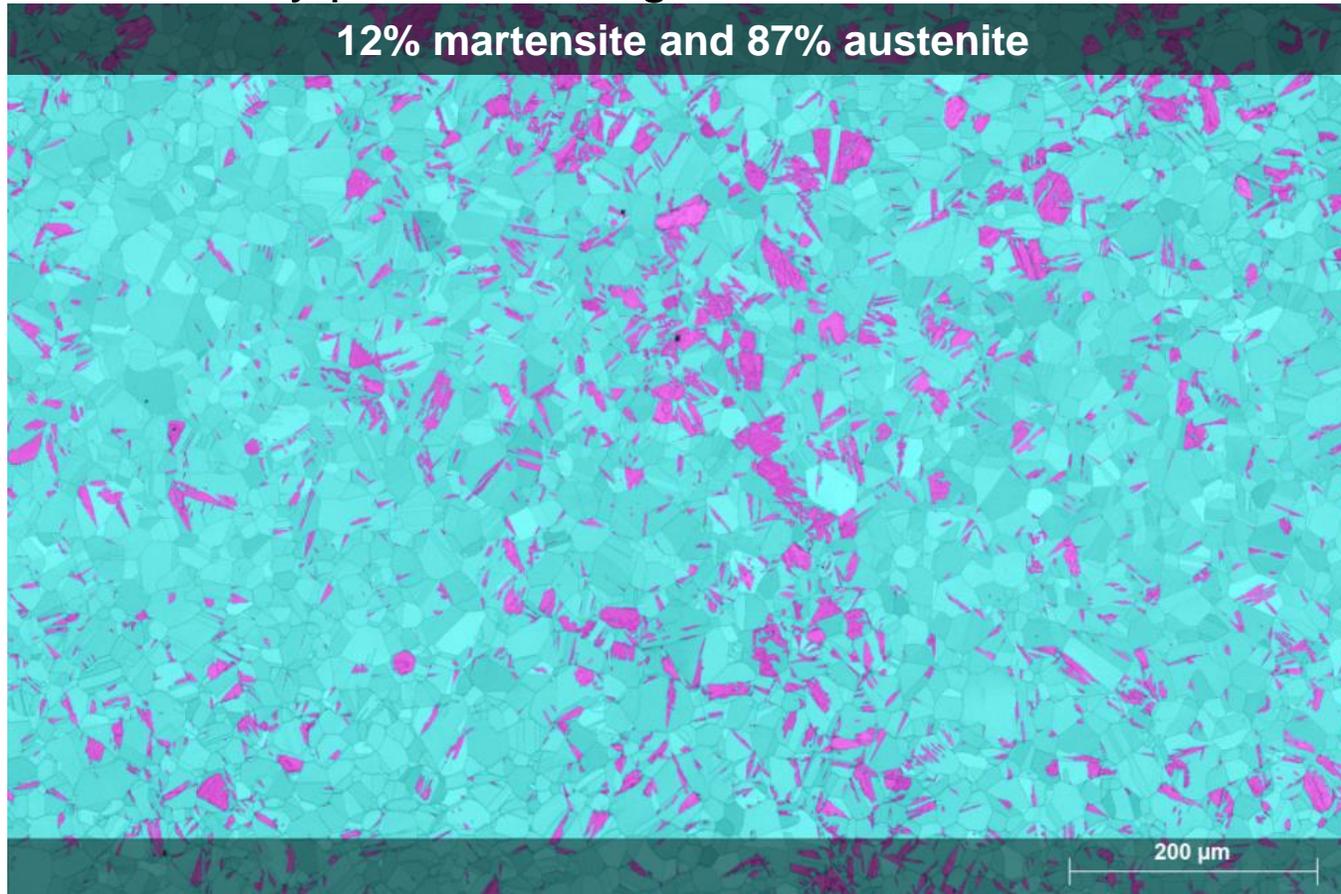

Phase map

Step size: 0.5 μm

Pattern resolution: 80x60 pixels

Acquisition speed: 950 Hz

Indexing rate: 99%

	Martensite	12%
	Austenite	87%


Dynamic strain-induced phase transformation of austenite to martensite during mechanical polishing

Austenitic 300 series

EBSD phase distribution map



Ion milled at 5 and 2 keV energy; 2° angle

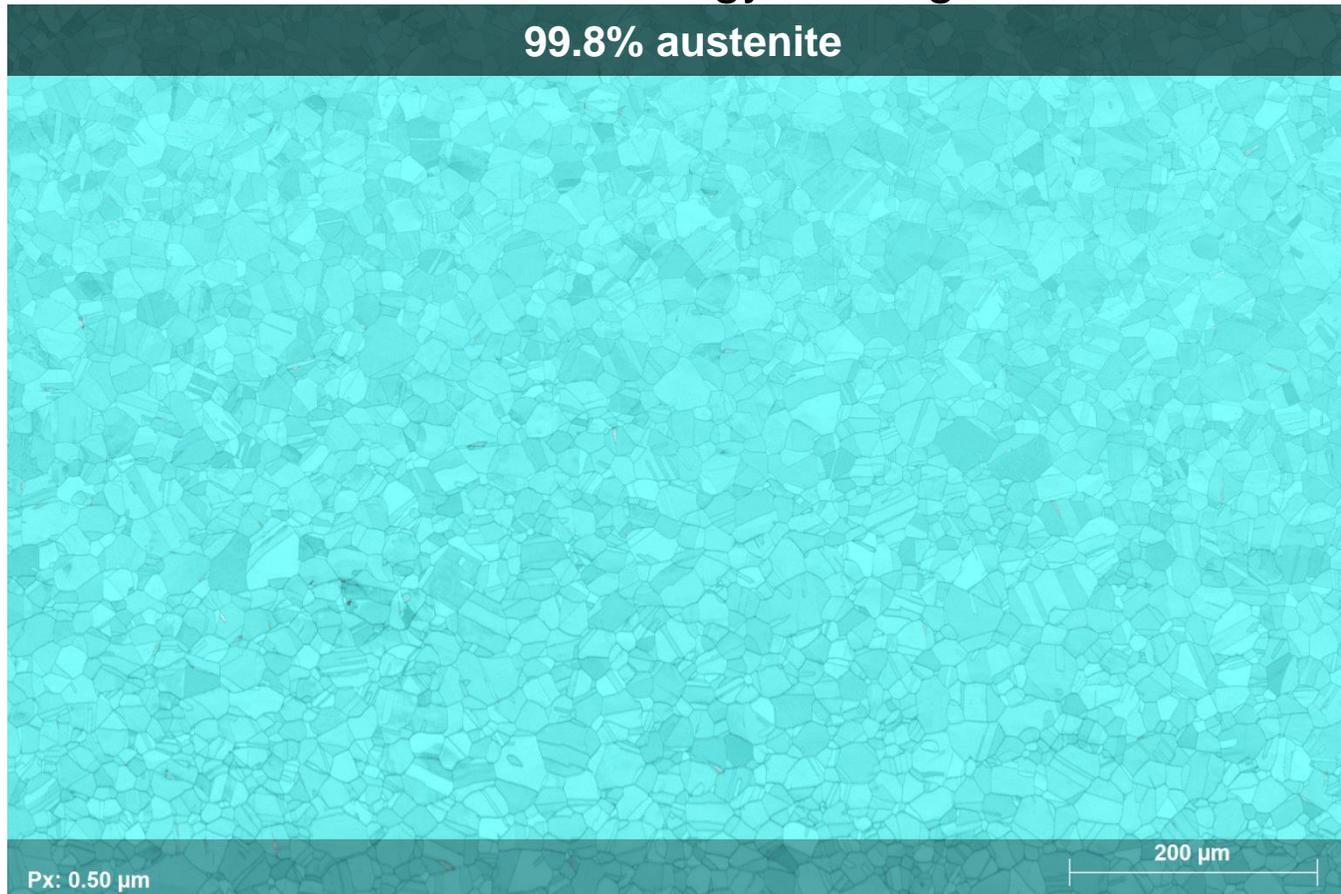

Phase map

Step size: 0.5 μm

Pattern resolution: 80x60 pixels

Acquisition speed: 950 Hz

Indexing rate: 99.8%

	Martensite	0%
	Austenite	99.8%


No dynamic strain-induced phase transformation

Austenitic 300 series


Texture



Mechanically polished: colloidal silica

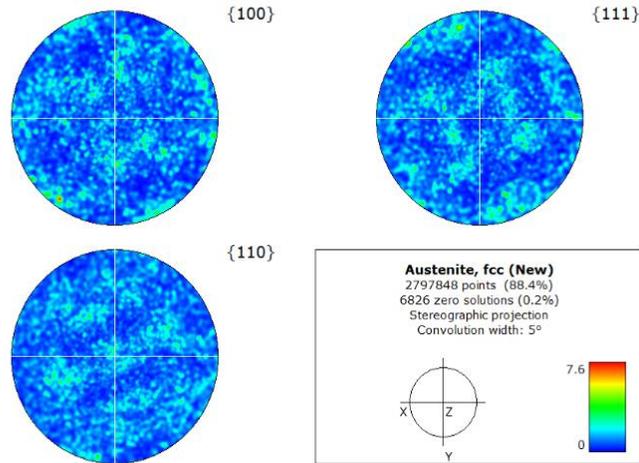

Ion milled: 5 and 2 keV energy; 2° angle

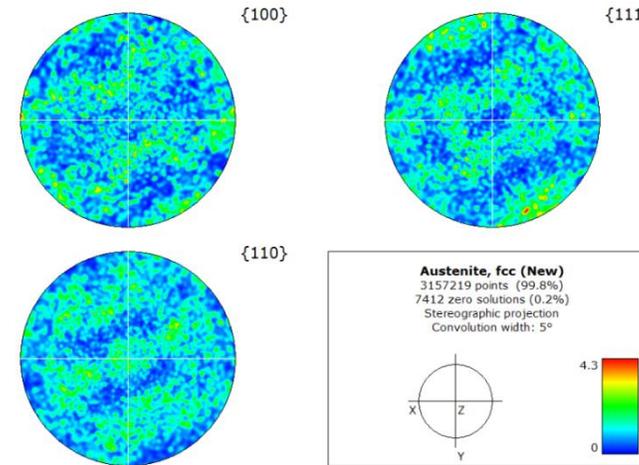


Austenitic 300 series



Texture



Mechanically polished: colloidal silica

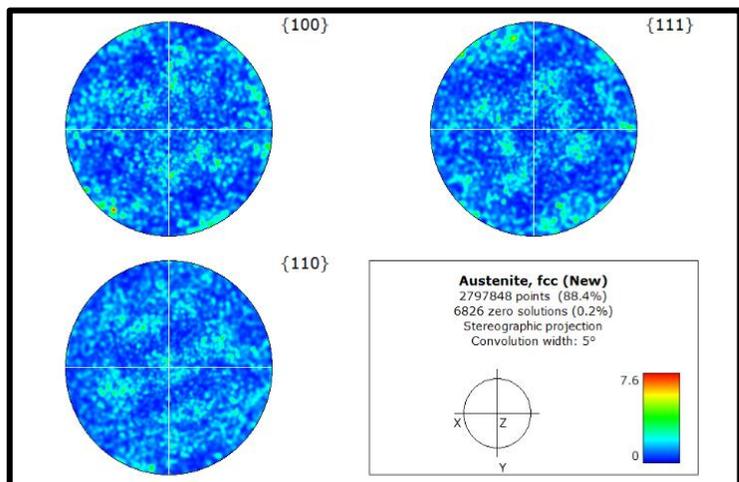


Ion milled: 5 and 2 keV energy; 2° angle

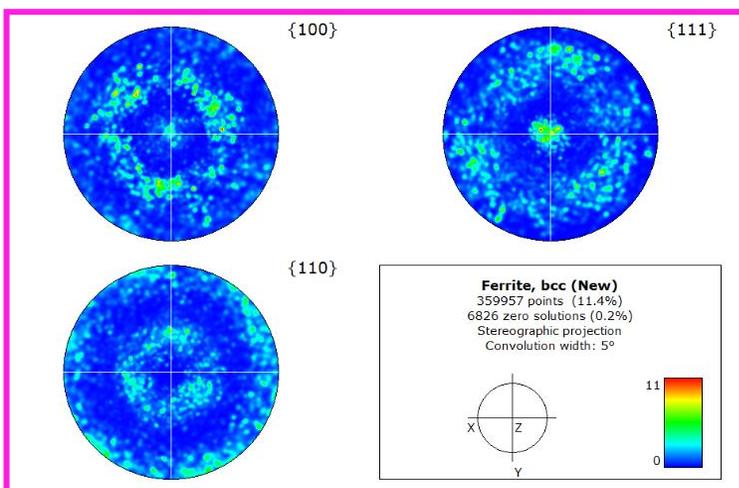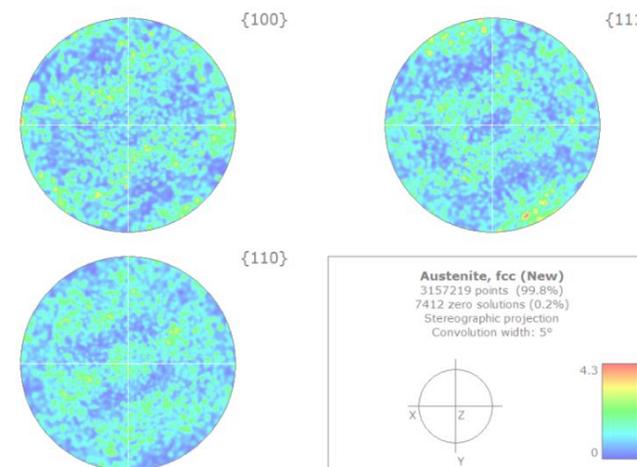



Austenitic 300 series

Texture

Mechanically polished: colloidal silica

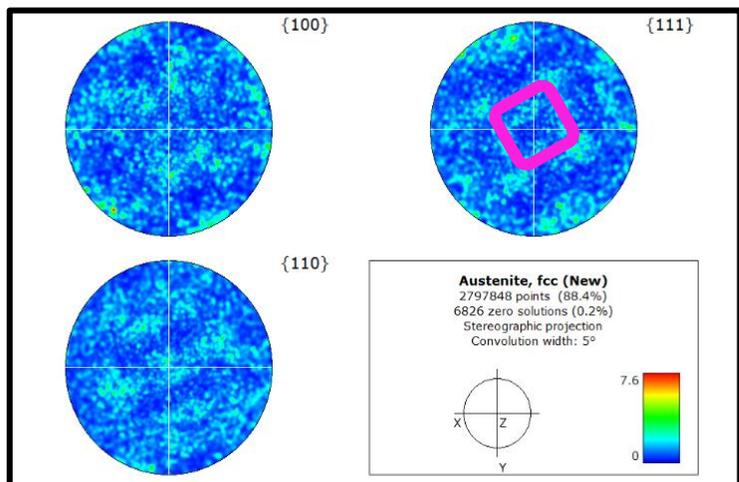

Ion milled: 5 and 2 keV energy; 2° angle

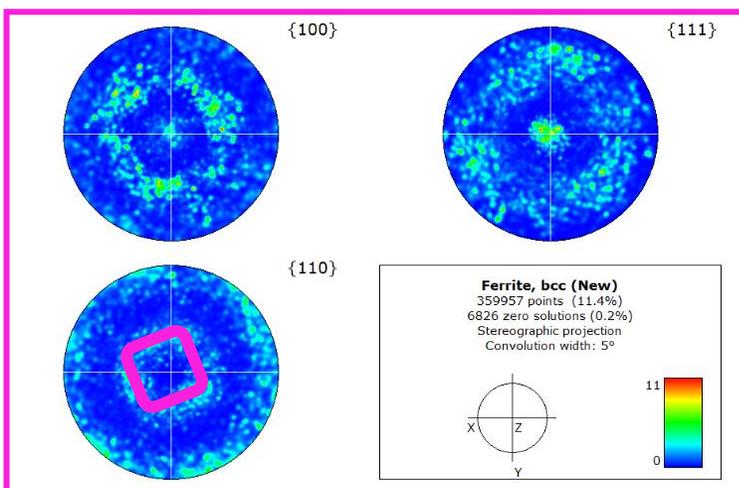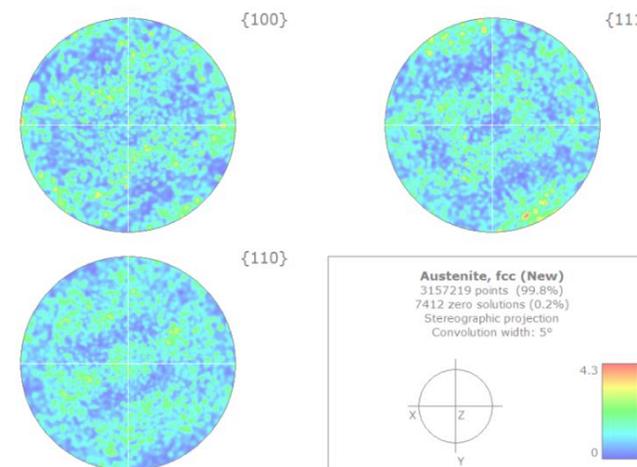

Kurdjumov–Sachs orientation relationship
 $\{111\}_\gamma \parallel \{110\}_{\alpha'}$ and $\langle 1-10 \rangle_\gamma \parallel \langle 11-1 \rangle_{\alpha'}$


Austenitic 300 series

EBSD phase distribution map



Mechanically polished: colloidal silica

12% martensite and 87% austenite

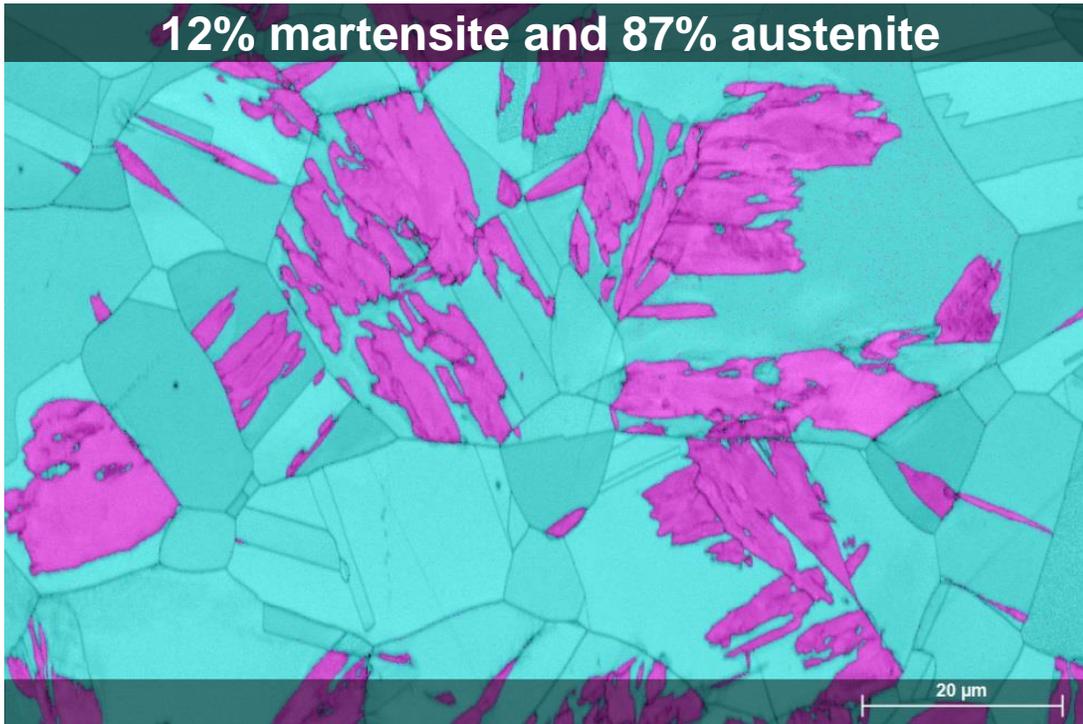

Ion milled: 5 and 2 keV energy; 2° angle

100% austenite



Step size: 100 nm
Pattern resolution: 160x120 pixels
Indexing rate: 99%




Austenitic 300 series

EBSD IPF map




Mechanically polished: colloidal silica

12% martensite and 87% austenite

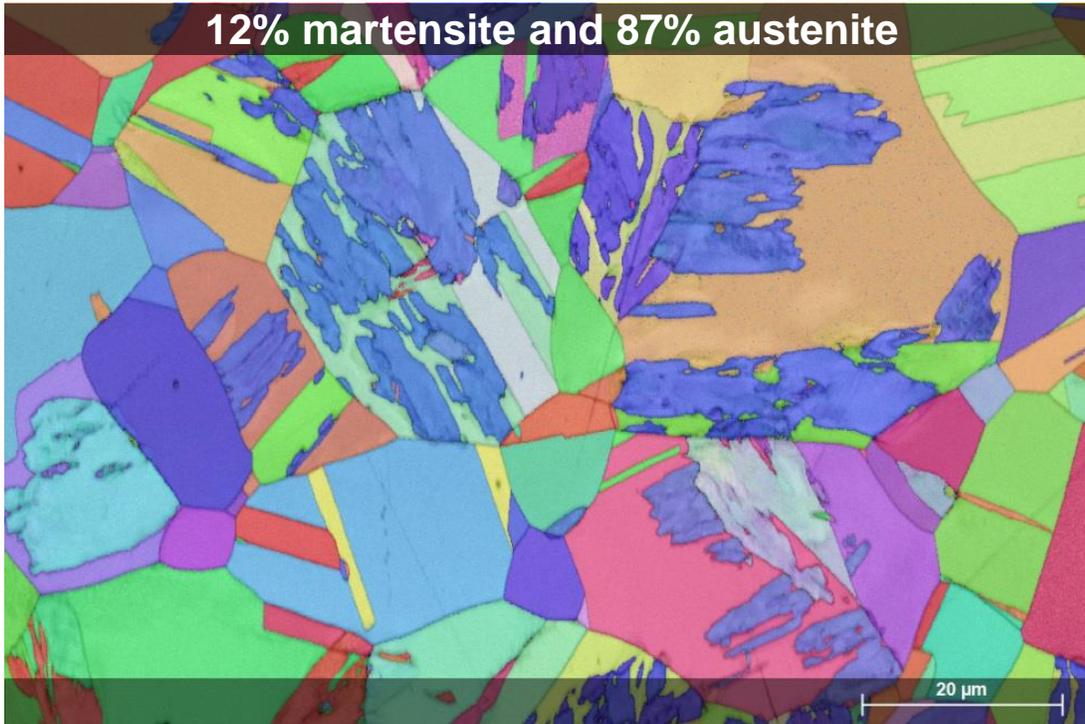


Ion milled: 5 and 2 keV energy; 2° angle

100% austenite

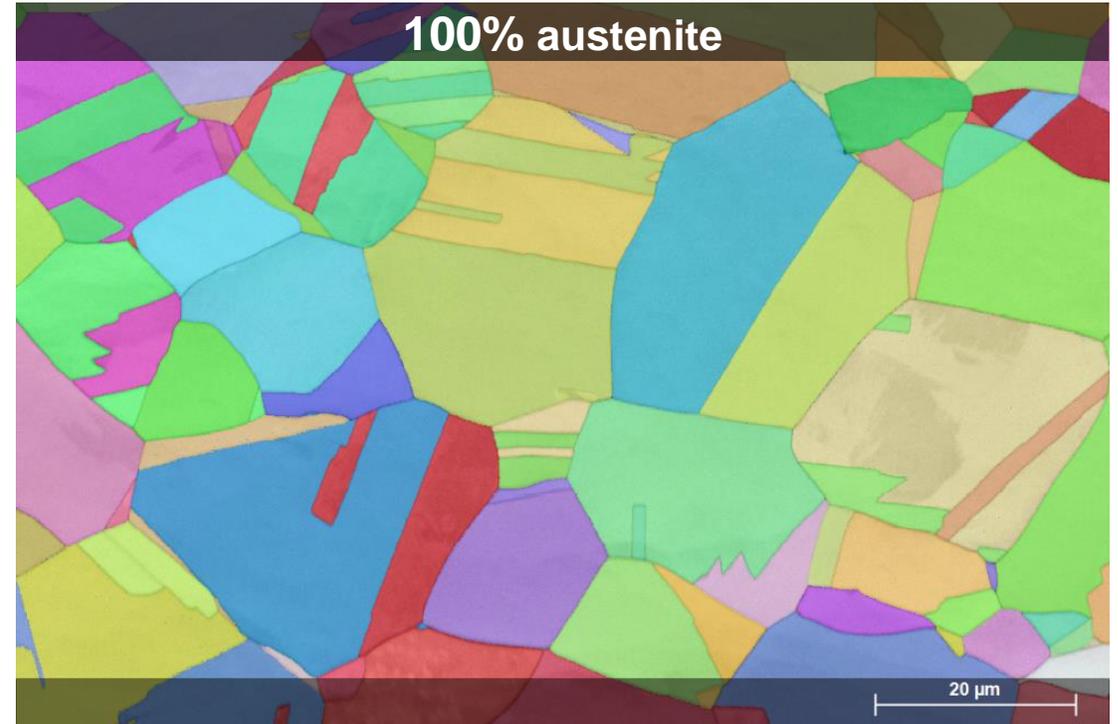

Step size: 100 nm
Pattern resolution: 160x120 pixels
Indexing rate: 99%

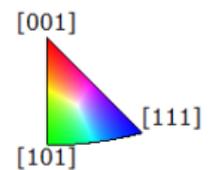


Austenitic 300 series

Kernel average misorientation map



Mechanically polished: colloidal silica


12% martensite and 87% austenite

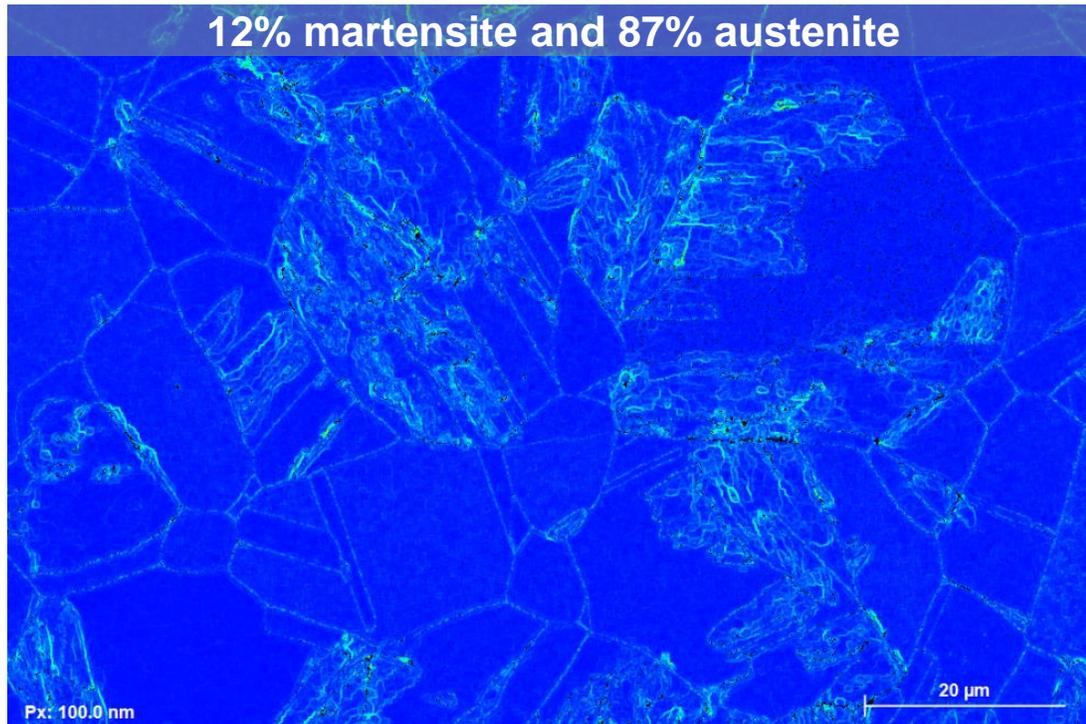

Step size: 100 nm
Pattern resolution: 160x120 pixels
Indexing rate: 99%

Ion milled: 5 and 2 keV energy; 2° angle

100% austenite

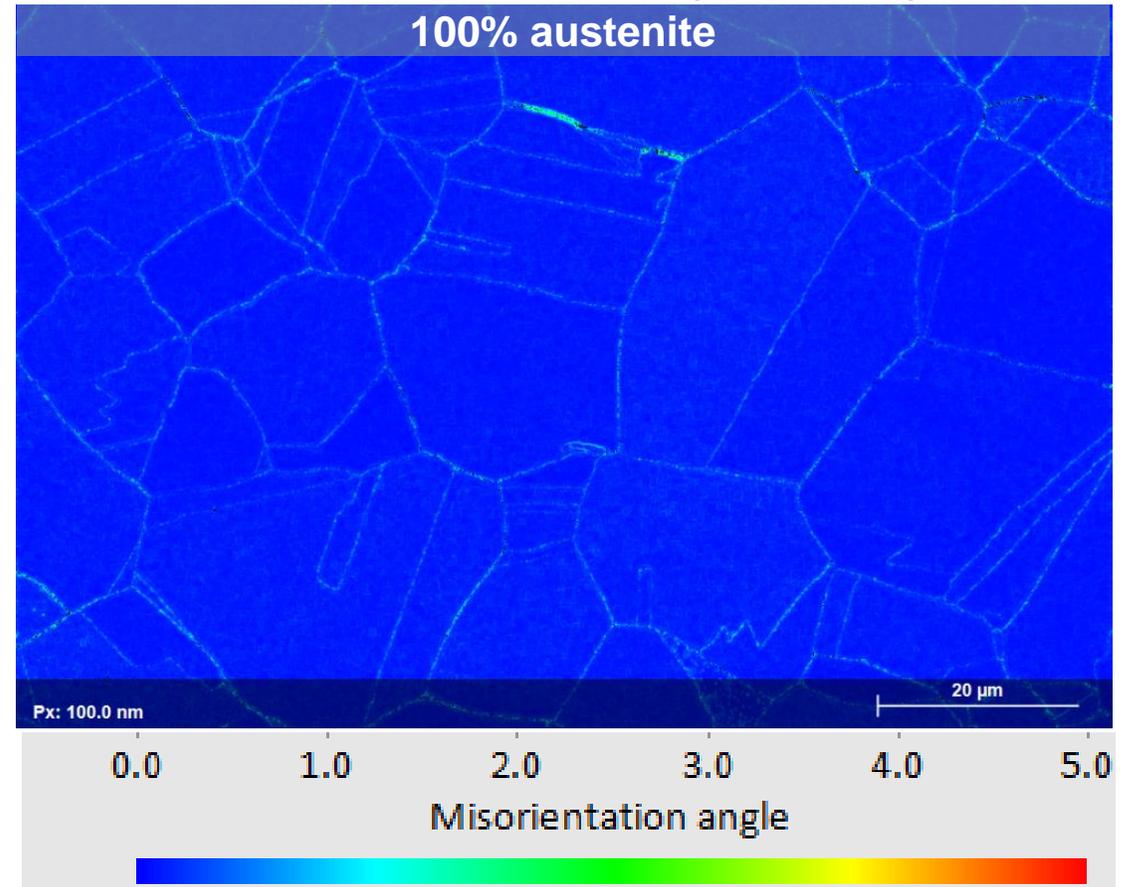




Zirconium alloy

Zirconium 702 alloy

Orientation contrast imaging with ARGUS™



Mechanically polished: colloidal silica

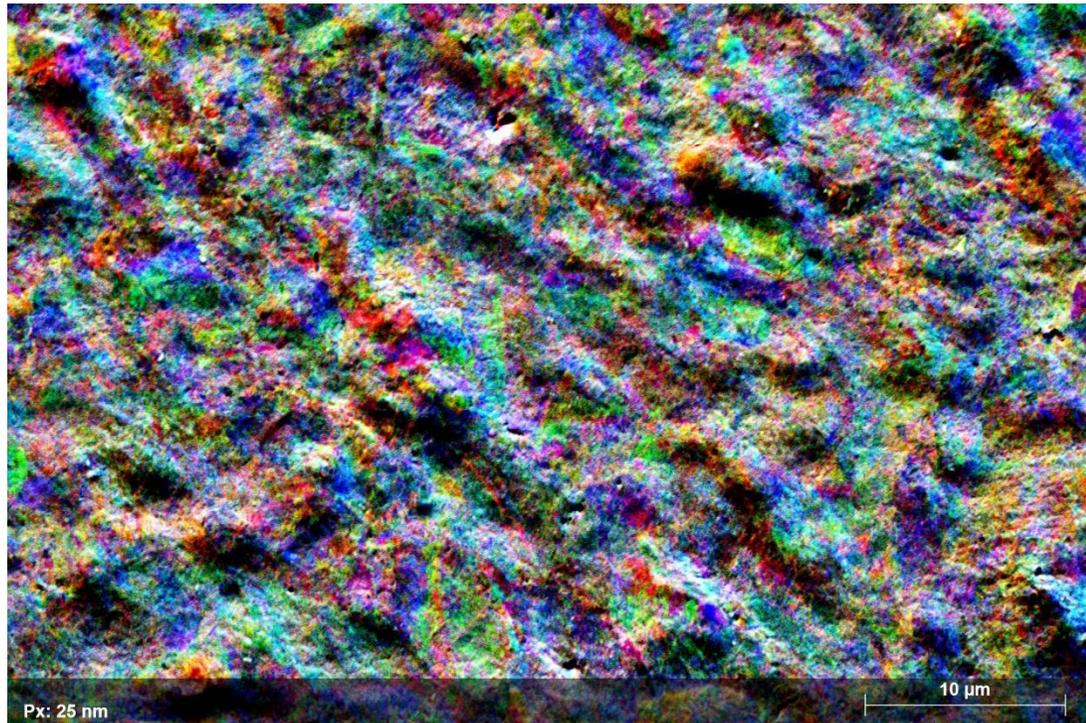

The microstructure is not visible

Ion milled: 8 and 2 keV energy; 3° angle




The fine microstructure is revealed
(inter- and intragranular)

Zirconium 702 alloy

EBSD patterns




Mechanically polished: colloidal silica

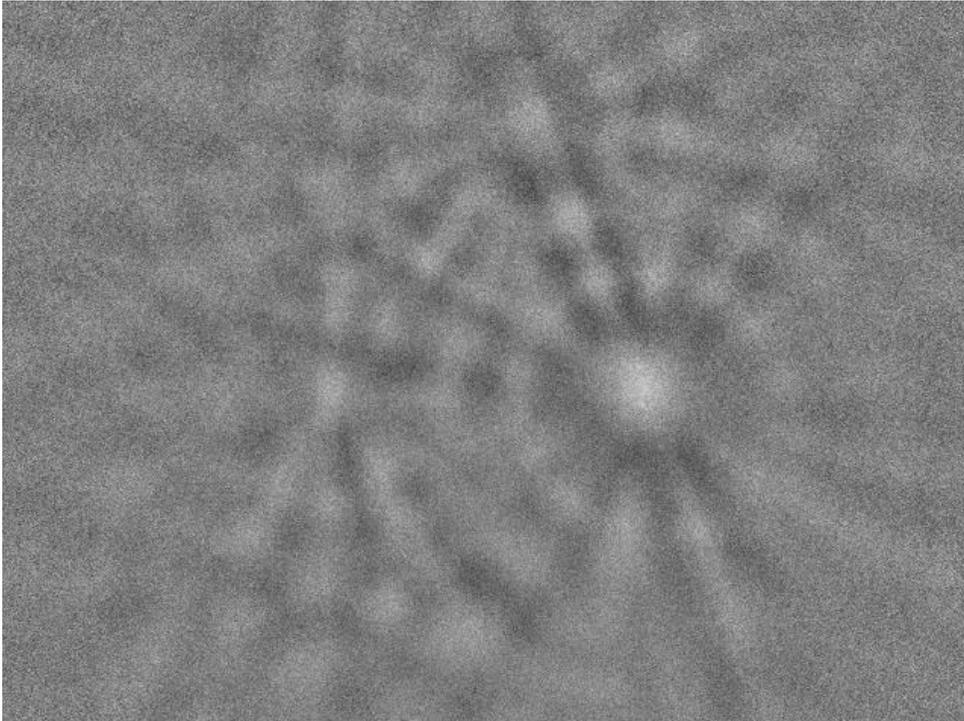

Very few patterns

Ion milled: 8 and 2 keV energy; 3° angle

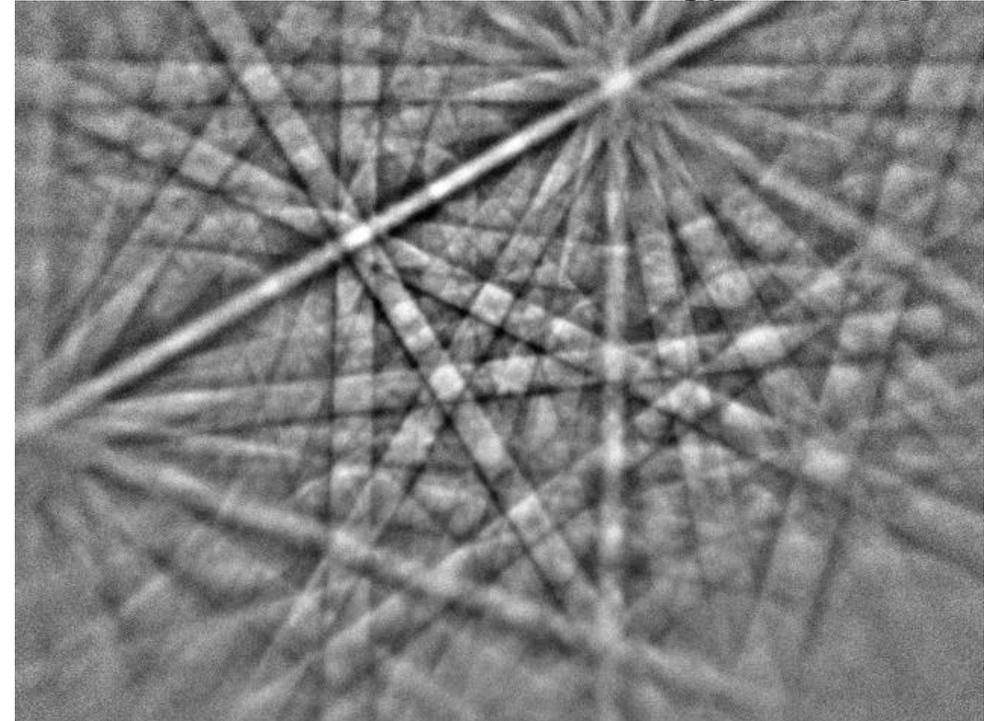


Many excellent quality patterns

Zirconium 702 alloy

Pattern quality EBSD maps




Mechanical polishing by colloidal silica

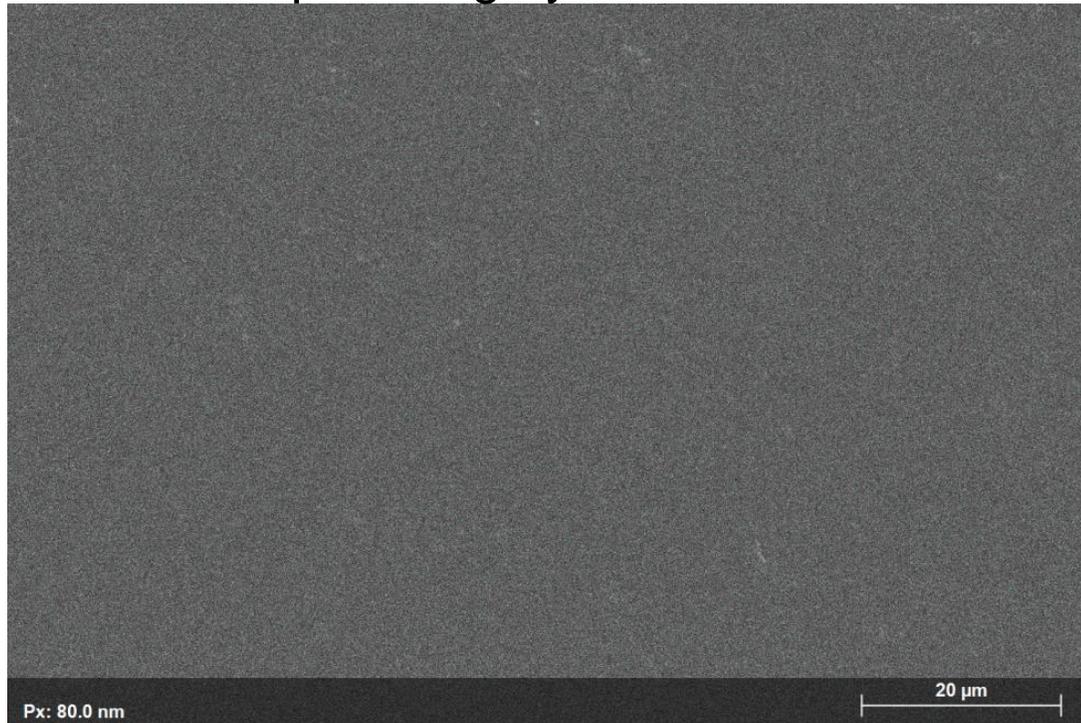

Indexing rate: < 1%

Ion milled: 8 and 2 keV energy; 3° angle

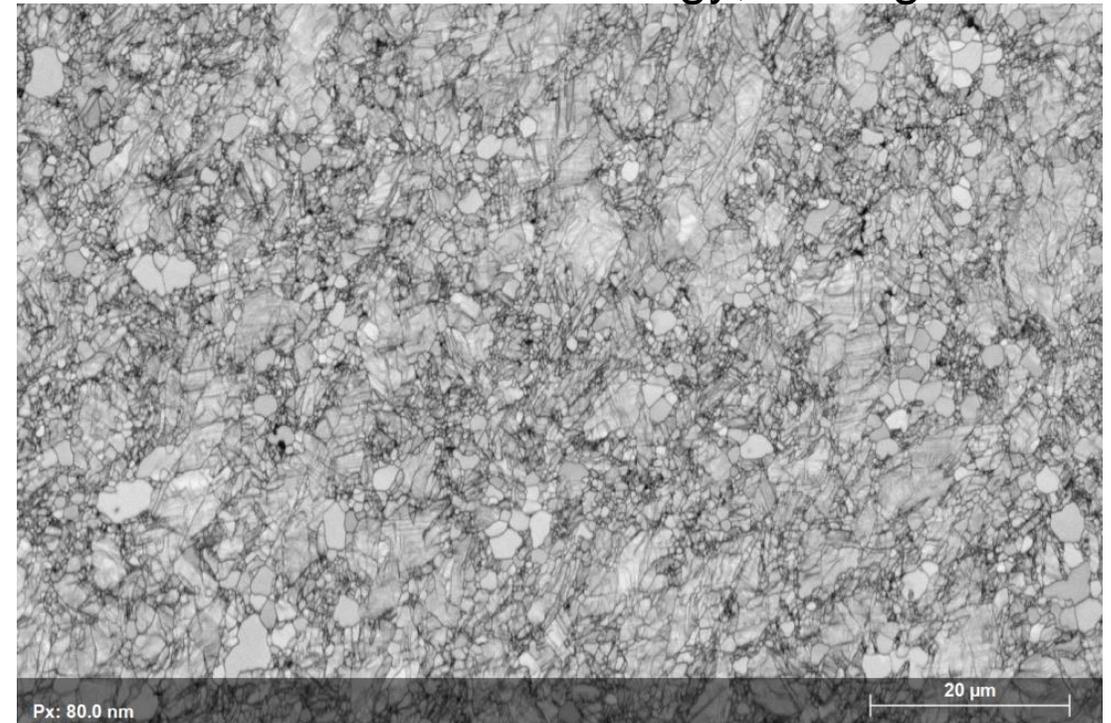


Indexing rate: 98.6%

Zirconium 702 alloy

EBSD IPF maps




Ion milled: 8 and 2 keV energy; 3° angle

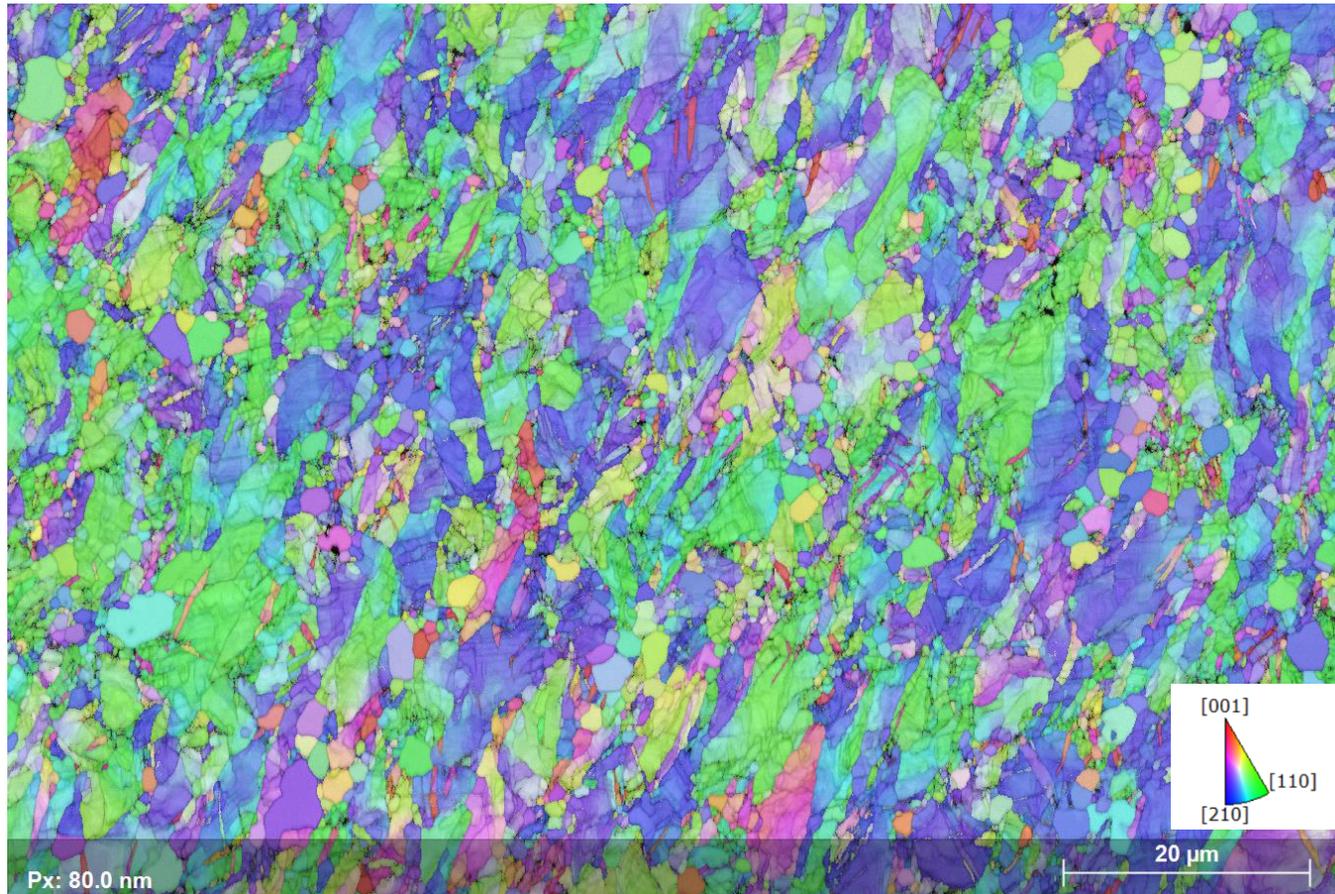

IPF Z map

Step size: 80 nm

Indexing rate: 98.6%

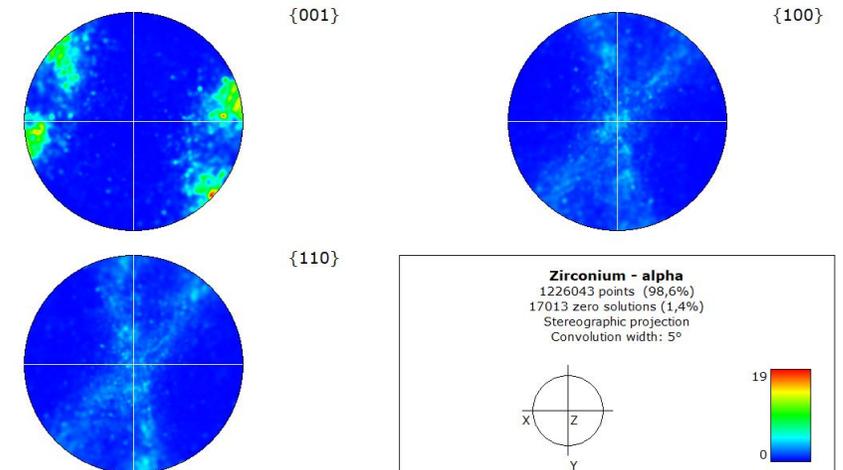


Zirconium 702 alloy


Grain average misorientation: Strain localization



Ion milled: 8 and 2 keV energy; 3° angle

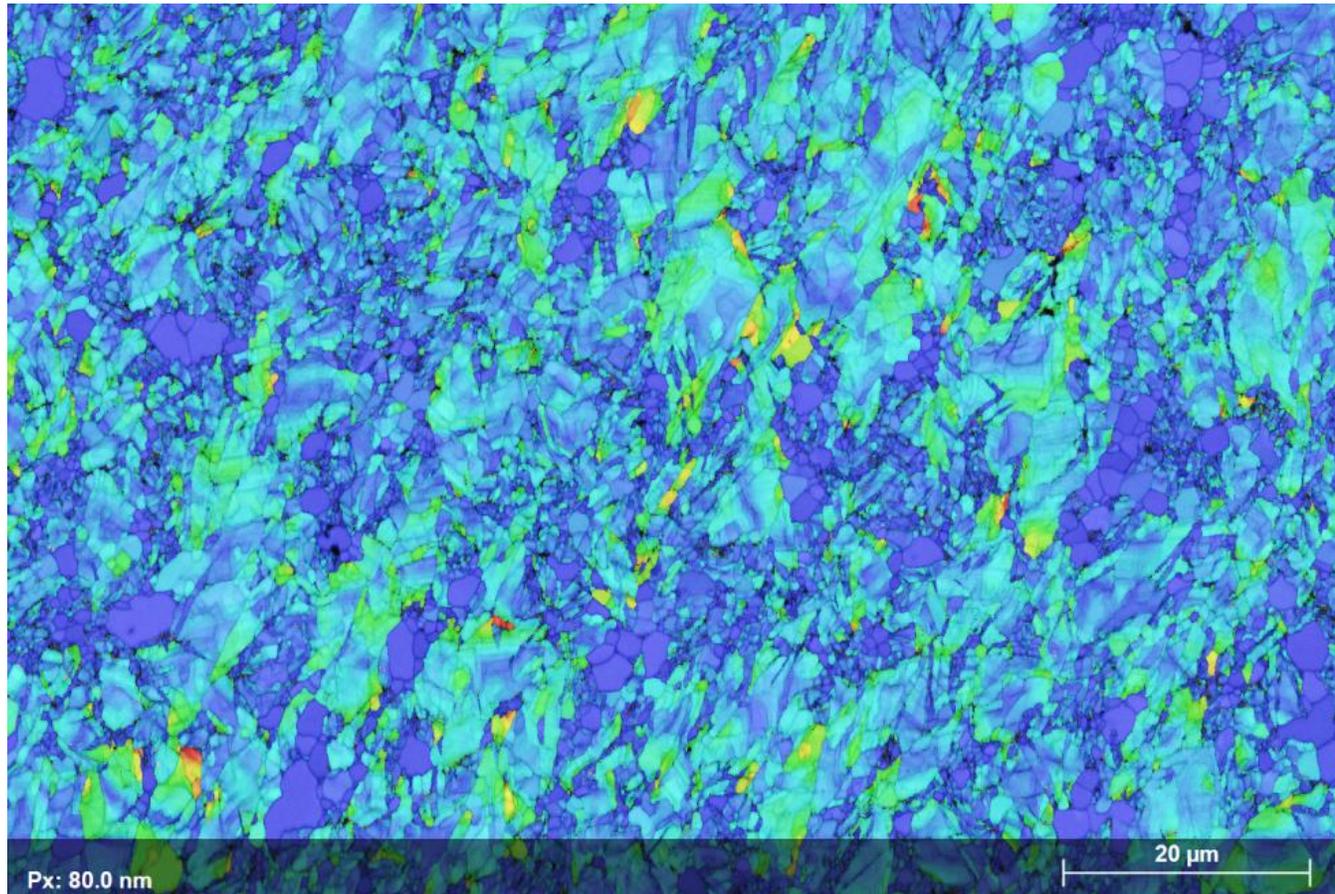

No artifact is observed after ion milling

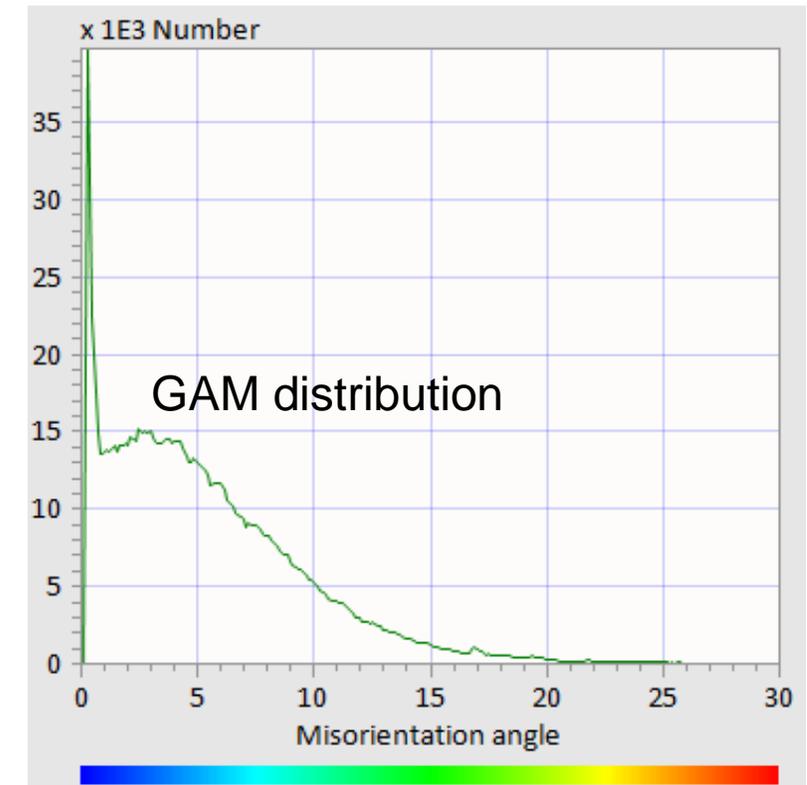


Zirconium 702 alloy


Grain average misorientation: Recrystallized grains



Ion milled: 8 and 2 keV energy; 3° angle

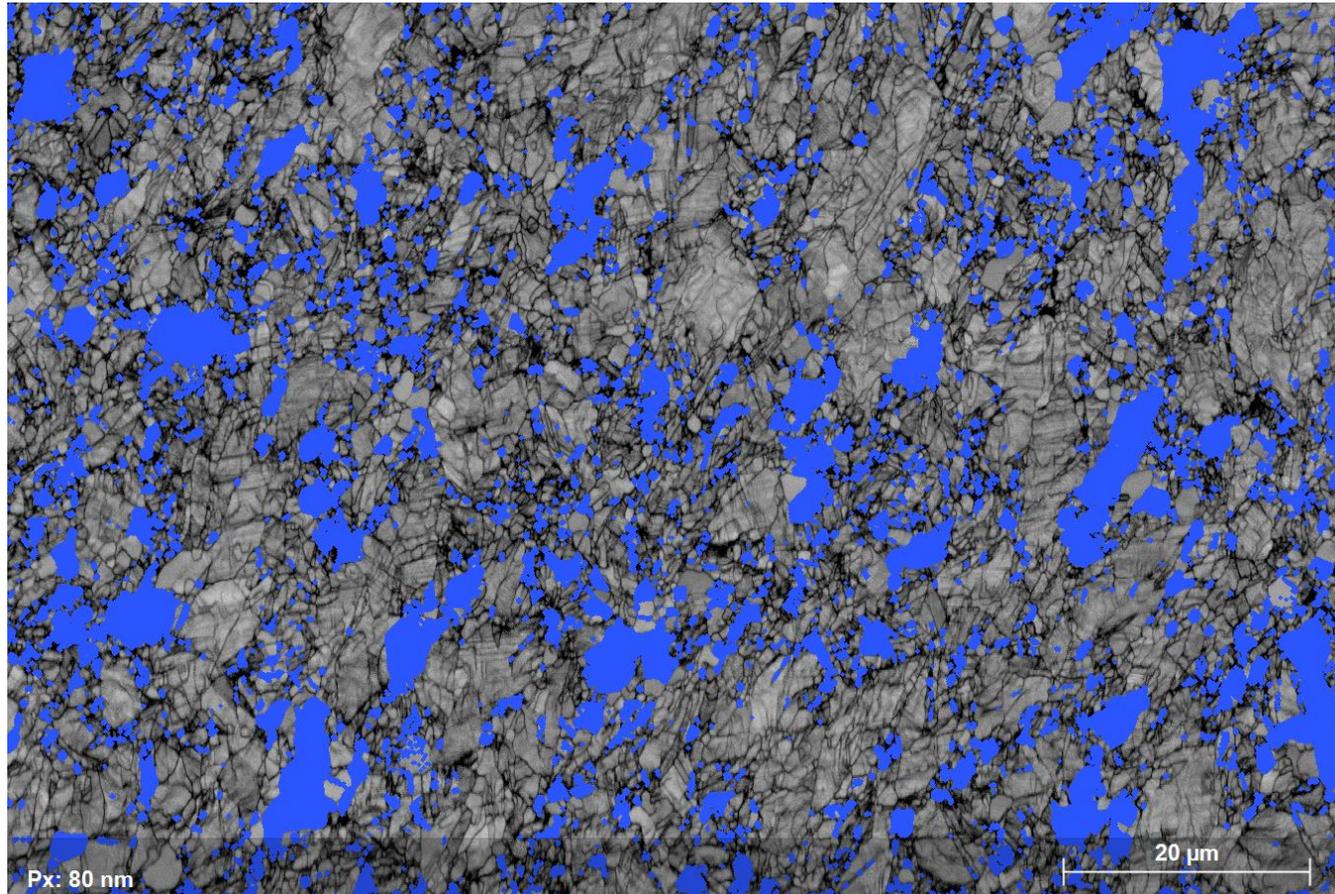

Recrystallized grains:
Misorientation angle from 0.01° to 0.6°

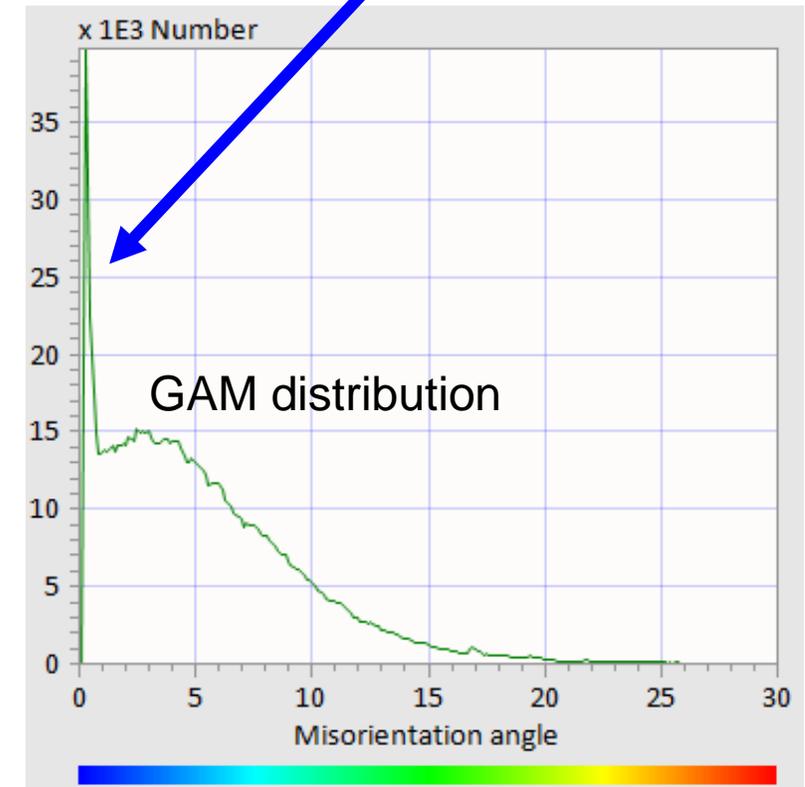


Zirconium 702 alloy


Kernel average misorientation



Ion milled: 8 and 2 keV energy; 3° angle

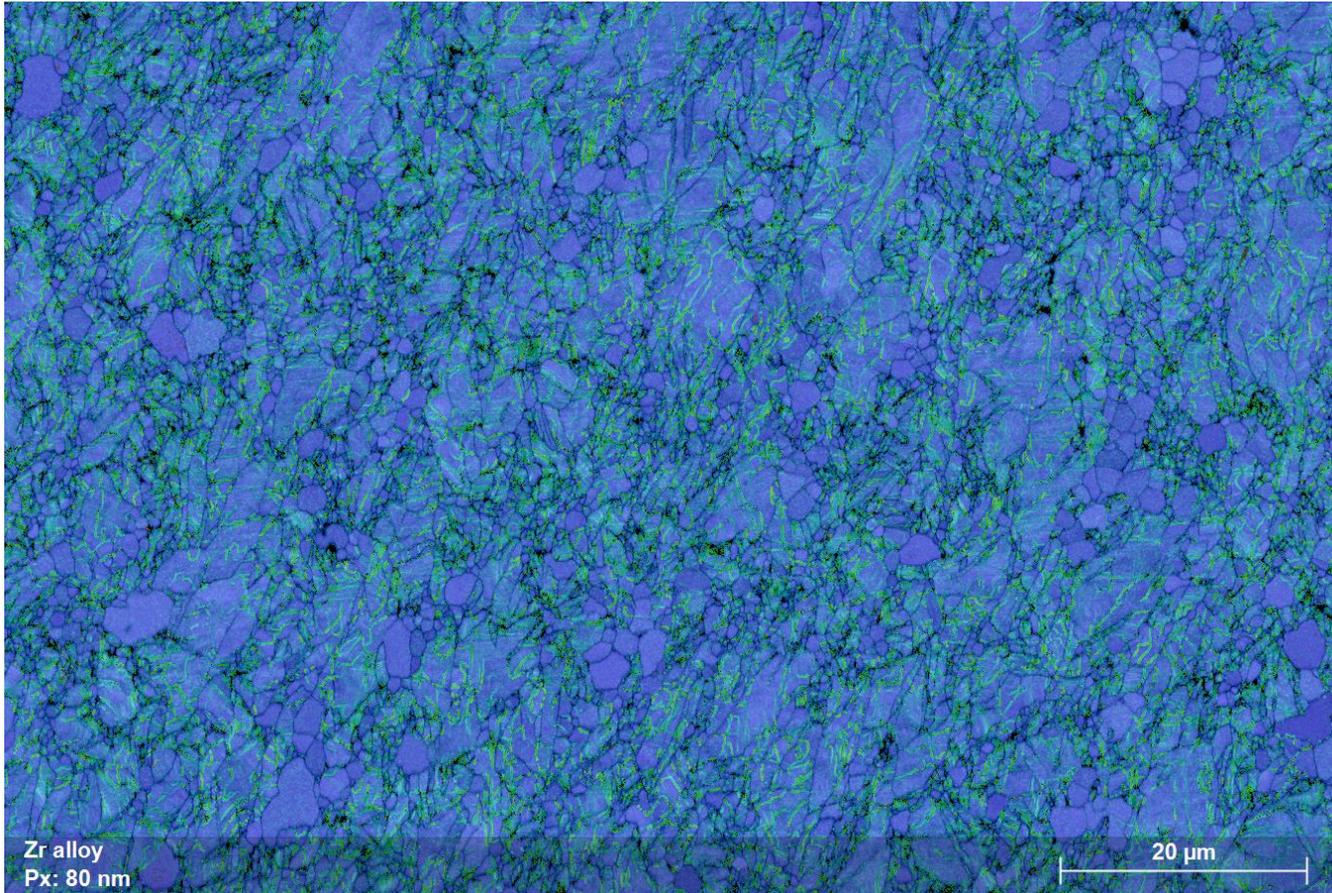

No artifact is observed after ion milling

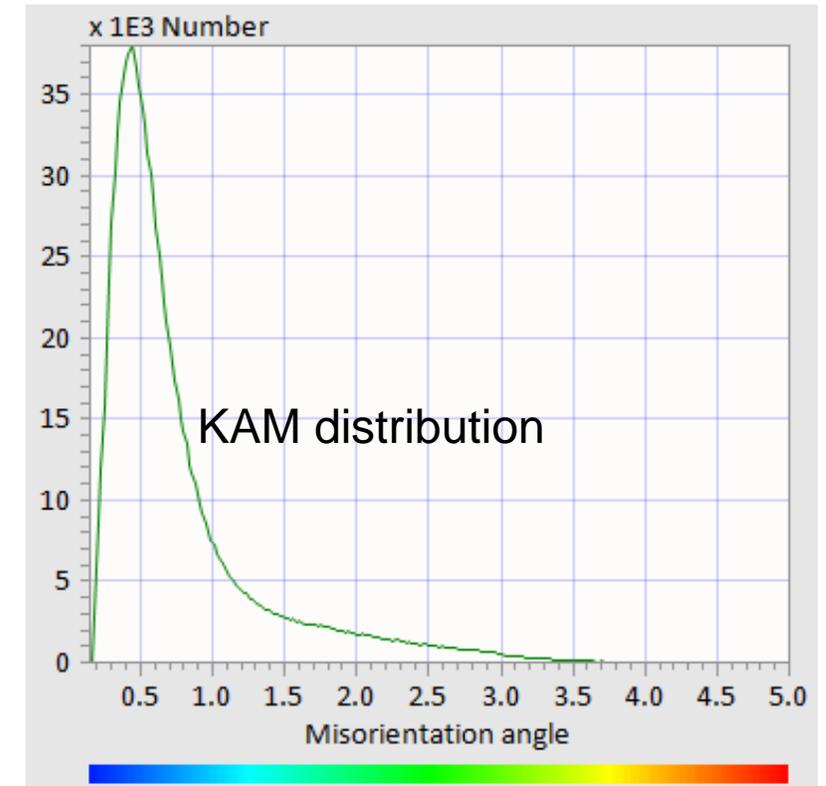



Titanium grade 2

Titanium grade 2

Orientation contrast imaging with ARGUS™



Ion milled: 8, 6, and 2 keV; 3° angle




Titanium grade 2

Orientation contrast imaging with ARGUS™



Ion milled: 8, 6, and 2 keV; 3° angle


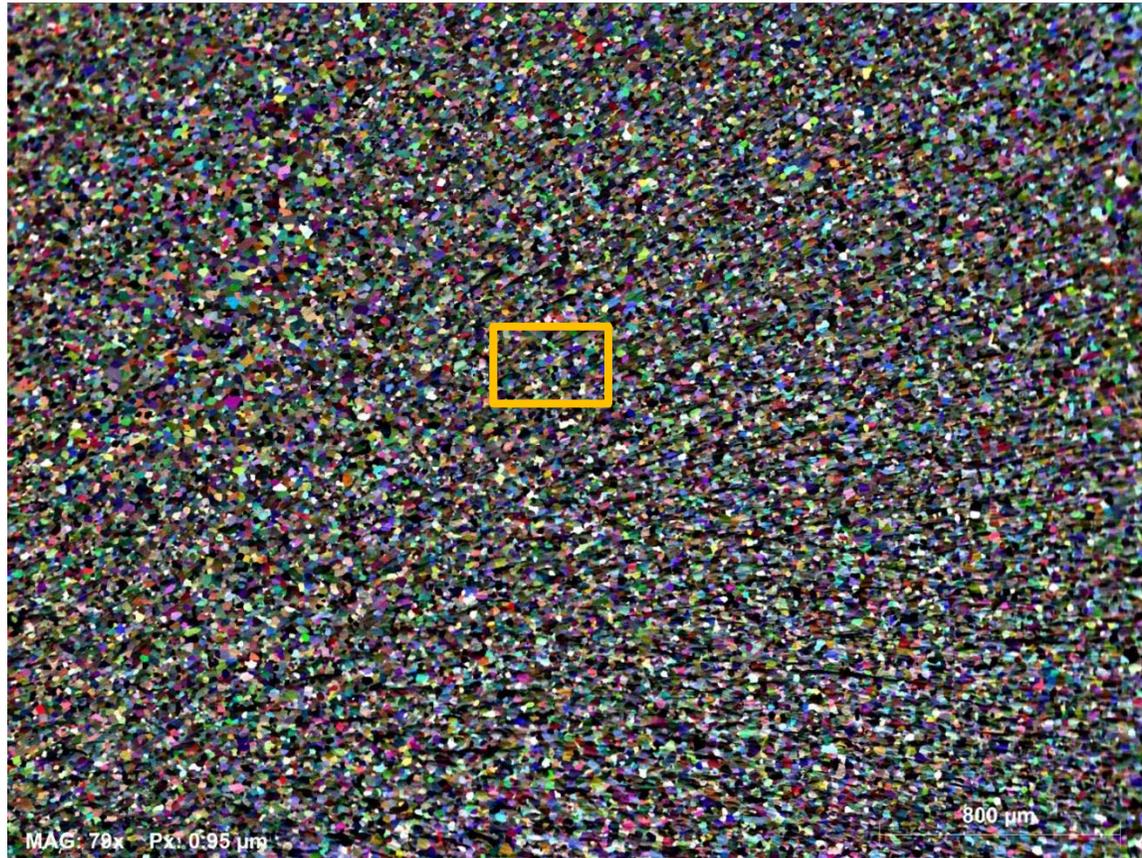


Titanium grade 2

Orientation contrast imaging with ARGUS™



Ion milled: 8, 6, and 2 keV; 3° angle

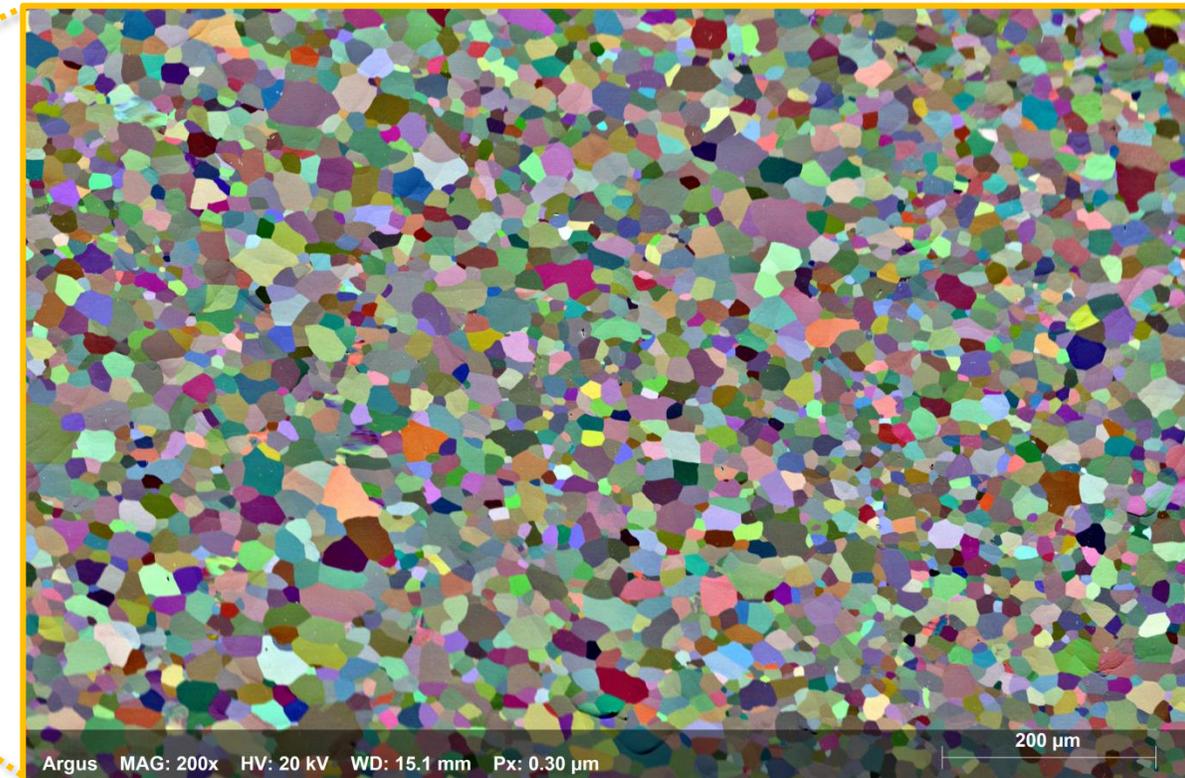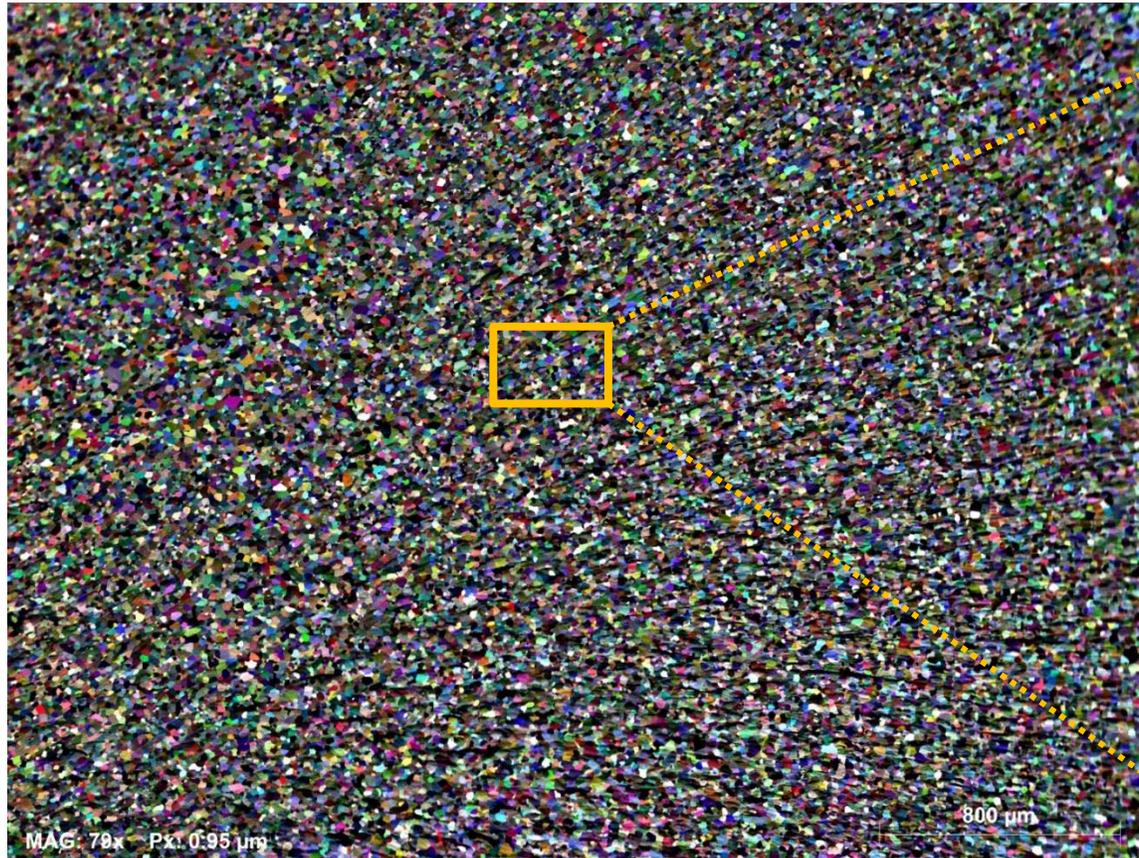

Titanium grade 2

EBSD pattern



Phase file

PHASE EDITOR

Titanium

Unit Cell Atoms

Crystal symmetry P 6₃/mmc

194 IT No.

P 6₃/m2/m2/c Space group selection

P 6₃/m2/m2/c Transform

6/m mm Laue group

Lattice parameters

2,951	a ₀ [Å]	90	α [°]
2,951	b ₀ [Å]	90	β [°]
4,685	c ₀ [Å]	120	γ [°]

Volume [Å³] : 35.333

Mass [u] : 95.76

Density [g/cm³] : 4.5

Titanium

Ti

ICSD-ID: 43733

Identifier


Chem. formula

Reference

OK Cancel

Replace old phase

Diffraction pattern

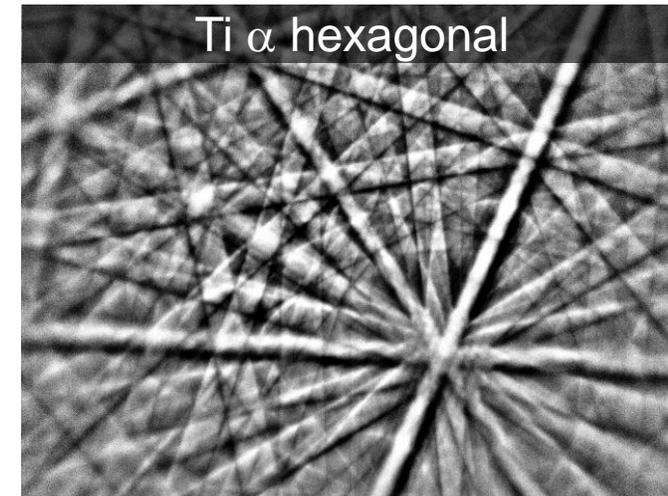


Titanium grade 2


Large-area milling and EBSD map



Ion milled: 8, 6, and 2 keV; 3° angle

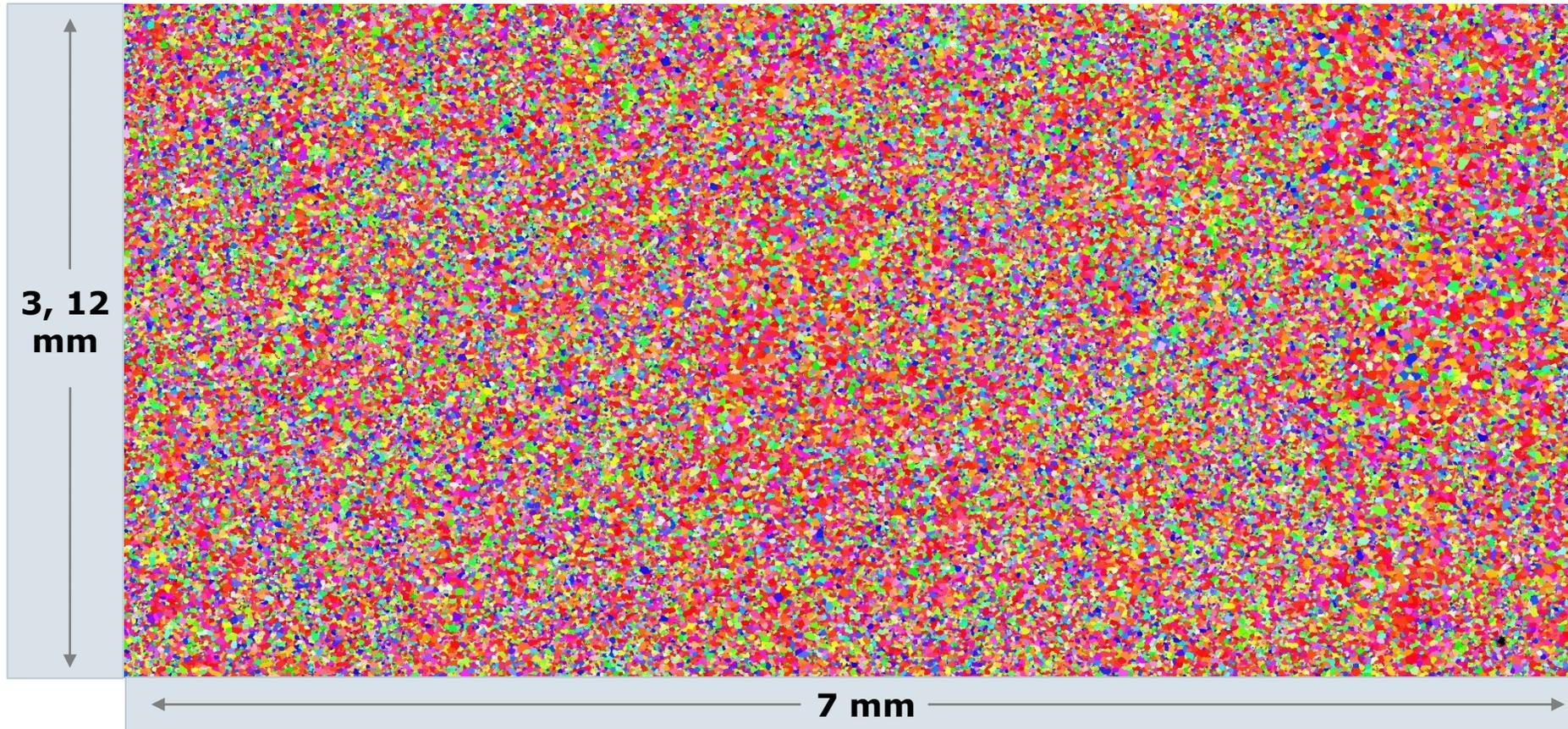

Step size: 1 μm
Indexing rate: 99%
EBSP resolution:
80x60 pixels
Measurement speed:
936 Hz

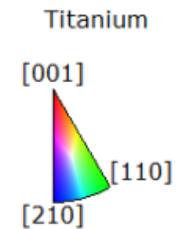


Titanium grade 2


Large-area milling and EBSD map



Ion milled: 8, 6, and 2 keV; 3° angle

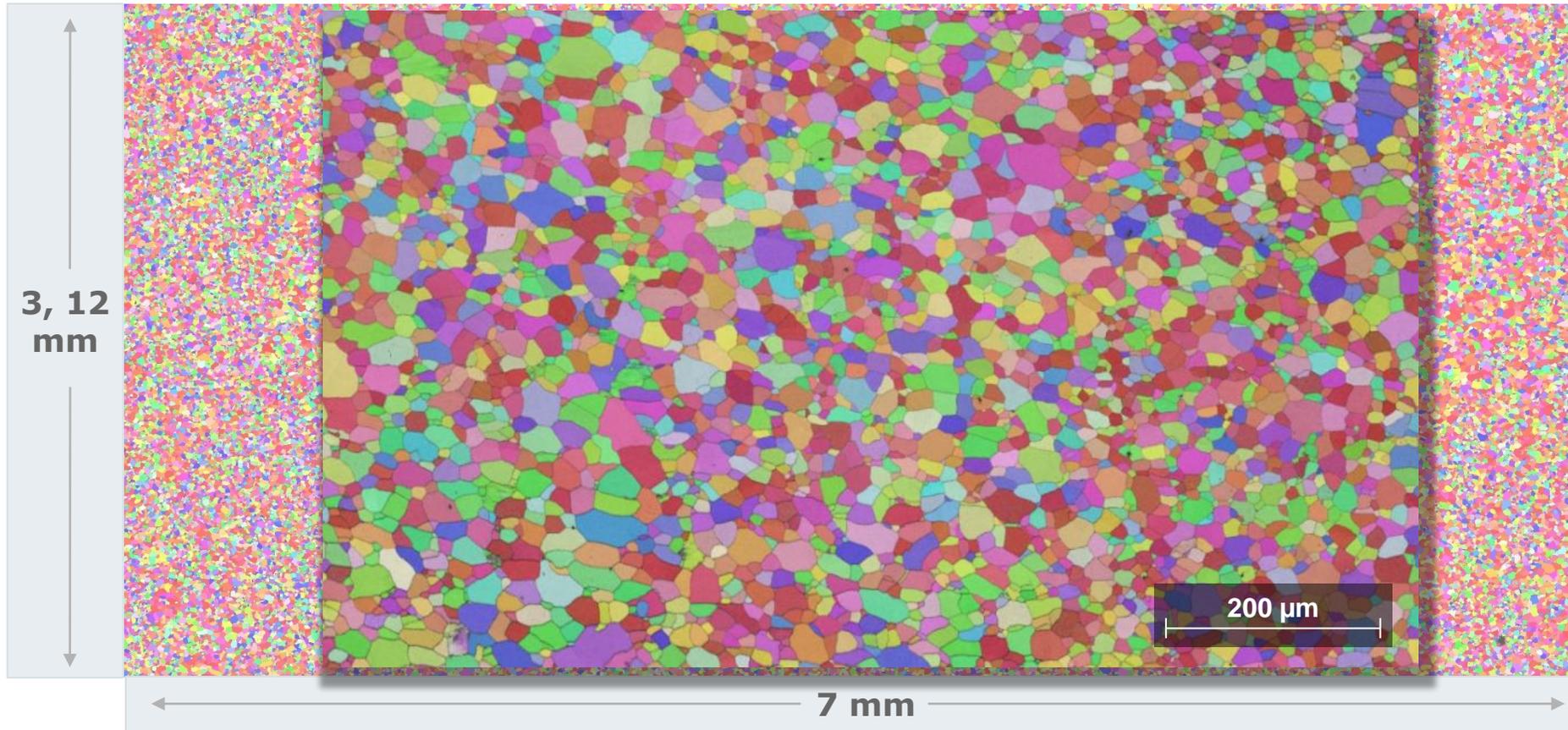

Step size: 1 μm
Indexing rate: 99%
6 min per map

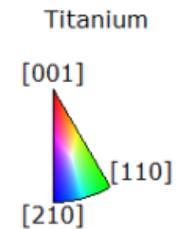


Titanium grade 2

Large-area milling and EBSD map



Grain size distribution

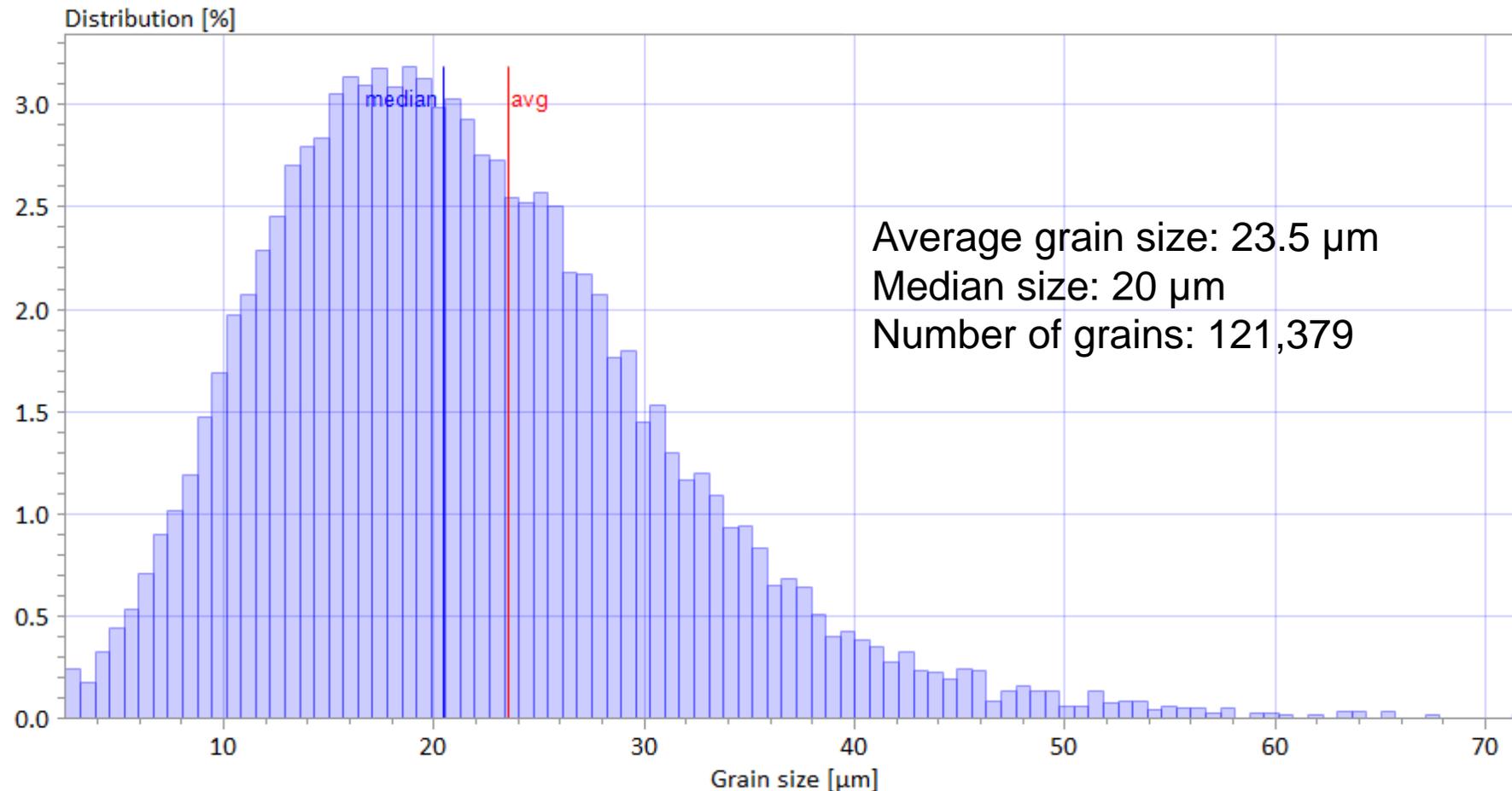




Titanium grade 5 (Ti-6Al-4V)

Titanium grade 5 (Ti-6Al-4V)

Orientation contrast imaging with ARGUS™



Ion milled: 6 and 2 keV; 3° angle

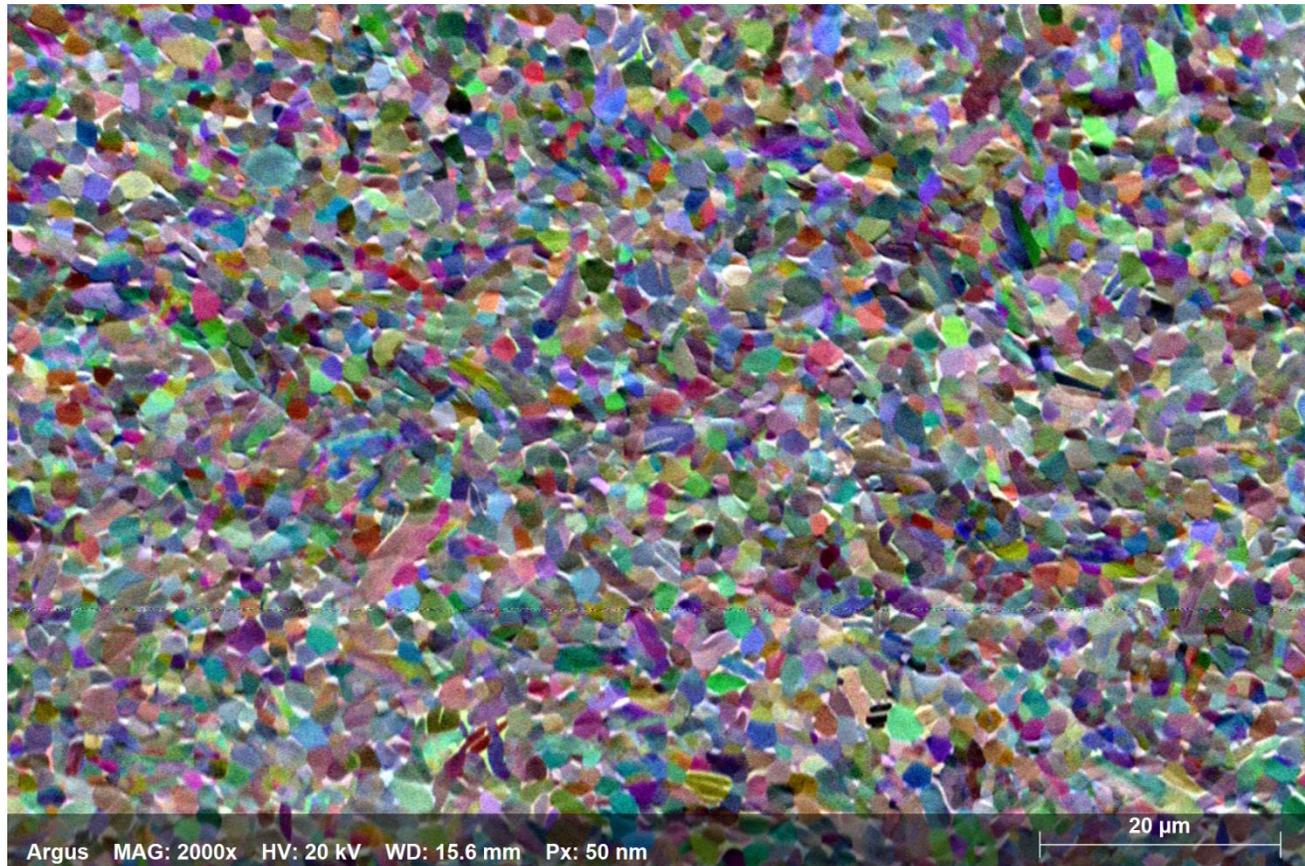

Titanium grade 5 (Ti-6Al-4V) EBSD patterns



PHASE EDITOR

Titanium

Unit Cell Atoms

Crystal symmetry $P 6_3/mmc$

IT No. 194

Space group selection $P 6_3/m2/m2/c$

Transform

Laue group $6/m\ mm$

Detect

Lattice parameters

2,951	a_0 [Å]	90	α [°]
2,951	b_0 [Å]	90	β [°]
4,685	c_0 [Å]	120	γ [°]


Volume [Å³] : 35.333
Mass [u] : 95.76
Density [g/cm³] : 4.5

Dynamical simulation

Titanium Identifier
Ti Chem. formula
ICSD-ID: 43733 Reference

OK Cancel

Replace old phase

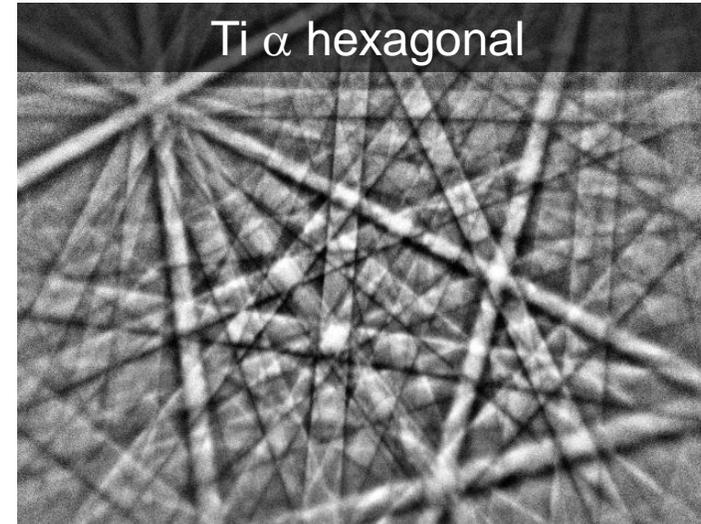

PHASE EDITOR

Titanium - Beta, Ht

Unit Cell Atoms

Crystal symmetry $I m\ 3m$

IT No. 229

Space group selection $I 4/m\ 32/m$

Transform

Laue group $m\ 3m$

Detect

Lattice parameters

3,311	a_0 [Å]	90	α [°]
3,311	b_0 [Å]	90	β [°]
3,311	c_0 [Å]	90	γ [°]


Volume [Å³] : 36.298
Mass [u] : 95.76
Density [g/cm³] : 4.381

Dynamical simulation

Titanium - Beta, Ht Identifier
Ti Chem. formula
ICSD-ID: 44391 Reference

OK Cancel

Replace old phase

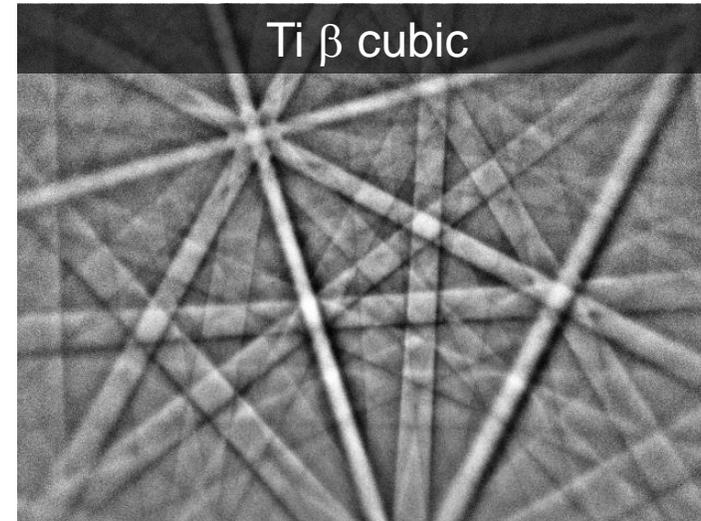


Titanium grade 5 (Ti-6Al-4V)

EBSD phase distribution map




Ion milled: 6 and 2 keV; 3° angle

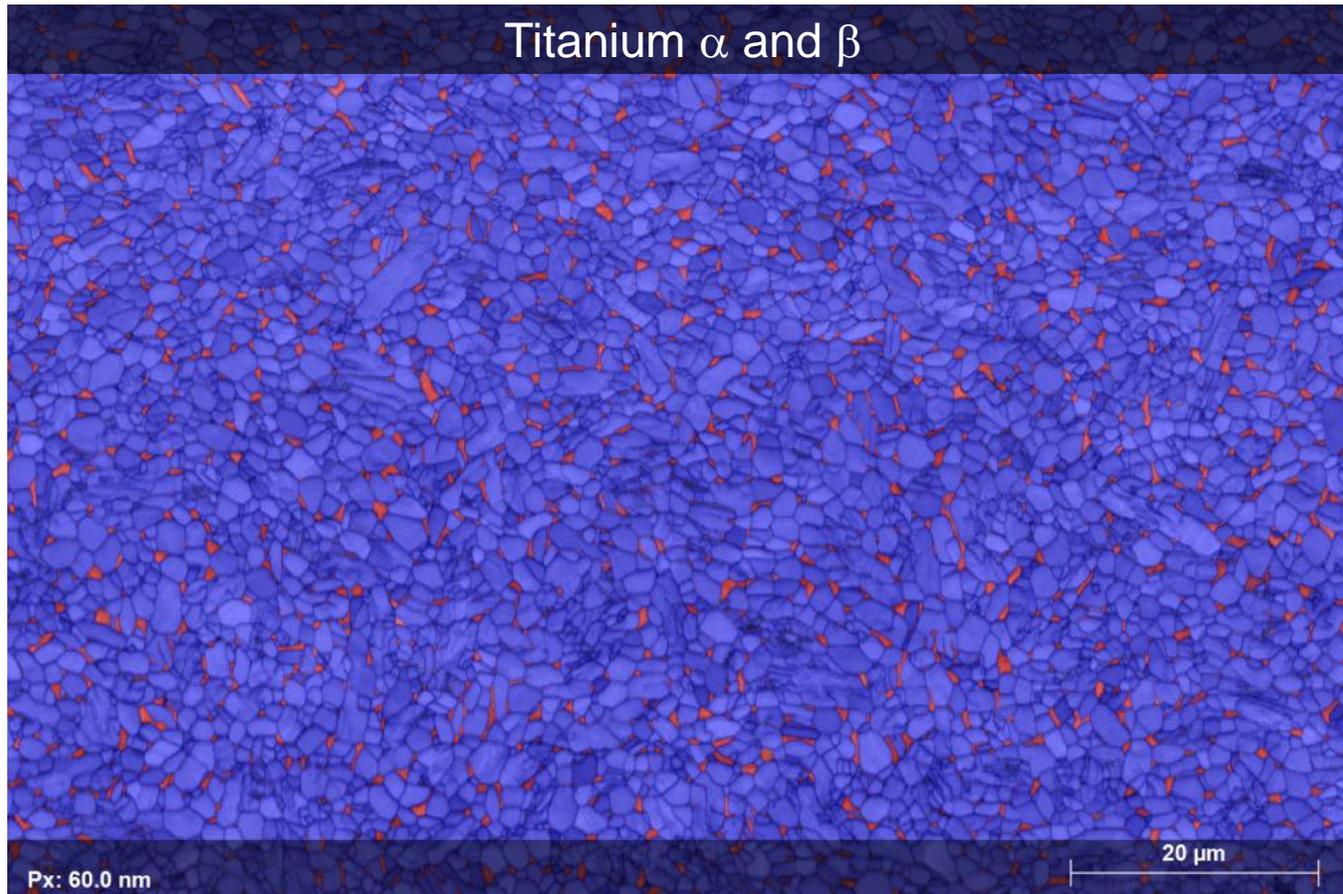

Phase map

Step size: 80 nm

Indexing rate: 98.6%


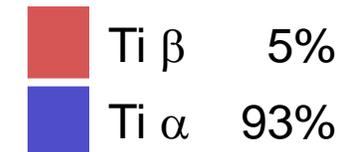


Titanium grade 5 (Ti-6Al-4V)

EBSD IPF map



Ion milled: 6 and 2 keV; 3° angle

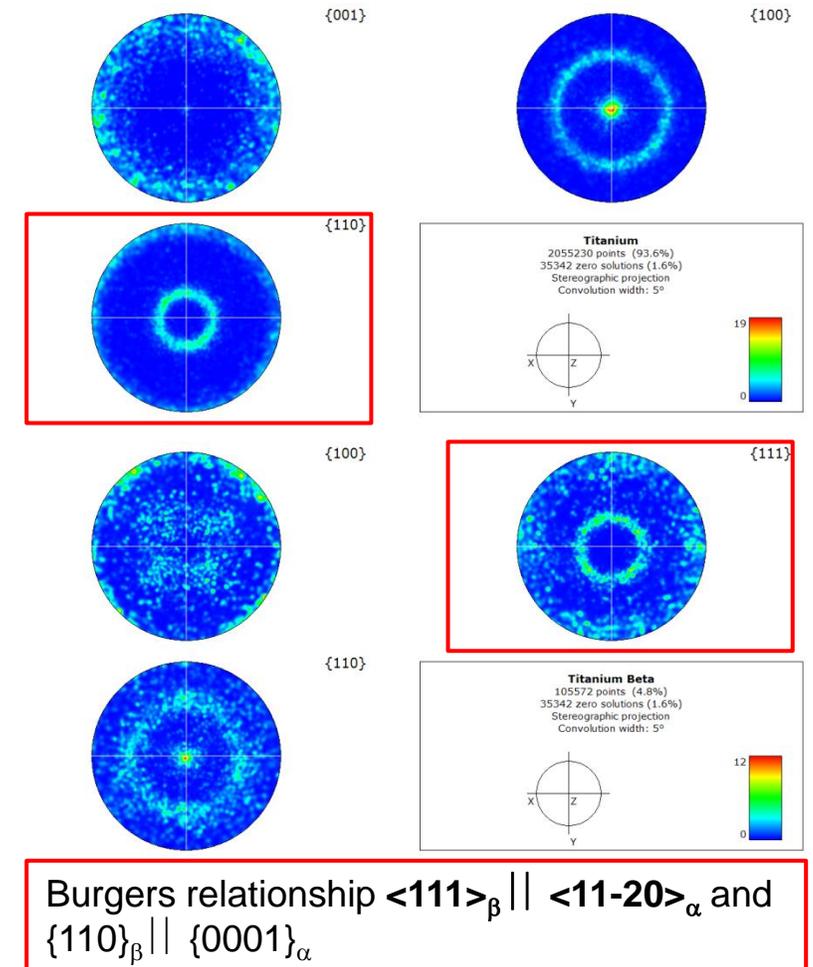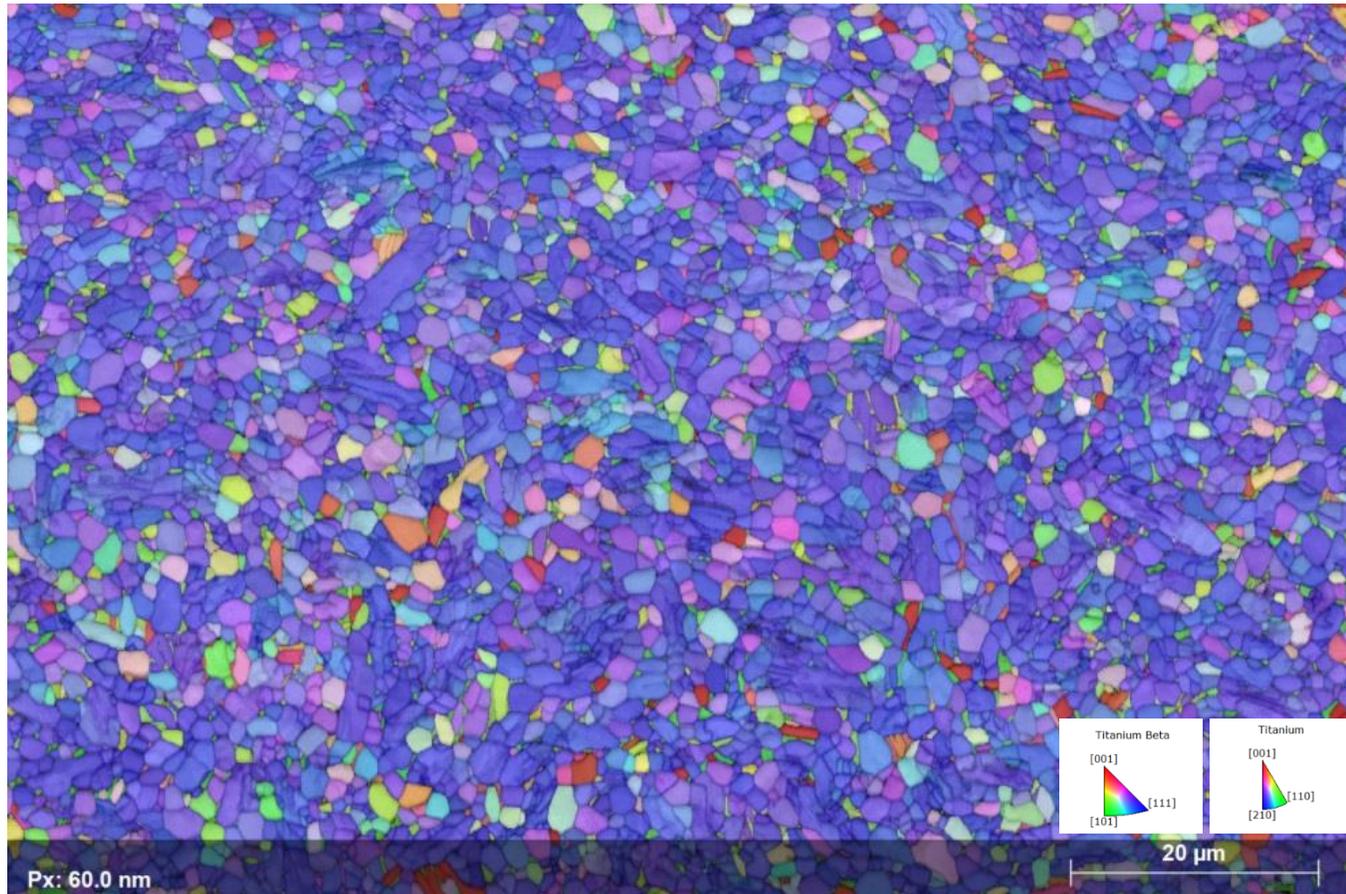




Solder bumps

Solder bumps

Cross-section sample schematic

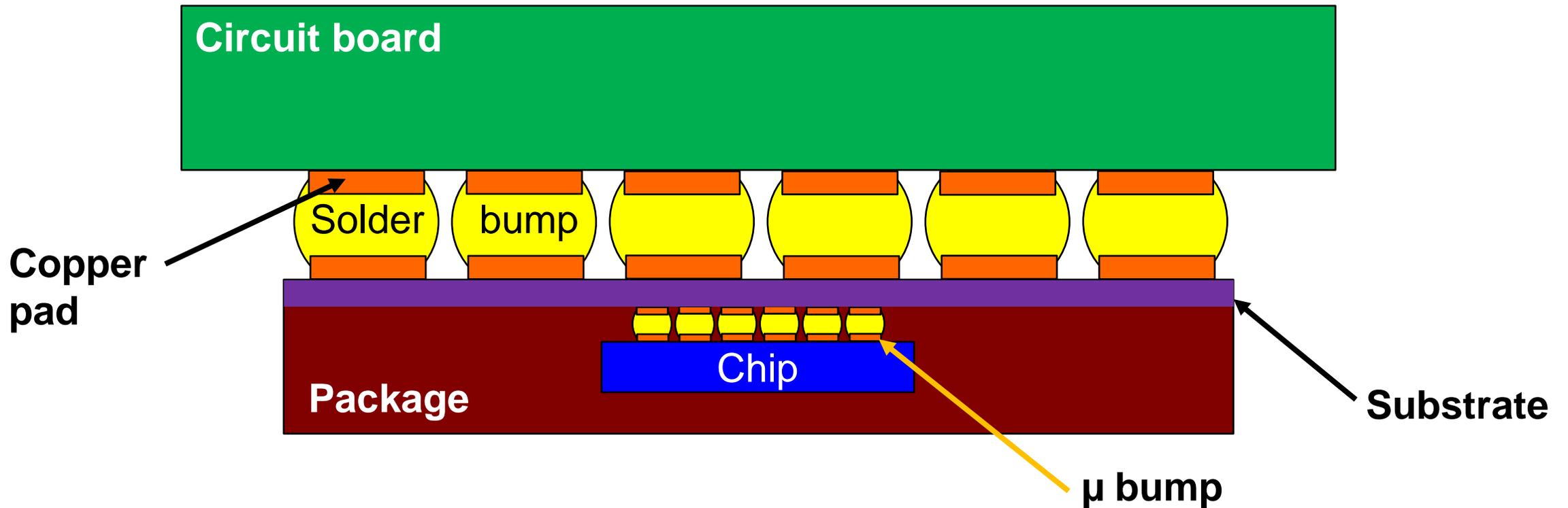


Structure is fragile

Solder bumps


EBSD IPF map

Mechanical polishing by colloidal silica

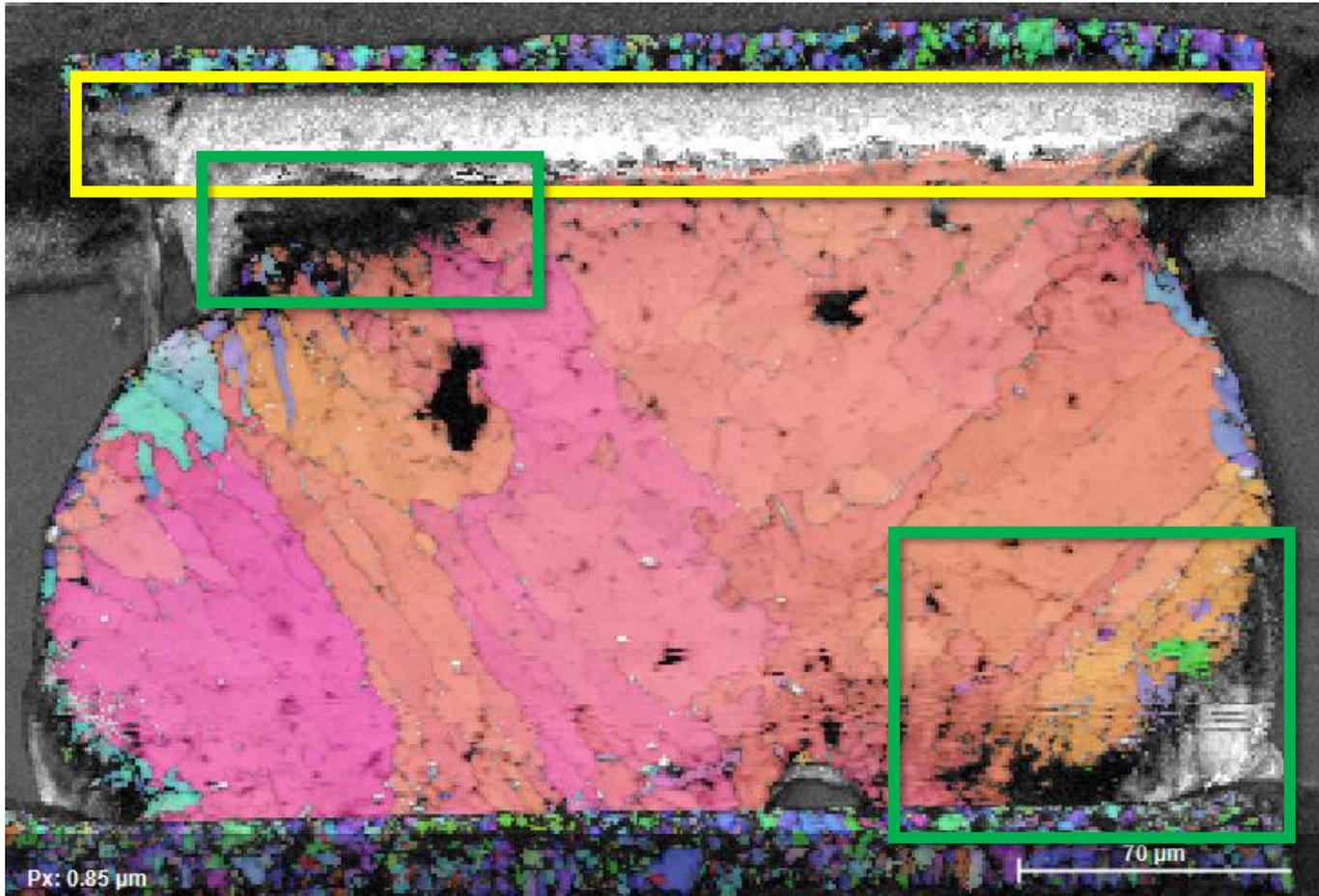

IPF Z map

- SiO_2 particles are difficult to remove
- Areas of the interface are not visible due to shadowing by the copper pad caused by differential material removal
- Corrosion of packaging components leaves surface contamination

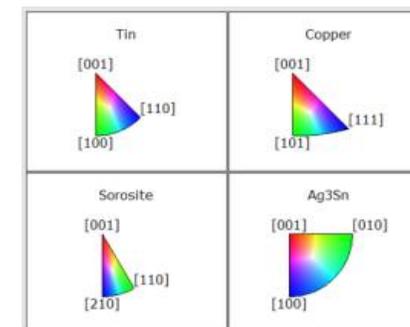


Solder bumps

Cross-section sample



Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

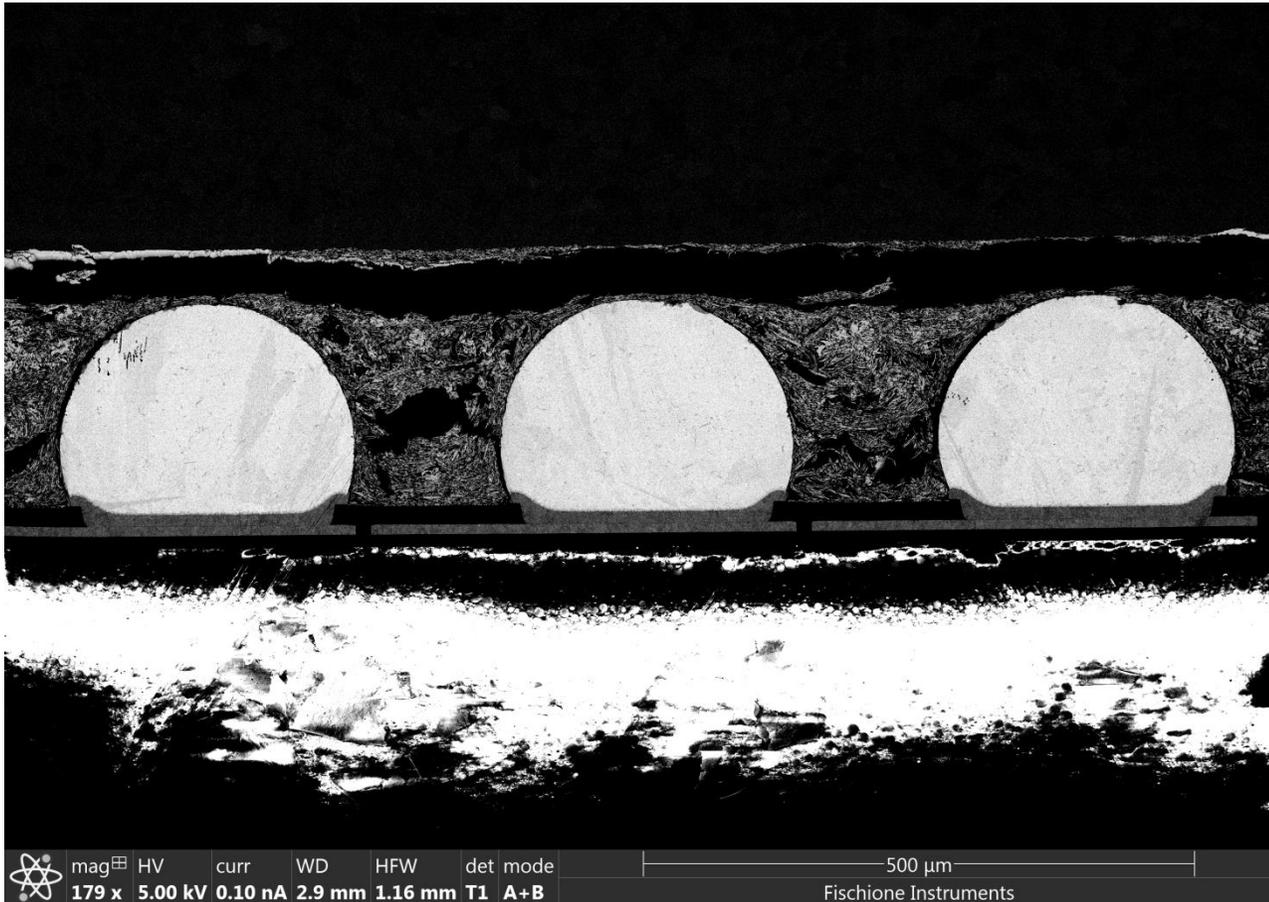


Solder bumps

Backscatter electron contrast imaging







Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

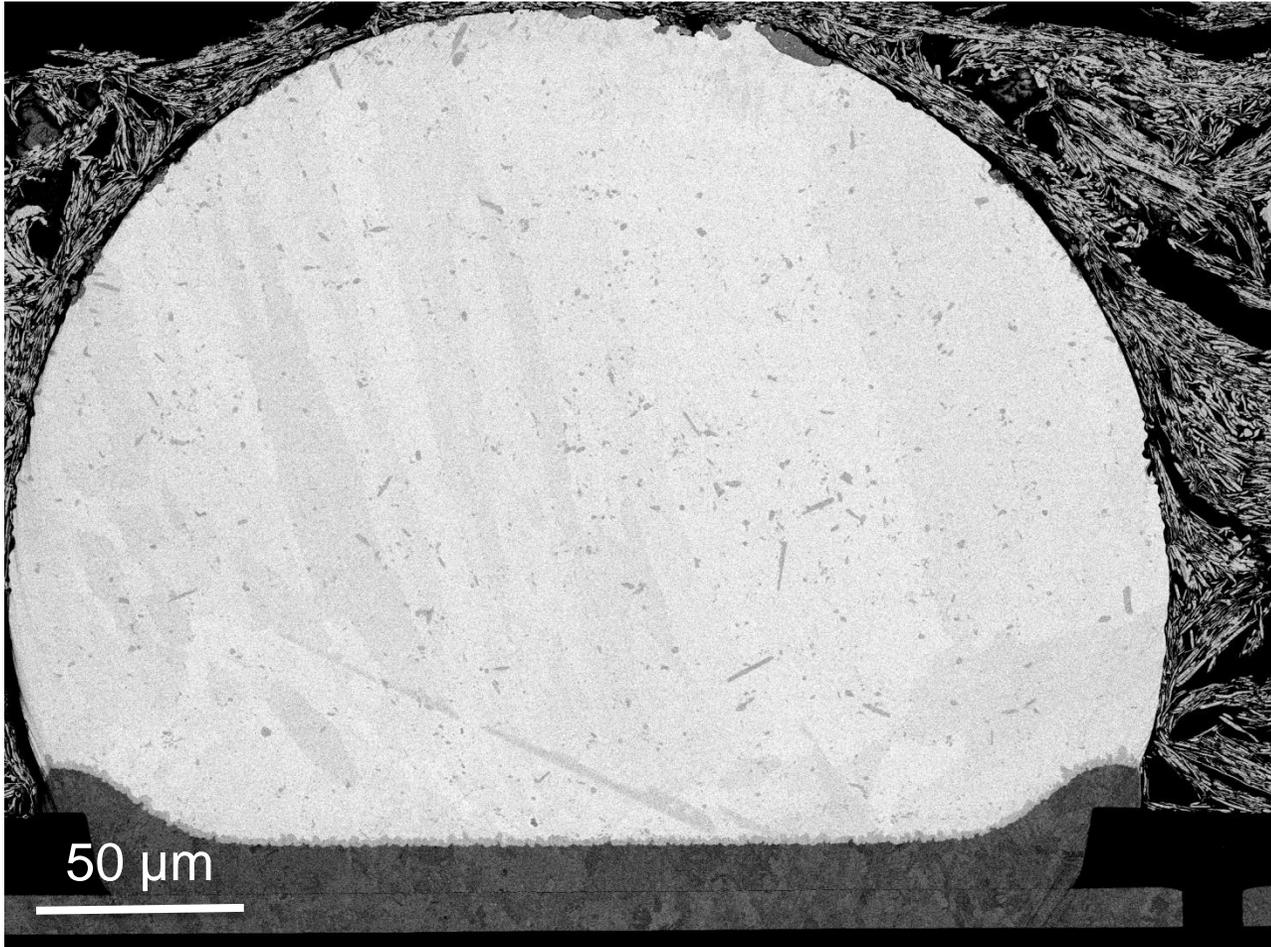

Solder bumps

EBSD patterns

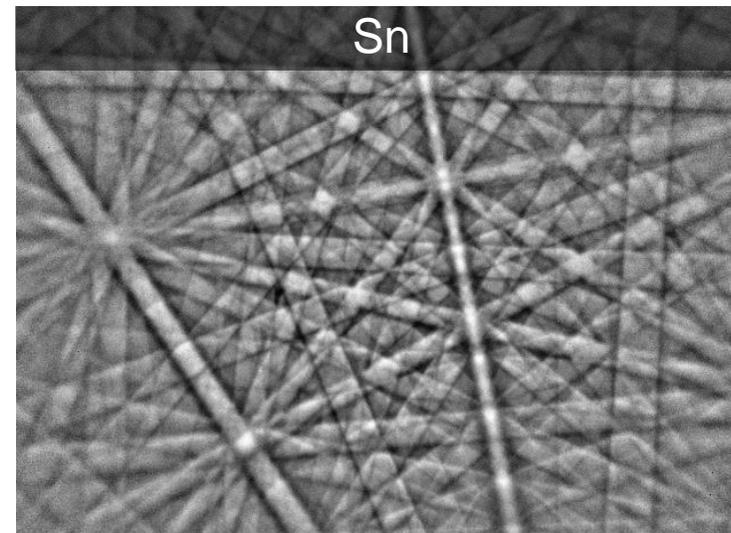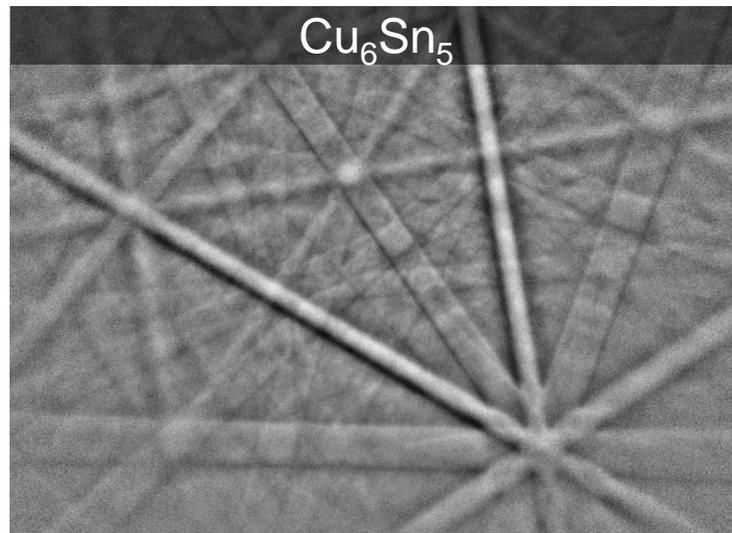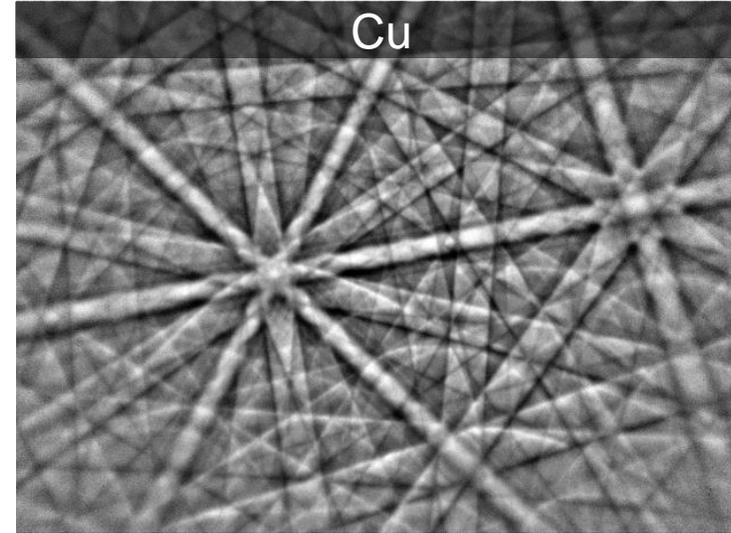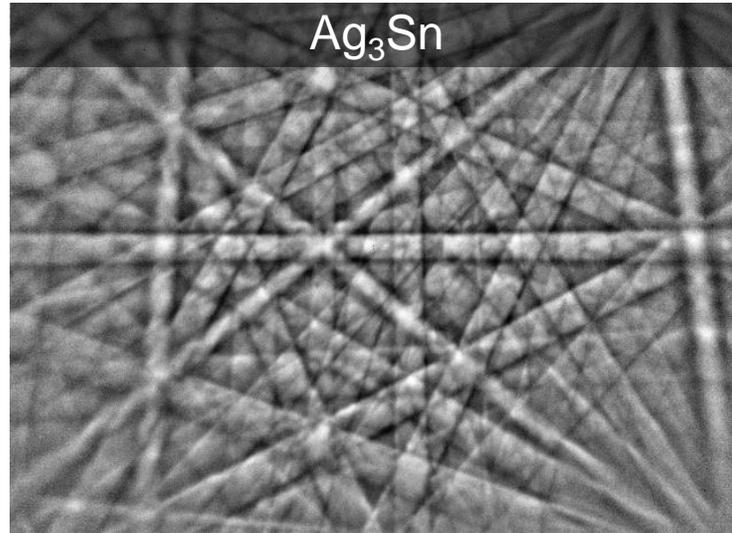


Solder bumps

EBSD phase distribution map



Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

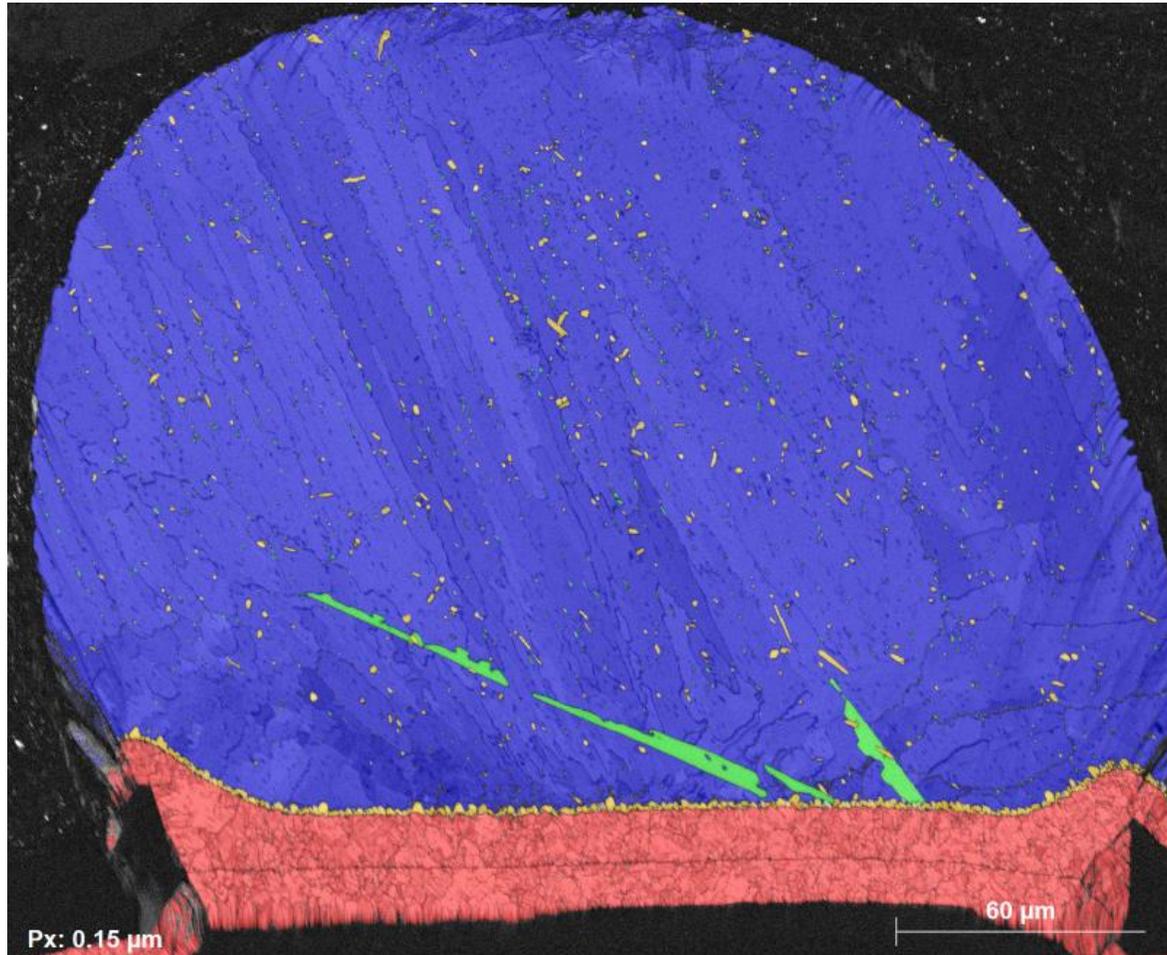

Phase map

Step size: 150 nm

Pattern resolution: 80x60 pixels


Indexing rate: 95%

Phase	Color	% normalized	Average grain size
Cu	Red	12.5	10 μm
Cu ₆ Sn ₅	Yellow	1.5	1,7μm
Ag ₃ Sn	Green	1	700 nm (fine grains only)
Sn	Blue	85	NA (only 4 grains)


Solder bumps

EBSD IPF map

Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

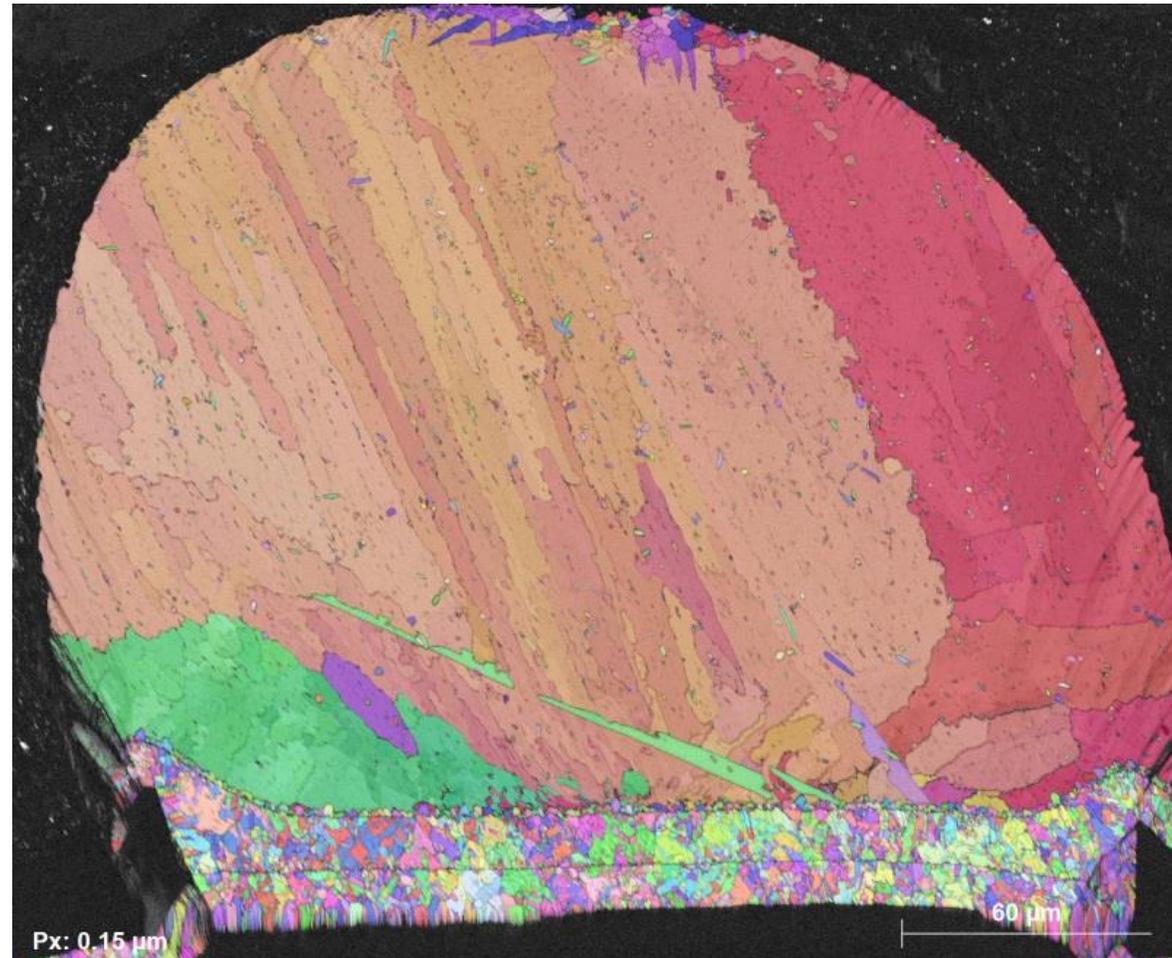

IPF Z map


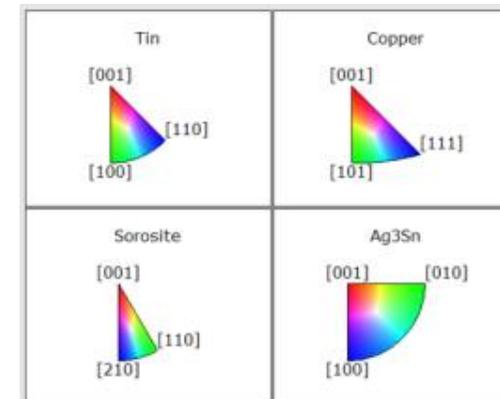


Solder bumps

Kernel average misorientation



Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

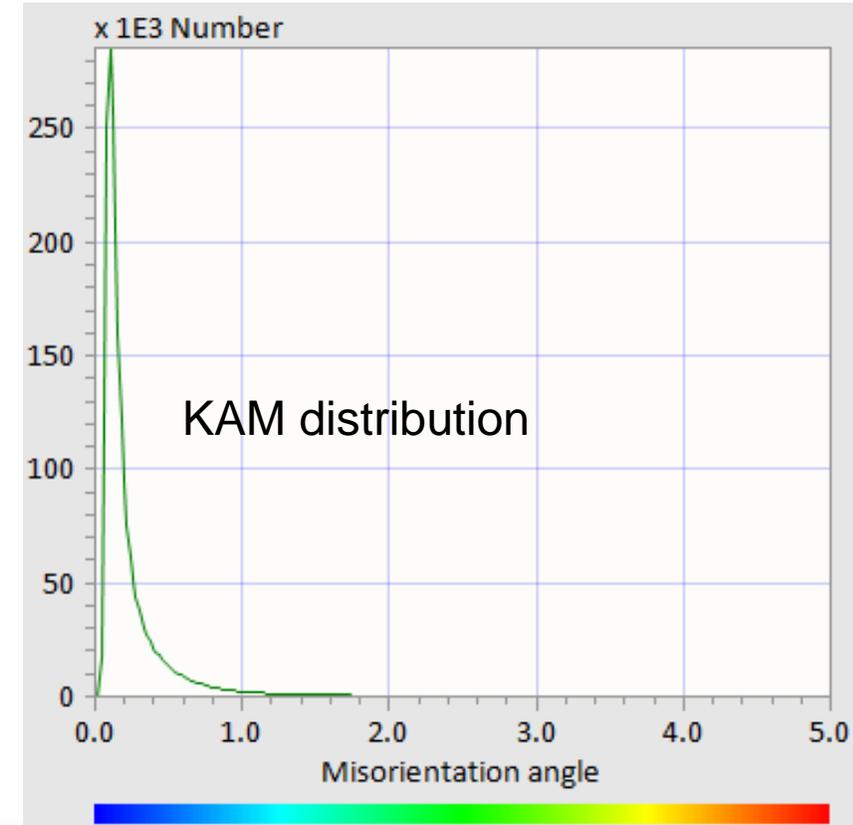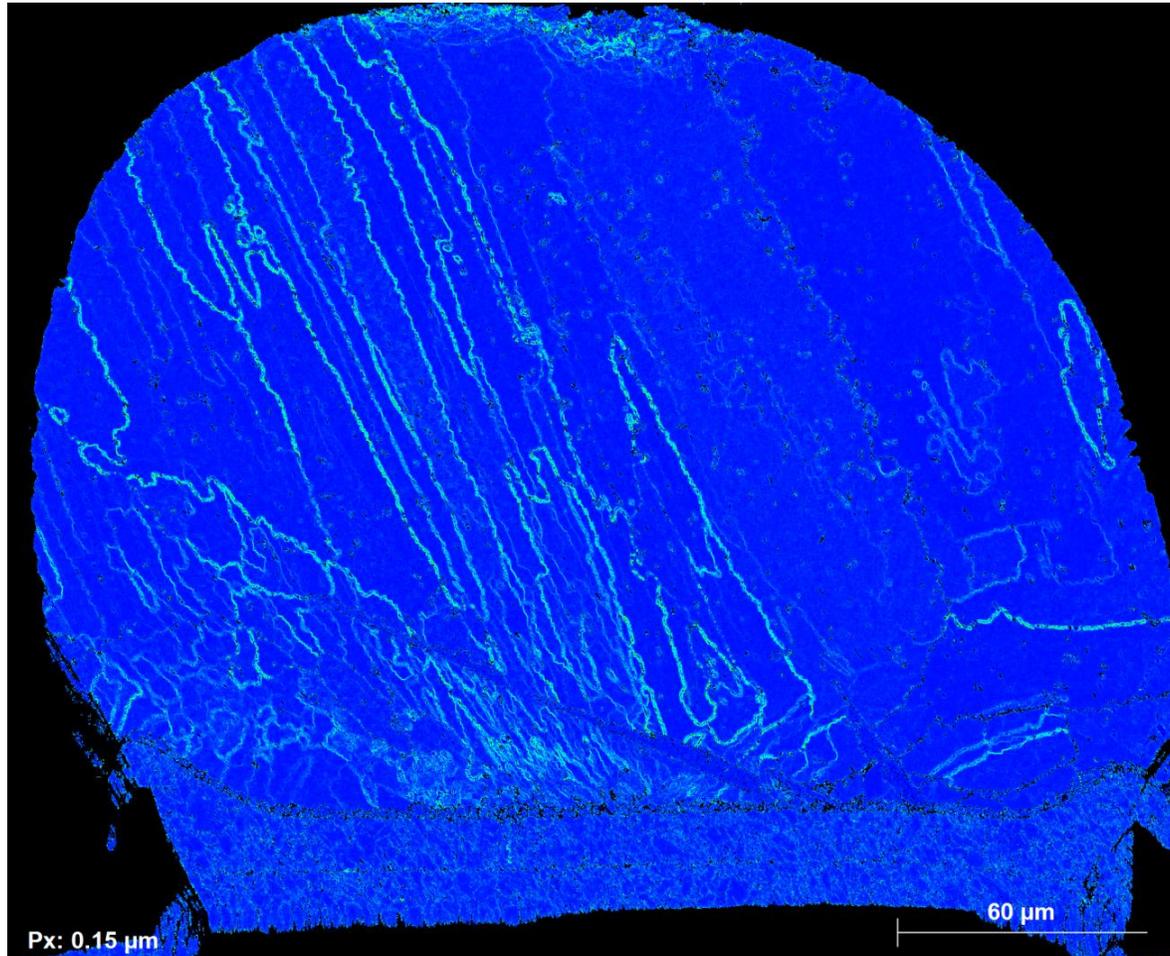


Solder bumps

Orientation contrast imaging with ARGUS™



Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

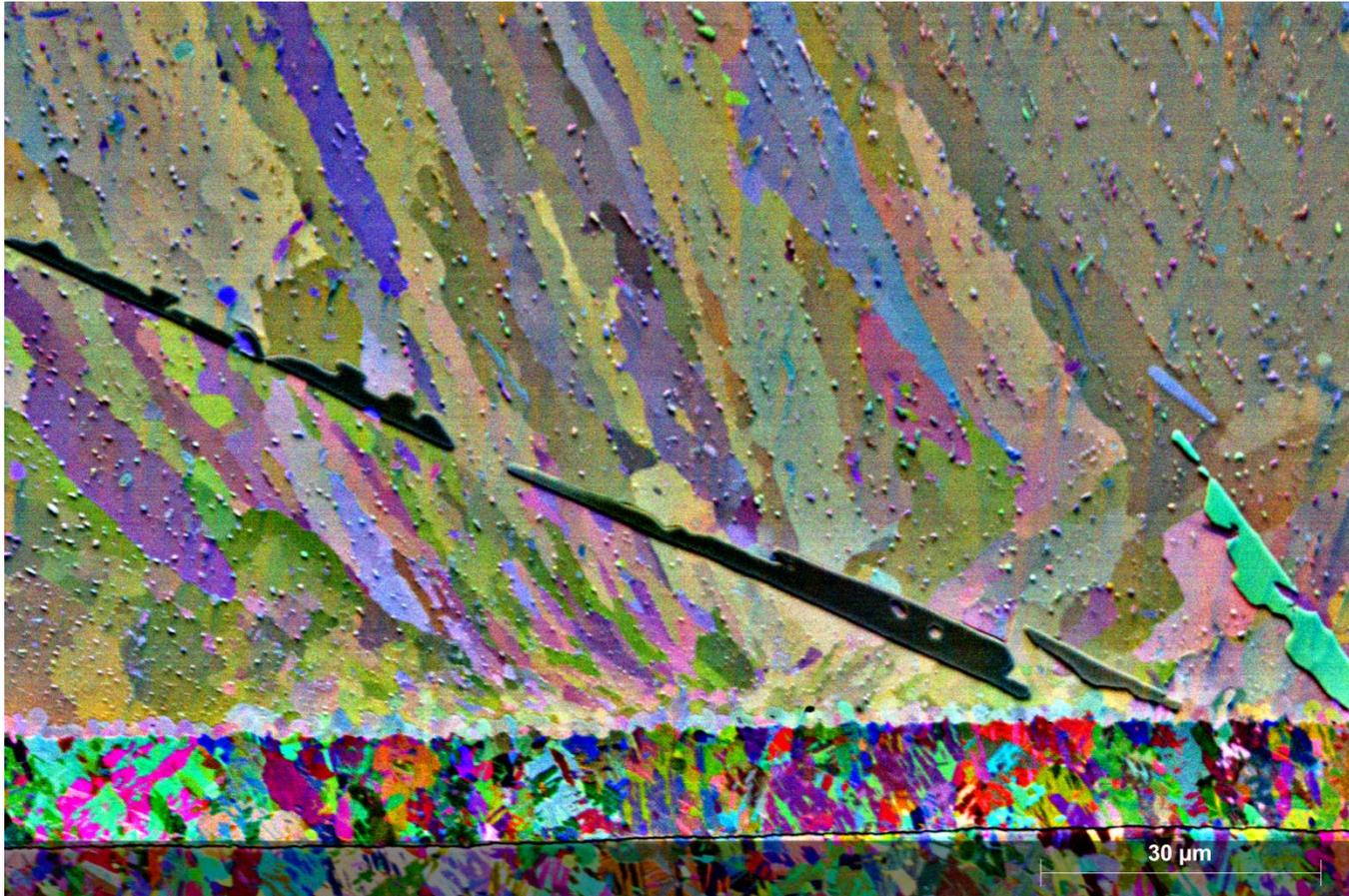


Solder bumps

EBSD phase distribution map



Ion milled at 6 and 2 keV; cross-section milling with mask under **cryo conditions**

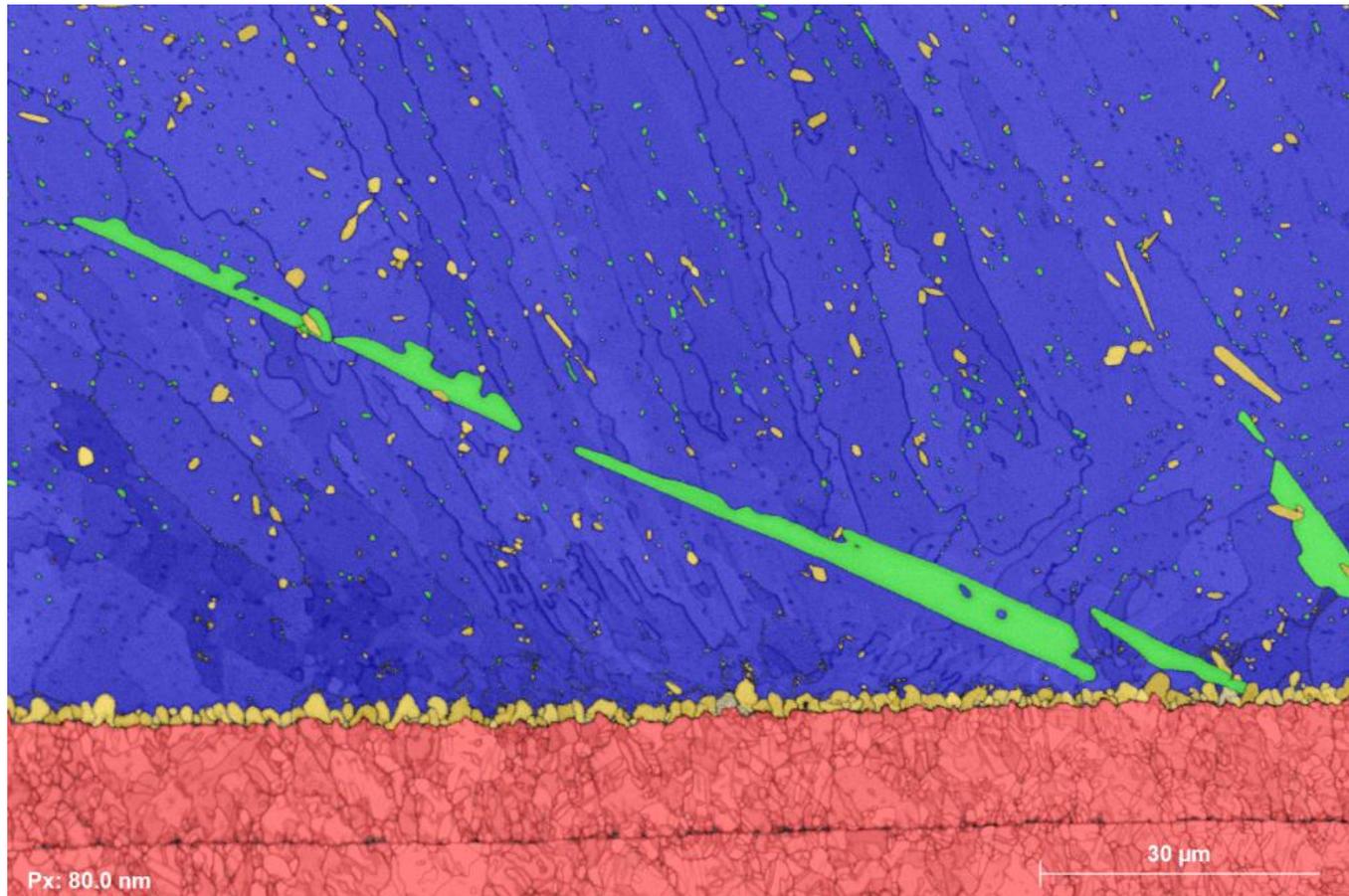


Phase map

Step size: 80 nm

Pattern resolution: 80x60 pixels

Indexing rate: 96%

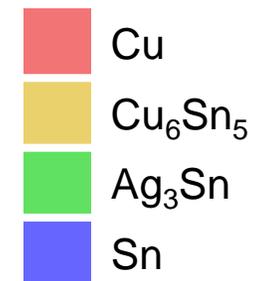

Achieve excellence in EBSD analyses

Summary




Combining Fischione Instruments and Bruker NanoAnalytics technologies allows all user levels to quantitatively characterize a sample's microstructure in a timely and accurate manner.

- **Fischione Instruments** provides fast and reliable sample preparation solutions for a wide range of applications and a wide range of challenging materials, such as thermally sensitive samples (solder bumps) and environmentally sensitive samples (lithium-ion batteries)
- **Bruker QUANTAX EBSD** delivers a complete solution for EBSD and TKD, with:
 - ARGUS™ imaging system with unmatched sensitivity and speed for sample preparation quality control and qualitative mapping of plastic and elastic strain in the specimen
 - high resolution EBSD measurement to characterize quantitatively the micro/nano-structure, phase distribution, texture,...., making it the ideal analytical tool for deformation studies



Are there any questions?

Please type your questions
in the Q&A box and press *Send*

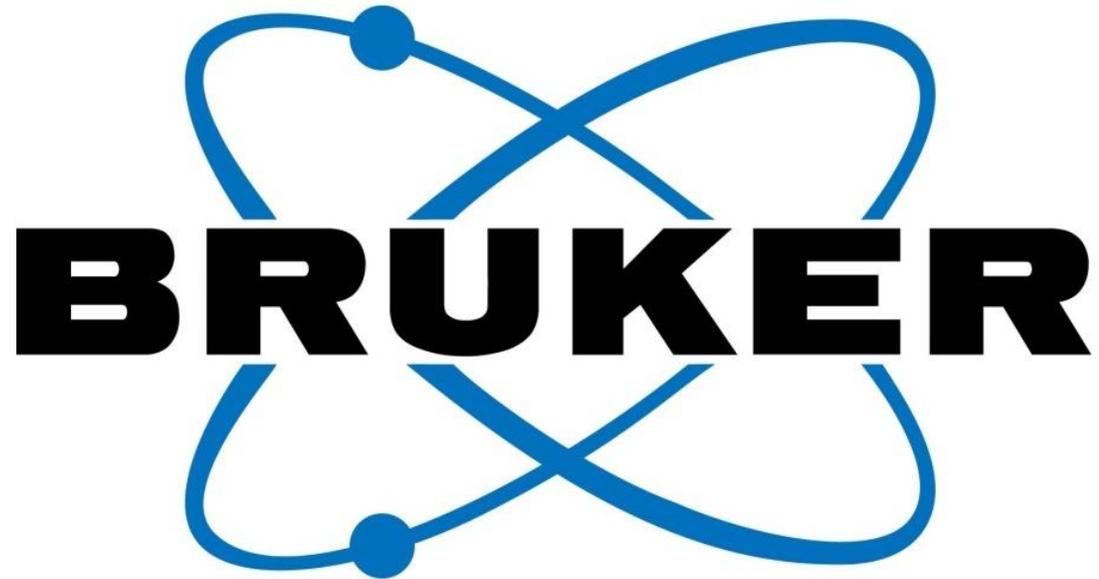

www.bruker.com



www.fischione.com