

Combination of targeted and non-targeted workflows for the identification of pollutants in river water using a passive sampling method



ASMS 2019 TP 199

Anthony Gravell¹; Melanie Schumacher¹; Bob Galvin²; Carsten Baessmann³
 Natural Resources Wales, Swansea University, Swansea, United Kingdom; ²Bruker UK Ltd., Coventry, United Kingdom; ³Bruker Daltonik GmbH, Bremen, Germany

Introduction

Many pollutants are ubiquitous in surface waters because of continuous discharges from municipal wastewater treatment plants and we still do not know which pollutants are reaching the environment, the size of the problem for exposed fauna, nor what effects, if any, of that exposure may be. Investigative monitoring of water bodies failing ecological standards as set out in the Water Framework Directive is now a requirement of all European Union member states. In this study we test the combination of a passive sampling device in combination with targeted and non-targeted workflows on a QTOF MS for monitoring surface water.

Methods

Chemcatcher[®] passive samplers were deployed for four weeks in the River Clun, Wales, UK, whose WFD status is classified as poor, to determine which chemicals may be responsible. The extracts obtained were analyzed using an impact II (Bruker) LC-QTOF-MS followed by targeted and non-targeted processing. The extracts were eluted using a 15 min gradient including a flow gradient (0.4-0.6 mL/min) on an UHPLC using a C18 (2.1 x 100 mm, 2.2 µm) column, with acidified water and methanol. For targeted screening and statistical analysis MS data were acquired using bbCID switching between MS and all ion fragmentation MS/MS data. For identification of unknowns some selected samples were analyzed using data dependent AutoMS/MS. Targeted data analysis for the identification of 'known unknowns' was performed with TASQ 2.1 using a combined database containing pesticides and drugs. Non-targeted data analysis for the identification of 'unknown unknowns' was performed with MetaboScape 4.0.

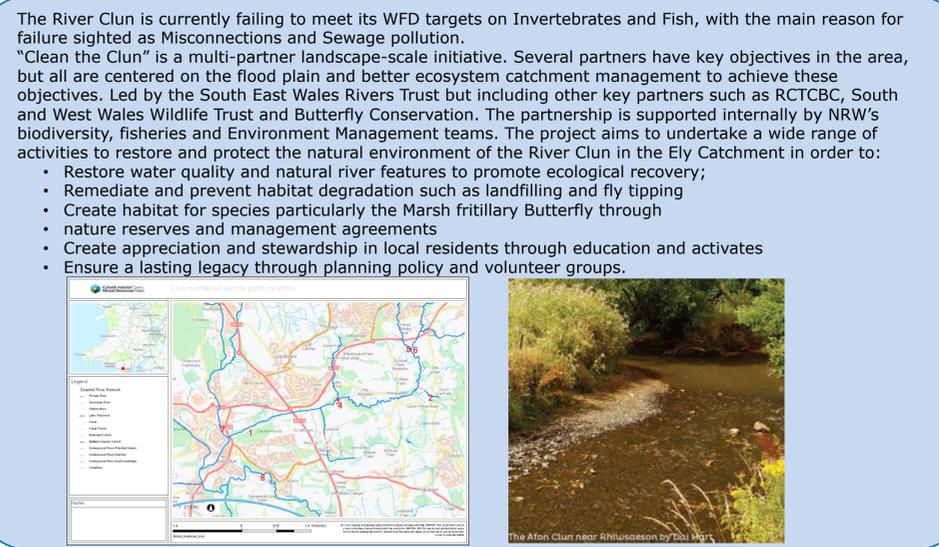


Fig. 1: Case study – Clean the Clun: Samples were taken over 10 months from 8 different sampling points using Chemcatcher passive samplers. The samplers stayed in the water for 4 weeks before analyzing the samples.

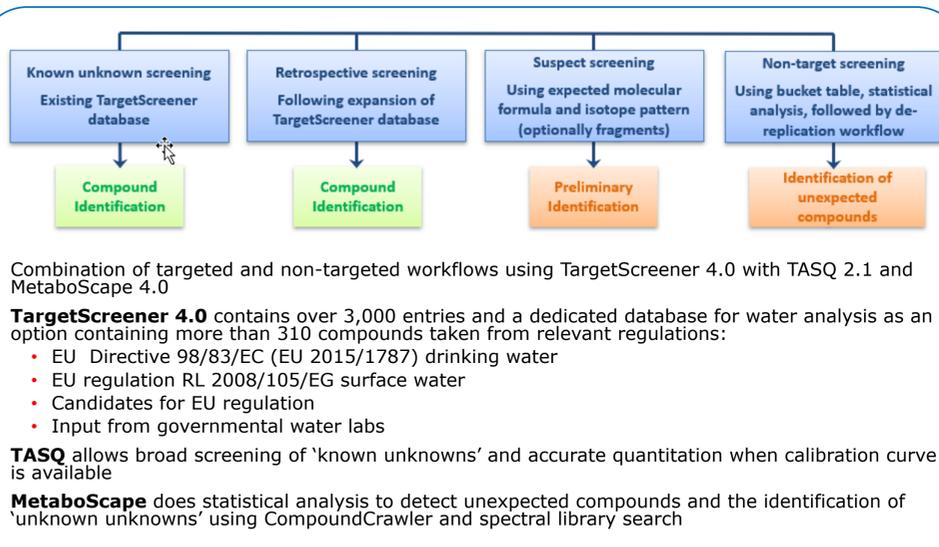


Fig. 2: Approach to screen for pollutants in the River Clun using targeted and non-targeted workflows.

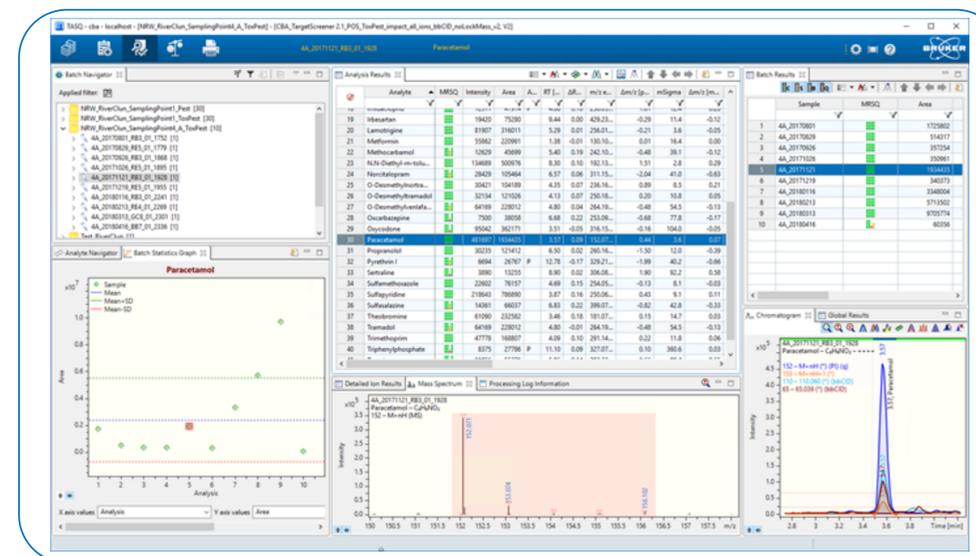


Fig. 3: TASQ analysis results from sampling point 4 and the sampling period 10/26/2017-11/21/2017. 6 pesticides and 36 drugs could be identified with a high MRSQ score and all mandatory qualifier ions found. The batch statistics shows a time course for 10 analyses from 7/4/2017-4/16/2018. Every data point represents a sampling duration of 4 weeks.

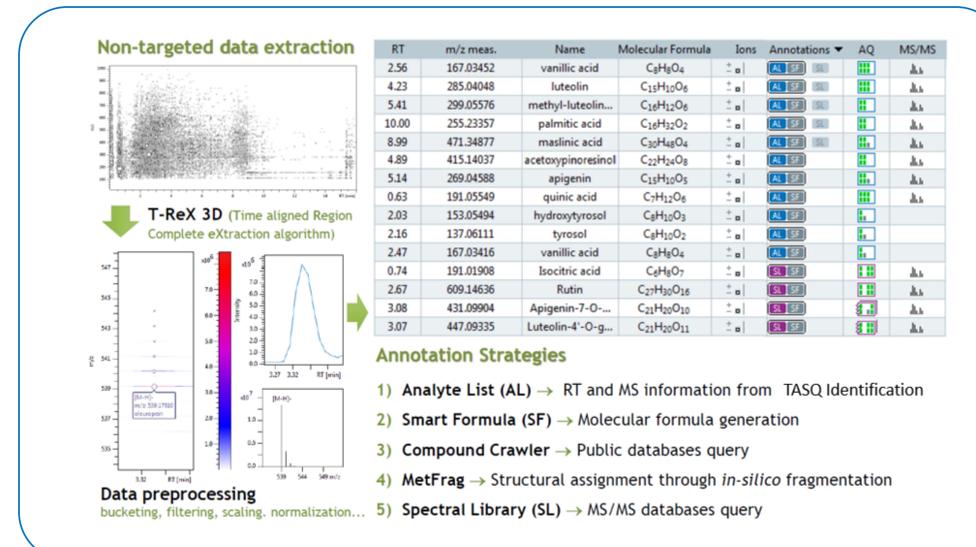


Fig. 4: Workflow for the identification of 'unknown unknowns' with MetaboScape

Results

Samples were acquired from July 2017 until April 2018 from 8 different places of the River Clun. For each site there was a time course of 10 data points covering 10 months. With the passive sampling unit every sample is an average of the pollution flowing past the sampling device during the 4 weeks it is submerged in the river. As expected, pollution in the river water due to discharges from municipal wastewater treatment plants varies significantly depending on the location and season. With the targeted approach between 5 and 93 compounds per sample could be identified with high confidence; in average about 50 compounds per sample. Most compounds were pharmaceuticals like Carbamazepine, Codeine, Citalopram, Paracetamol or Tramadol. For statistical analysis (PCA) and the identification of unknowns MetaboScape 4.0 (Bruker) is used. This is work in progress. We found that the passive sampling device can be used for monitoring pollutants in water and to obtain time-weighted average (TWA) or equilibrium concentrations of a wide range (2-3 orders) of pollutants in water. Most other sampling techniques involve the periodic collection of spot samples of water which is just snapshot of the conditions at the time of sampling. As pollutant concentrations fluctuate over time and may only be present at trace concentrations, the passive sampling device has many advantages compared to spot sampling techniques.

Conclusions

- The passive sampling device allows the end user to obtain a more representative picture of pollutants that may be present in the aquatic environment than other sampling techniques.
- TASQ allows rapid screening and quantitation of 'known unknowns' using databases with more than 3000 compounds.
- MetaboScape enables a de-replication workflow for the identification of 'unknown unknowns' using an Analyte List from TASQ, Smart Formula, Compound Crawler, MetFrag and Spectral Library Search.

QTOF / Environmental